

Help Manual

Help Manual	1
1 How to start?.....	1
1.1 Start	1
1.2 Create the camera video window.....	1
2 Video window GUI.....	2
2.1 Video window GUI	2
2.2 How to close the video window?.....	3
3 Image window GUI.....	4
3.1 Image window GUI	4
3.2 How to close the image window?.....	4
4 UI toolbar.....	7
5 Camera Sidebar.....	8
5.1 Capture & Resolution group.....	8
5.2 Exposure & Gain group	8
5.3 White Balance group	9
5.4 Color Adjustment group.....	9
5.5 Frame Rate group	10
5.6 Flip group	10
5.7 Color/Gray group	10
5.8 Power Frequency group	10
5.9 Sampling group	10
5.10 Histogram group	11
6 File.....	12
6.1 Open Image•••  Ctrl+O	12
6.2 Open Video••• 	14
6.3 Camera List	15
6.4 Twain: Select Device•••.....	16
6.5 Twain: Acquire•••	16
6.6 Save  Ctrl+S	18
6.7 Save As•••	18

Contents

6.8	Batch Save•••	22
6.9	Paste as New File	24
6.10	Print Setup•••	24
6.11	Print Preview••• Ctrl+Shift+P	24
6.12	Print••• Ctrl+P	24
6.13	Recent Files	25
6.14	Exit	25
7	Edit	26
7.1	Undo/Redo Ctrl+Z	26
7.2	Forward	27
7.3	Backward	28
7.4	Cut Ctrl+X	29
7.5	Copy Ctrl+C	29
7.6	Paste Ctrl+V	31
7.7	Image Select 	31
7.8	Select All Ctrl+A	32
7.8.1	Select all on the background layer	32
7.8.2	Select all objects over the background layer	32
7.8.3	Select all files in the browse window	33
7.9	Select None Ctrl+D	33
7.10	Inverse Selection	33
7.11	Delete File Delete	33
8	View	35
8.1	Browse  Ctrl+B	35
8.1.1	Open the browse window	35
8.1.2	Browse window right mouse button context menu	35
8.2	Sort>Sort by Names	36
8.3	Sort>Sort by Type	36
8.4	Sort>Sort by Size	36
8.5	Sort>Sort by Width	36
8.6	Sort>Sort by Height	36
8.7	Sort>Ascending	36
8.8	Sort>Descending	37

Contents

8.9	Icon>Large Icons	37
8.10	Icon>Small Icons	37
8.11	Refresh F5.....	37
8.12	Measurement Sheet	37
8.12.1	File Import•••	38
8.12.2	File Save•••	38
8.12.3	Export>To Html File	38
8.12.4	Export to Excel	39
8.12.5	Auto Highlight	39
8.12.6	Settings•••	39
8.13	Sidebar	40
8.13.1	Sidebar overview	40
8.13.2	Sidebar>Camera	41
8.13.3	Sidebar>Folders	41
8.13.4	Sidebar>Undo/Redo	42
8.13.5	Sidebar>Layer	42
8.13.6	Sidebar>Measurement	43
8.14	Ruler and Grid	43
8.14.1	Grids>No Grids	43
8.14.2	Grids>Manual Grids	43
8.14.3	Grids>Auto Grids	44
8.14.4	Grids>Remove All Grids.....	44
8.14.5	Settings•••	45
8.15	Best Fit	46
8.16	Actual Size	46
8.17	Track 	46
8.18	Properties	46
9	Setup.....	48
9.1	Start/Pause Pause	48
9.2	Full Screen ESC.....	48
9.3	View Properties•••	48
9.4	Video Overlay•••	49
9.4.1	Video Overlay: Overlay.....	49

Contents

9.4.2	Video Overlay: Marker•••	50
9.5	Video Watermark•••	51
9.6	Move Watermark	53
9.6.1	Move to•••	53
9.6.2	Move to zero	53
9.7	Rotate Watermark	53
9.7.1	Rotate to•••	53
9.7.2	Rotate to zero	54
9.8	Gray Calibration•••	54
9.9	Manual Fusion••• 	55
9.10	Video Source Property•••	56
9.11	Video Stream Format•••	56
9.12	Still Image Options•••	57
10	Capture	59
10.1	Capture Image F8	59
10.2	Time-lapse (Auto Capture)•••	59
10.3	Start Record••• F9	60
11	Image	64
11.1	Mode	64
11.1.1	Color Quantize•••	64
11.1.2	Gray Scale	64
11.2	Adjust	64
11.2.1	Curve•••	64
11.2.2	Auto Level	66
11.2.3	Auto Contrast	67
11.2.4	Histogram Equalization	68
11.2.5	Brightness/Contrast•••	68
11.2.6	Color•••	69
11.2.7	HMS•••	71
11.2.8	Gamma•••	71
11.2.9	Filter Color•••	71
11.2.10	Extract Color•••	72
11.2.11	Invert	72

Contents

11.3	Rotate	72
11.3.1	90 (CW)	72
11.3.2	180 (CW)	72
11.3.3	270 (CW)	72
11.3.4	Arbitrary***	73
11.3.5	Flip Horizontal	73
11.3.6	Flip Vertical	73
11.4	Crop Shift+C	73
11.5	Image Scale***	74
11.6	Histogram***	75
11.7	Resolution***	77
11.8	Overlay Scale Bar***	77
12	Process.....	79
12.1	Stitch	79
12.1.1	Video Stitch	79
12.1.2	Image Stitch	79
12.2	Filter*** Shift+F.....	81
12.2.1	Image Enhance	82
12.2.2	Edge Enhance	85
12.2.3	Morphological	86
12.2.4	Kernel	88
12.3	Range*** Shift+R.....	90
12.4	Segmentation*** Shift+S.....	91
12.5	Binary*** Shift+B.....	92
12.6	Emboss*** Shift+E.....	92
12.7	Pseudo Color***	93
12.8	Surface Plot***	93
12.9	Line Profile***	94
12.10	Diffuse*** Shift+D.....	95
12.11	Granulate*** Shift+G.....	96
12.12	Mosaic***	96
12.13	Fusion***	98
12.14	Color Composite***	99

Contents

13	Layer.....	106
13.1	About layer	106
13.2	Organizing layers	106
13.3	Layers for non-destructive measurement and label	106
13.4	Layer Sidebar	106
13.5	Layer menu and layer sidebar page context menu... ..	107
13.6	New•••	107
13.7	Remove•••	108
13.8	Current•••	108
13.9	Show/Hide•••	108
13.10	Rename•••	108
13.11	Export to Image	108
13.12	Export to Microsoft Excel.....	108
14	Measurements.....	109
14.1	Object Select 	109
14.2	Angle 	110
14.3	Point 	111
14.4	Line	111
14.4.1	Line>Arbitrary Line 	111
14.4.2	Line> Horizontal Line 	112
14.4.3	Line> Vertical Line 	112
14.5	Parallel 	113
14.6	Vertical 	114
14.6.1	Vertical>Four Points. 	114
14.6.2	Vertical>Three Points 	114
14.7	Rectangle 	115
14.8	RoundRect 	115
14.9	Ellipse 	116
14.10	Circle 	117
14.10.1	Circle>Center+Radius 	117
14.10.2	Circle>Two Points 	117
14.10.3	Circle>Three Points 	118
14.11	Annulus 	118

Contents

14.12	Two Circles 	119
14.12.1	Two Circle>Center+Radius.....	119
14.12.2	Two Circle>Three Points.....	119
14.13	Arc 	120
14.14	Text 	120
14.15	Polygon 	121
14.16	Z Order	121
15	Options.....	123
15.1	Preferences•••	123
15.1.1	File	123
15.1.2	Plugin	123
15.1.3	Print	124
15.1.4	Rulers and Grids	126
15.1.5	Cursor	127
15.1.6	Capture	128
15.1.7	Misc	128
15.2	Measurements•••	129
15.2.1	General>General	129
15.2.2	General>Length Unit	131
15.2.3	General>Angle Unit	133
15.2.4	General>Measurement Sheet.....	133
15.2.5	Object>Point	134
15.2.6	Object>Line	135
15.2.7	Object>Other Objects	136
15.2.8	Magnifications•••	136
15.3	Calibrate••• 	137
15.4	Edit Dye List•••	139
15.5	Auto Correction•••	141
16	Windows.....	142
16.1	Close All	142
16.2	Windows•••	143
17	Help.....	144
17.1	Help Contents	144

Contents

17.2

About 144

1 How to start?

1.1 Start

1. Double click on the desktop icon “” to start ToupView;
2. Click the **Start** button (At your screen bottom left corner) and a **Start** menu will bring up. Move your mouse over the menu and try to locate ToupView, click to start it.

1.2 Create the camera video window

ToupView will detect all of the cameras that your computer has installed (Here, it is **UCMOS03100KPA**, a 3.1M pixel CMOS camera) and will append all the camera names as submenu to the **File>Camera List** menu (Here, the submenu name is “**UCMOS03100KPA**”).


Choosing the **File>Camera List>UCMOS03100KPA** will create a video window and start the video stream. The video window will be associated with the name of “**Video [UCMOS03100KPA]**” (i.e., its title bar name will be “**Video [UCMOS03100KPA]**”).

There are 3 methods to start the video stream, they are:

File>Camera List	Choose the File>Camera List>UCMOS03100KPA (Here, a 3.1M pixel camera is installed) command to create the video window;
Camera Sidebar	Click the Camera Sidebar (If it is not activated) and the Camera List to expand the Camera List group (if not expanded). Click the camera name (Here it is UCMOS03100KPA) to create the video window; 
Toolbar button	Click the button  ’s down arrow on the toolbar to expand the camera list and choose the right camera (Here it is UCMOS03100KPA) to create the video window. 

2 Video window GUI

2.1 Video window GUI

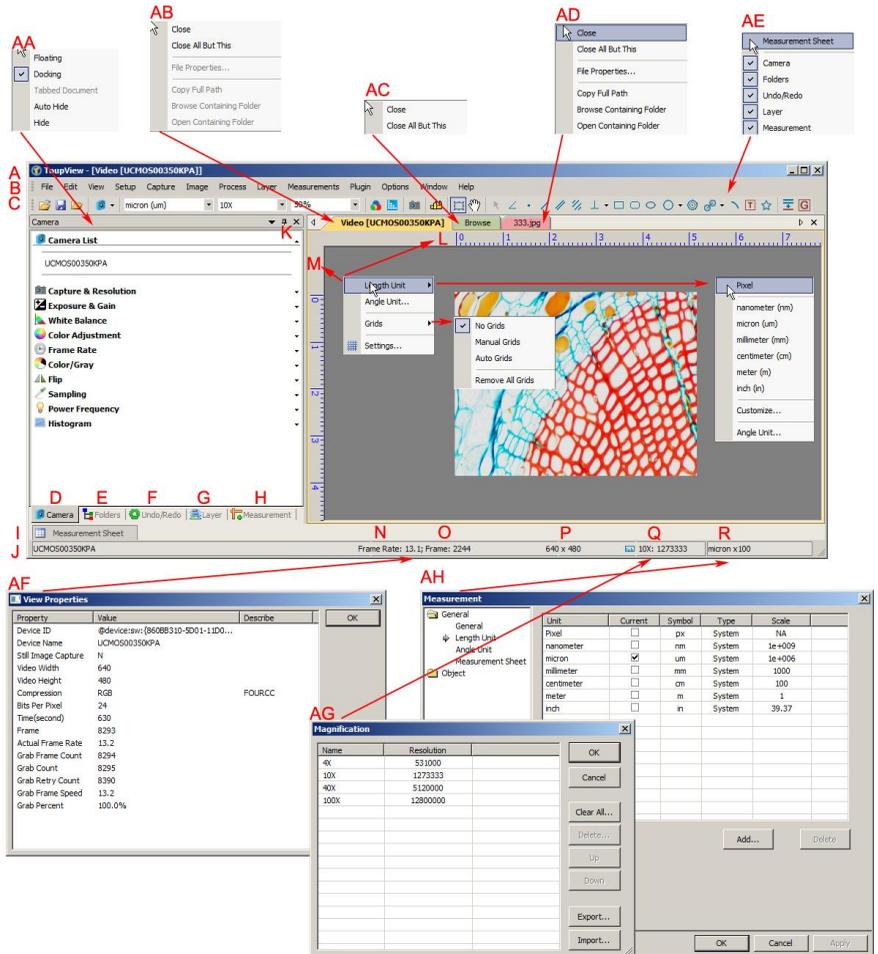

A:ToupView; B: Menu; C:ToupView toolbar D:Camera Sidebar ;
 E: Folders Sidebar ; F: Undo/Redo Sidebar ; G: Layer Sidebar ;
 H: Measurement Sidebar ; I: Measurement Sheet; J:Statusbar;
 K: Auto Hide button L: Horizontal ruler; M:Vertical ruler
 N:Frame Rate O:Frames captured P:Current Video sizes


- Q: Selected microscope **Magnification** R: Current **Unit**;
- AA: **Sidebar** right mouse button context menu;
- AB: **Video** window right mouse button context menu;
- AC: **Browse** window right mouse button context menu;
- AD: **Image** window right mouse button context menu;
- AE: **Frame** window right mouse button context menu;
- AF: Double-click bring up **Video Properties** dialog;
- AG: Double-click bring up **Magnification** dialog;
- AH: Double-click bring up **Measurement** dialog;
- AI: **Horizontal Ruler** or **Vertical Ruler** right mouse button context menu

2.2 How to close the video window?

- 1 Double-clicking the tabbed video window title or clicking **x** on  will close the video window directly;
- 2 For MDI video window, click **x** on  to close the video window;
3. Choose the **Windows>Close All** command to close the video window.

3 Image window GUI

3.1 Image window GUI

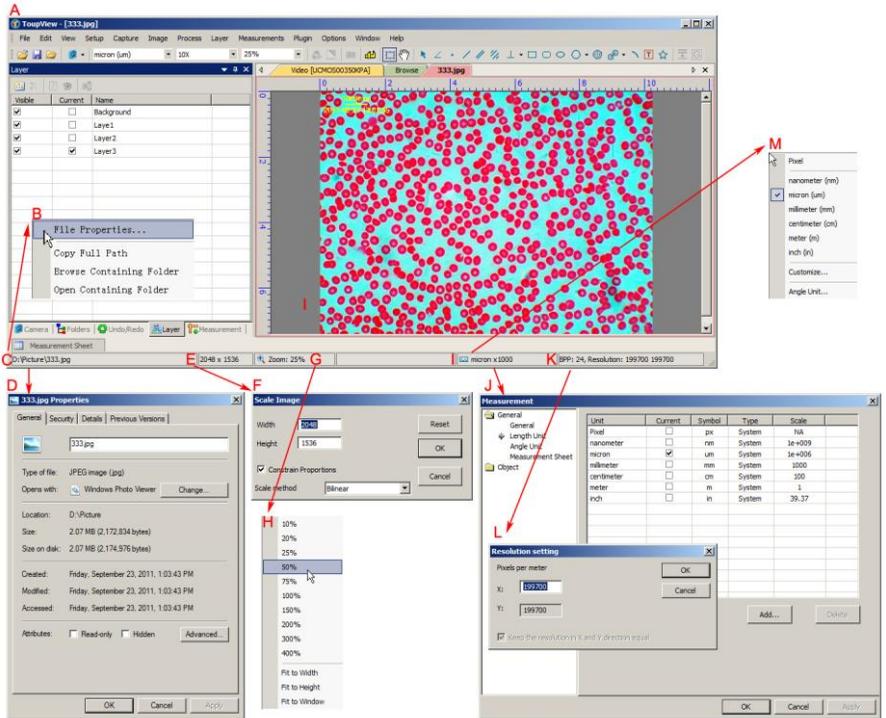

A:TopView; B:The opened file right mouse button context menu on status bar;

C:The opened file name and directory;

D:Double-click bring up opened file **Properties** dialog;

E:Current image width and height;

F:Double-click bring up **Scale Image** dialog;

G:Image **Zoom** ratio, double-clicking will zoom the image to 100%;

H:**Zoom** ratio right mouse button context menu;

I:Currently selected **Unit** J: Double-click bring up **Measurement** dialog;

K:Image **BPP & Resolution**;

L:Double-click bring up **Resolution Setting** dialog;


M: **Unit** right mouse button context menu.

3.2 How to close the image window?

Help Manual


1. Tabbed window

a) If you have modified an image before attempting to close it, double-clicking the tabbed image window title or clicking **x** on  will bring up a **ToupView** dialog:

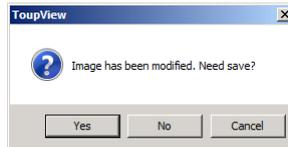


Clicking **Yes** will save the changes with its old name and close the window quickly, **No** will close the file immediately with no changes and no warning, or **Cancel** will cancel the **Close** command and leave the window there with no changes;

b) If the **Image** window is snapped from the video window and with number as its title, double-clicking the tabbed image window title or clicking **x** on  will bring up a warning **ToupView** dialog:

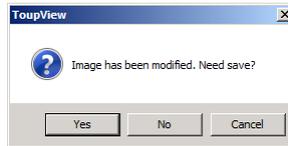

Clicking **Yes** will bring up the **Save As** dialog:

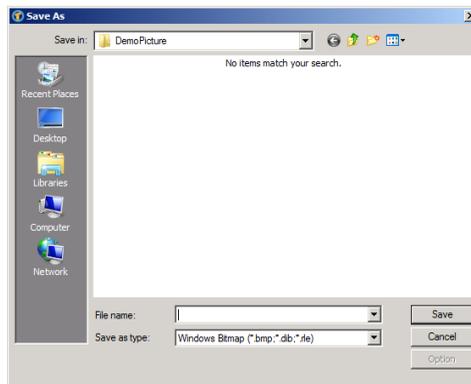

Select the driver and folder to which you want your image file saved in the **Save in** combobox and enter the file name in the **File name** edit box.

Click **Save** to save the captured image with the specified directory and file name, or **Cancel** to close the **Save As** dialog and return to image window.

Clicking **No** on the **ToupView** dialog will close the file immediately with no changes and no warning or **Cancel** on the **ToupView** dialog will cancel the **Close** command and return to image window.

Note: Choosing the **Windows>Close All** command can also close the tabbed image window. Check the **Windows>Close All** for detail.

2. MDI window

a) For an opened and modified image window, click **X** on  to bring up the same **ToupView** dialog. The next operations are just the same as the tabbed window in step 1a;

b). For a **Video** captured image with number as its title, clicking **X** on  will bring up the same **ToupView** dialog. The next operations are just the same as step 1b in the tabbed window

Note: Choosing the **Windows>Close All** command will also close the MDI image windows in a quick way. Check the **Windows>Close All** for detail.

4 UI toolbar

When the camera is started or the image is opened, most of the icons on the toolbar will be available for the quick setup of the video or image characteristic.



- 1: **Open** (Ctrl + O)
- 2: **Camera List**
- 3: **Save** (Ctrl + S)
- 4: **Unit**
- 5: **Magnification**
- 6: **Zoom**
- 7: **Video Source Properties**(will be enabled only for the camera that support directshow interface)
- 8: **Video Stream Format**(will be enabled only for the camera that support directshow interface)
- 9: **Video/Image Select**
- 10: **Track**(enabled only when the video/image sizes large than the window size)
- 11: **Object Select**(will be enabled only when an object is existed on the layer above the background layer)
- 12: **Angle**
- 13: **Point**
- 14: **Line**
- 15: **Parallel**
- 16: **Two Parallel**
- 17: **Vertical**
- 18: **Rectangle**
- 19: **RoundRect**
- 20: **Ellipse**
- 21: **Circle**
- 22: **Annulus**
- 23: **Two Circles**
- 24: **Arc**
- 25: **Text**
- 26: **Polygon**
- 27: **Calibration** (for both video/image window)
- 28: **Gray Calibration** (will be enabled only when a rectangle video area is selected)
- 29: **Manual Fusion**.
- 30: **Manual Fusion**.

5 Camera Sidebar

The **Camera Sidebar** is used for the control of ToupCam camera, it includes 11 groups. The group can be expanded by clicking the group name or clicking the **Down Arrow** at the right of the group name.



The down arrow to expanded the group



5.1 Capture & Resolution group



Snap: Continuously Snap images by clicking it;
Record: Record video stream in wmv/asf or avi format;
Live: Set the video resolution;

Snap: Set the Snap resolution for the image capture process.

5.2 Exposure & Gain group

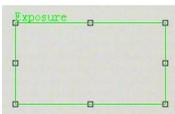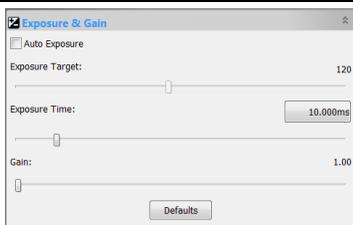

Rectangle for Auto Exposure


1. When the **Exposure & Gain** group is expanded, a green rectangle marked with "Exposure" will be overlaid on the video. This marked region is a reference region for judging if the image brightness reach to the **Exposure Target** value. Drag or resize the **Exposure ROI** to the dark area will increase the video brightness and drag it to the brighter area will decrease the video brightness;
2. Uncheck the **Auto Exposure** box to switch the **Auto Exposure** mode to the **Manual Exposure**.

Help Manual

The **Exposure Target** slider will be disabled in this mode;

3. Tune the microscope light source to a bright state, and then drag the **Exposure Time** slider left or right until the image brightness is normal;
4. If and only if the microscope light intensity is too low to meet the **Exposure Time** up limit, drag the **Gain** slide right until the video brightness is normal;
5. The exact **Exposure Time** can also be entered by clicking the edit box at the right of the **Exposure Time**:. This will bring up a dialog called **Exposure Time**. You can type the number in the field to set the exact **Exposure Time** value.

5.3 White Balance group

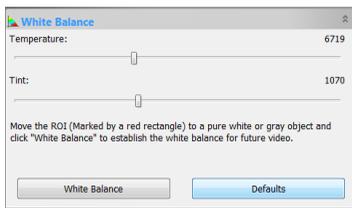


1. Click the **White Balance** bar to expand the **White Balance** group and a red rectangular marked with **White Balance** will be overlaid on the video window;
2. Drag or resize the red rectangle to a pure white or gray area and click the **White Balance** button to establish the video **White Balance** for

future video streaming process;

3. If the automatic setting and the actual result still has deviation, drag the **Temperature** and **Tint** slides to left or right to manually correct the **White Balance**.

5.4 Color Adjustment group

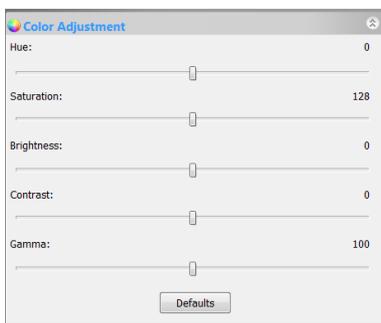

1. **Hue**: Adjusts the **Hue** of the video. Drag the slider to the right to increase or drag to the left to decrease hue;
2. **Saturation**: Adjusts the **Saturation** of the video. Drag the slider to the right to increase or drag to the left to decrease video saturation;
3. **Brightness**: Adjusts the video **Brightness**. Drag the slider to the right to increase or drag to the left to decrease the video brightness;


4. **Contrast**: Adjusts the video **Contrast**. Drag the slider to the right to increase or drag to the left to decrease the video contrast;

5. **Gamma**: Adjusts the video **Gamma**. Drag the slider to the right to increase or drag to the left to decrease the video gamma;

Help Manual

6. **Defaults**: Click the **Defaults** to clear all the changes and reset them to default ones;
7. All of your settings will be saved for future **Color Adjustment** application.

5.5 Frame Rate group

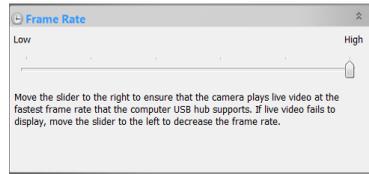

Drag the slider to the right (**High**) to ensure that the camera can run the video at the fastest **Frame Rate** that your computer USB hub supports. If the video fails to display, drag the slider to the left (**Low**) to reduce the **Frame Rate** and enable the video in a low speed mode.

5.6 Flip group



to set the video direction to the right one.

If the video on the screen appears in different directions from what is viewed under the camera, check the "**Horizontal**" or "**Vertical**" box

5.7 Color/Gray group



If you wish to preview **Color** video, select the "**Color**" button;

If you wish to preview **Gray** video, check the "**Gray**" button.


5.8 Power Frequency group



1. A CMOS sensor captures each row of pixels (from top to bottom) in sequential order, creating a rolling effect, hence the name "Rolling Shutter". Instead of being relatively constant, so for example, as the commercial main frequency in Europe is 50Hz, fluorescent lights in Europe flicker at 100 times per second and as the main frequency in US is 60Hz, so in the USA they flicker at 120 times per second;
2. This flickering problem is solved by capture row pixels in over the duration of integer number of (n) flicker periods;
3. check **50HZ** will delete the rolling dark band for the **50HZ** light fluctuation;
4. check **60HZ** will delete the rolling dark band for the **60HZ** light fluctuation;
5. For **DC** power, no light fluctuation is existed and no compensation is needed.

5.9 Sampling group


Help Manual

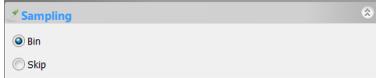

1. **Bin**: Pixel binning refers to the method of combining (averaging) pixels of blocks of neighboring same color pixels to resize the video the lower resolution;

2. **Skip**: Also called "Decimation", means that a certain amount of pixels are not read out but skipped (horizontally, vertically or in both axes). This reduces resolution of the resulting video but introduces subsampling artifacts.

5.10 Histogram group

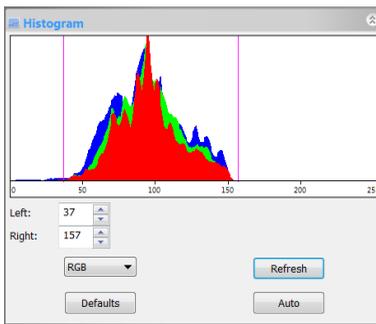

1. A **Histogram** illustrates how pixels in an image are distributed by graphing the number of pixels at each color intensity level. The **Histogram** shows detail in the shadows (shown in the left part of the histogram), midtones (shown in the middle), and highlights (shown in the right part). A **Histogram** can help you determine whether an image has enough detail to make a good correction;

2. This group shows the **Histogram** of current active video. Two vertical line markers show the upper and lower limits of the intensity levels. These markers can be dragged with mouse. If you are looking at a color image, the **Histogram** will reflect the **RGB**(red, green and blue channels histogram at the same time) **R**(red), **G**(green), and **B**(blue) values with lines of the same color;

3. You can also enter directly the desired values in the **Left** or **Right** boxes below the **Histogram** chart for both **Left** and **Right Histogram** boundaries;

4. Click the "Refresh" button to update the **Histogram** display if the sample under observation is moved or changed;

5. Clicking the **Defaults** button will return the **Left** and **Right Histogram** boundaries to its original ones;


6. Click the **Auto** button to locate the two boundaries automatically to get the best video quality.

6 File

6.1 Open Image●●● Ctrl+O

Choose the **File>Open Image●●●**  command to open an existing image file. The **Open Image●●●** command can also be used to preview an image in small size, or to view its statistics and information without actually opening the image itself. These capabilities can be used to quickly locate a particular image.

ToupView supports and can open many image formats. These are identified in the **Files of type** list box. You may also open an image file type with **ToupView File Type (*.tft)** format which has **Measurement Objects** overlaid on the image.

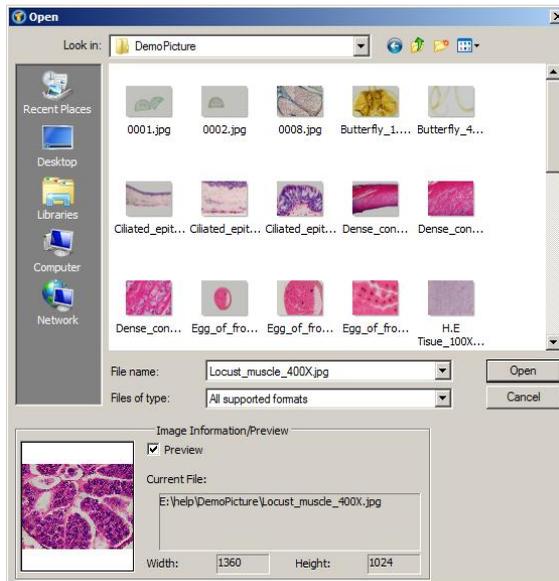

More than one image can be opened within ToupView simultaneously by a) with Ctrl + left mouse button method and b) Shift + left mouse button method to highlight the files to be opened;

When opening an image, ToupView places it into a new image window. It then becomes the active image.

Note: ToupView maintains, at the **File>Recent Files** submenu, a list of the last 4(Can be 1-8) opened files. Any of these files can be accessed by simply clicking on file

Help Manual

name. If no files are listed in the **File>Recent Files** submenu, the **Open Image...** command must be used to open the file.

Also, the **View>Browse** can be used to view images in an **Icon** format under any selected directory. Brief information is given in the **View>Browse** menu.

File name	<p>From this list box, select the name of the file want to open. Either the type of the file name (with its entire path, if it is not in the current folder), or selecting Files of type to obtain a list of file names. Double-clicking a file name in the large combobox (where both folder and file names are listed) will automatically open it.</p> <p>Note: If just type in the file name, be sure that the Files of type field correctly identify the format of the file to open. Otherwise error messages will bring up when ToupView tries to open the file.</p>
Files of type	<p>In this list box, select the image format of the file to open. If one selects All supported formats, ToupView uses the file's extension to identify its format. ToupView supports the following file formats:</p> <ul style="list-style-type: none">Window Bitmap(*.bmp,*.dib,*.rie)JPEG(*.jpg,*.jpeg,*.jpe,*.jif,*.jfif)Portable Network Graphics(*.png)Tag Image File Format(*.tif, *.tiff)Compuserve GIF (*.gif)Targa(*.tga)PhotoShop(*.psd)ICON(*.ico)Enhanced Window Metafile(*.emf)Window Metafile(*.wmf)JBIG(*.jbg)Wireless Bitmap(*.wbmp)ToupView File Type(*.tft) <p>If the image file does not use standard format-identifying extensions, the file in the File name field must be typed, and then select its format from the Files of type list box. Otherwise, ToupView will select a format based on the file name extension.</p>

Help Manual


Preview	Click this button to preview image in small size. In preview mode, statistics and information about the image (i.e. image Width , Height and image location) will be displayed. The default is no Preview .
Current File	Current File location on your computer
Width	Current File image width;
Height	Current File image height;

6.2 Open Video●●●



The **File>Open Video●●●** menu will be enabled only when there is no video file opened in the ToupView frame or no camera start up.

1. Choose the **File>Open Video●●●** command to open an existing video file;

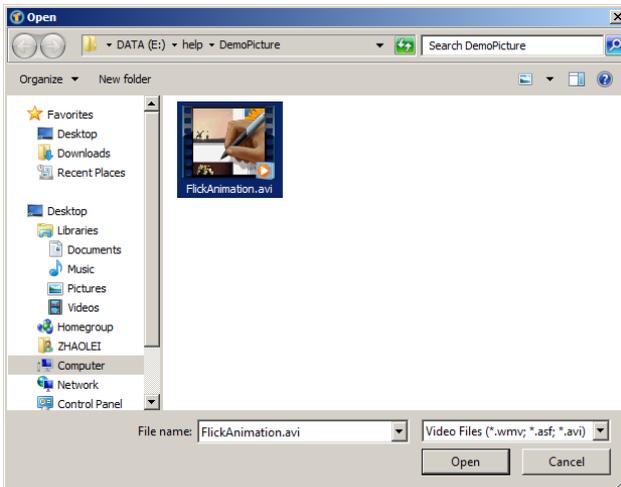

2. Select the name of the file you want to open. If the file does not appear, select the option for showing all files from the **Files of Type** combobox. The video file type can be ***.wmv***; ***.asf*** or ***.avi*** format.

3. Click **Open** to open a video file, this will create a video window and begin to start the video stream. The video window will be associated a name called **"Video [XXX.XXX]"** (i.e., its title bar will display **"Video [XXX.XXX]"**, here, **XXX.XXX** is the video file name).

4. Click **Cancel** to return to the application.

Note: Only a single video can be opened at a time. ToupView takes camera as an special video file, if the camera is started, this menu will be disabled and the video file cannot be opened anymore.

6.3 Camera List

The **File> Camera List** submenu will be disabled when there is a video file opened in the ToupView frame or a camera start up.

ToupView will detect all of the cameras that your computer has installed (Here, it is **UCMOS03100KPA**, a 3.1M pixel CMOS camera) and will append all the camera names as submenu to the **File>Camera List** menu (Here, the submenu name is "**UCMOS03100KPA**").

Choosing the **File>Camera List>UCMOS03100KPA** will create a video window and start the video stream. The video window will be associated with the name of "**Video [UCMOS03100KPA]**" (i.e., its title bar name will be "**Video [UCMOS03100KPA]**").

There are 3 methods to start the video stream, they are:

File>Camera List	Choose the File>Camera List>UCMOS03100KPA (Here, a 3.1M pixel camera is installed) command to create the video window;
Camera Sidebar	Click the Camera Sidebar (If it is not activated) and the Camera List group to expand it (if not expanded). Click the camera name (Here it is UCMOS03100KPA) to create the video window; 
Toolbar button	Click the Camera List button  's down arrow on the toolbar to expand the camera list and choose the right camera (Here it is UCMOS03100KPA) to create the video window. 

Note: a) Only a single video can be opened at a time. ToupView takes camera as a special video file, if the camera video window is opened, the **File>Open Video...** menu will be disabled and the video file cannot be opened anymore; b) If a video file is opened, the camera cannot be started.


6.4 Twain: Select Device●●●

Twain is a cross-platform interface for acquiring images captured by certain scanners, digital cameras, or frame grabbers. The manufacturer of the **Twain Device** must provide a **Source Manager** and **Twain Data Source** to work with ToupView.

Select the active device for the **Twain: Acquire●●●** menu from all devices available in the device combobox which are enumerated by the application.

One must install the **Twain Device** hardware and its driver first. See the documentations provided by the device manufacturer for the installation instructions.

Before begin to start the **Twain: Acquire** at the first time with ToupView, choose the **File>Twain: Select Device●●●** command first, this will invoke a dialog called Select Source as shown below:

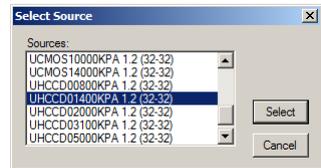

Source: Choose the right device from the **Source** window (Highlight it)

Select: Click **Select** to select the device. User does not need to repeat this step for subsequent choosing of the **Twain: Acquire●●●** command.

6.5 Twain: Acquire●●●

Introduction

There are basically two techniques used to capture the video images from video devices such as a PC camera, digital camera, and scanner. They are the **Twain: Acquire●●●** technique and the **DirectShow** technique (previously called **VFW**).


The most obvious characteristics of the **Twain** technique is that it previews the video in smaller resolution but captures the image with higher resolution. The **USCMOS** and **UHCCD** series cameras support all of these two video preview techniques.

Steps for Twain Acquire

Here we illustrate how to capture the image using a **UHCCD01400KPA** (1.4M pixels, USB2.0) camera as an example.

Help Manual

1. Install the camera driver (for example driver for **UHCCD01400KPA** hardware);
2. Install ToupView;
3. Plug the cameras **UHCCD01400KPA** (USB2.0) into the computer;
4. Start ToupView;
5. Choose the **File>Twain: Select Device...** command to select the device **UHCCD01400KPA** from the **Select Source** dialog (If never selected before).
6. Choose the **File>Twain: Acquire...** command. There should be a dialog box like below:

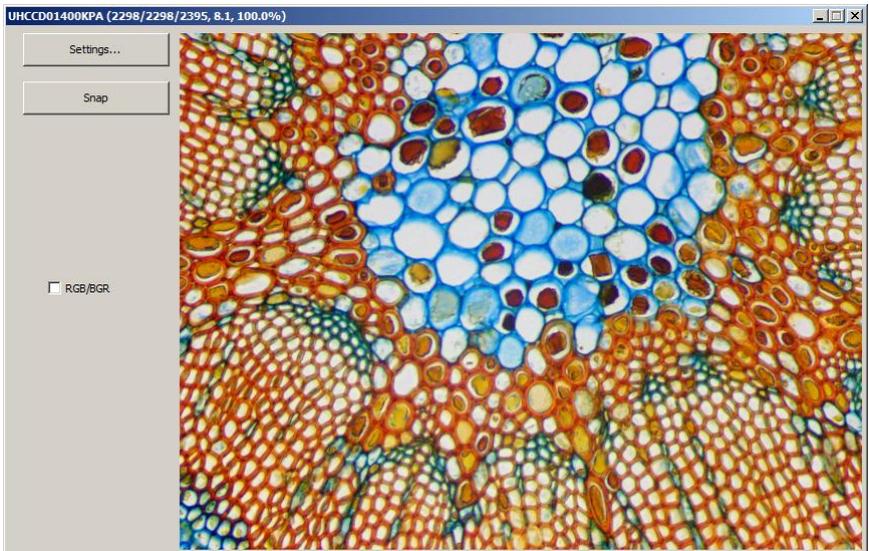

In this dialog, video **Resolution** can be selected (Combobox if it has). The **Video Source Property** can be set by clicking the **Setting...** button. Click the **Snap** button to capture an image. This will create a new window and its title bar will be assigned a digital as the image window name;

Check the **RGB/BGR** box to ensure the correct color encoding format compatible with your application.

Click **X** on  to close the **Twain: Acquire** dialog.

6.6 Save **Ctrl+S**

Choose the **File>Save**  command to immediately store the current window image to its file (the file name is listed on the window's title bar) while leaving the image still active in its window.

If the image is untitled or titled with a digit, ToupView will issue the **File>Save As...** dialog automatically. The default "Save as type" will be "Window Bitmap (*.bmp,*.dib,*.rle)". If there are **Measurement Objects** overlaid on the image, the default **Save as type** will be "ToupView File Type (*.tft)".

The **File>Save** command can be used to save the most recent changes to disk. It is often performed as a precautionary measure during lengthy or involved processes to reduce the amount of reprocessing that might be required in the event of a system failure or operational error. When an image is closed and not to save its changes is chosen, ToupView discards all changes made since the last **File>Save** operation.

Note: a).The **File>Save** command always saves the contents of the entire window, even if there is an **AOI (Area of Interest)** defined on it; b).The **File>Save** command will be disabled if the file is not changed or the changes have been saved.

6.7 Save As...

Choose the **File>Save As...** command to store the contents of the current window to a specified file format. At the end of a **File>Save As...** operation, the image window will be associated with the new file and the new format (i.e., its title bar will display the new file name). ToupView supported file save formats are:

Window Bitmap (*.bmp,*.dib,*.rle)

JPEG (*.jpg,*.jpeg,*.jpe,*.jif,*.jfif)

Portable Network Graphics (*.png)

Tag Image File Format (*.tif, *.tiff)

Compuserve GIF (*.gif)

PCX (*.pcx)


Targa (*.tga)

JBIG (*.jbg)

ToupView File Type(*.tft)

Save in: Find the folder where the file wishes to be saved. A new folder may be created using the **New Folders** button.

File name: Enter to be saved file name. To specify the file's location, either enter its entire path (disk and folder), or specify its location using the **Save in** list box.

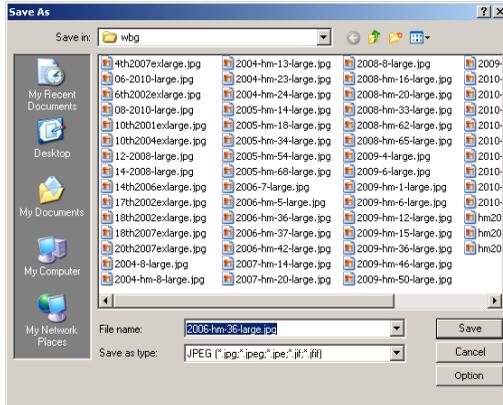


Save as type: In this list box, select the format in which the image wants to be saved. **Save As** is also used to convert a single image from one format to another. For example, if a **TIFF** file needs to convert to **PCX** format, open the **TIFF** image first, then choose **Save As** command with the **PCX** format option to save it to a new file.

The default "**Save as type**" will be "**Window Bitmap (*.bmp;*.dib;*.rle)**". If there are **Measurement Objects** overlaid on the image, the default **Save as type** will be "**ToupView File Type (*.tft)**".

The **Save As** command has several important uses beyond simply storing an image to a new file name. Click **Option** to select the different parameters to encode the file.


Help Manual

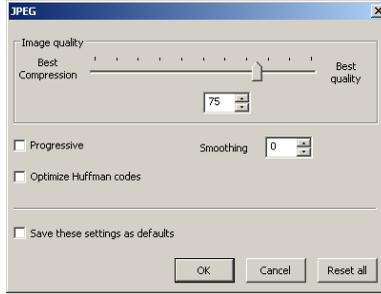

For JPEG (*.jpg,*.jpeg,*.jpe,*.jif,*.jfif), Option has the following items:

Image quality	If one save an image in JPEG format (*.jpg), one may adjust image quality in the edit box. The values range from 0 to 100. Default value: 75.
Progressive	The default is unchecked.
Optimize Huffman codes	The default is unchecked.
Smoothing	The values range between 0 and 100. Default value: 0.
Save these setting as defaults	When saving a file, the current settings will be saved as defaults for the next file save operation.


For Portable Network Graphics (*.png), Option has the following items:

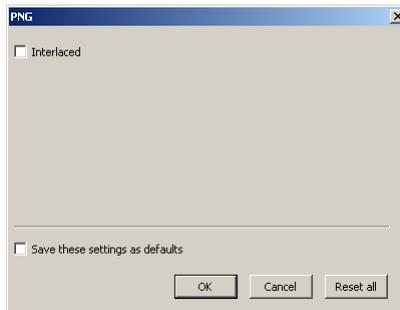

Interlaced	The default is unchecked.
Save these setting as defaults	When saving a file, the current settings will be saved as defaults for the next file save operation.

For Tag Image File Format (*.tif, *.tiff), Option has the following items:


Help Manual

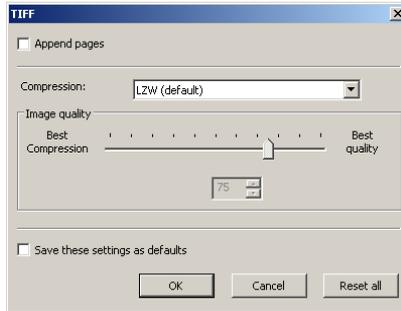

Appended pages	Determine whether the current image will be saved in multiple pages style or not.
Compressions	Specifies a method for compressing the composite image data. For saving a 32-bit TIFF file, one can specify that the file be saved with predictor compression, but have no option to use JPEG compression. Predictor compression offers improved compression by rearranging floating point values, and works with both LZW and ZIP compression.
Image quality	If choosing Compressions as "JPEG", the Image quality can be adjusted by the slider bar. The values range between 0 and 100. Default value: 75.
Save these setting as defaults	When saving a file, the current settings will be saved as defaults for the next file save operation.

For **Compuserve GIF (*.gif)**; **PCX(*.pcx)**; **Targa(*.tga)**; **JBIG(*.jbg)**; **ToupView File Type(*.tft)**, There is no **Option**.

Note: a) Detailed information of the above academic terminologies can be found in books about image processing and image compression or internet; b) The file saved directory can be restored for future application. To keep the directory unchanged when ToupView is started again, choose **Options>Preferences...** command, click **Misc** page and check **Restore the current directory when startup** under the **Privacy** item.

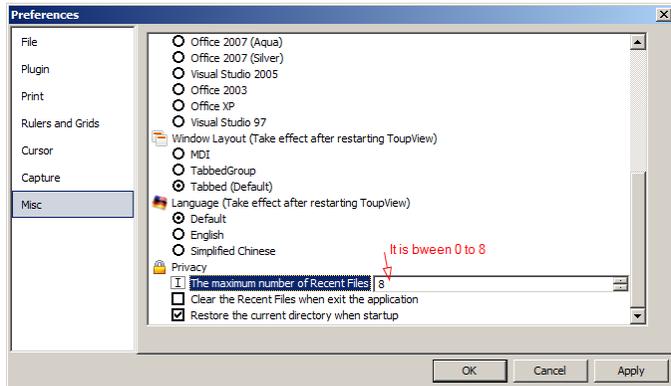

6.8 Batch Save●●●

The **File>Batch Save●●●** menu will be enabled when a) an image is opened; b) an image is captured from the camera; c) an image window is created by choosing the **File>Paste as New File** menu; d) an image window is created copied from the **Undo/Redo Sidebar**.

Batch Save for the a) style image window

1. If the image is modified, choosing the **File>Batch Save●●●** command will save the file with its opened file name and file extension;
2. If the image is modified because of the **Measurement Objects**, choosing the **File>Batch Save●●●** menu will save the file with its opened file name but with "ToupView File Type (*.tft)" file extension format;


Batch Save for the b), c) or d) style image window

If the image window has been created with b), c) or d) style, choosing the **File>Save As●●●** command to realize the file saving task will be time-consuming. The **Batch Save●●●** command will runs **File>Save As●●●** command with the name automatically specified according to the paradigm specified in the **Batch Save** dialog

To start the **File>Batch Save●●●** command, you have to

1. Start the camera;
2. Snap at least an image first;

3. Choosing the **File>Batch Save...** command will bring up a **Batch Save** dialog:

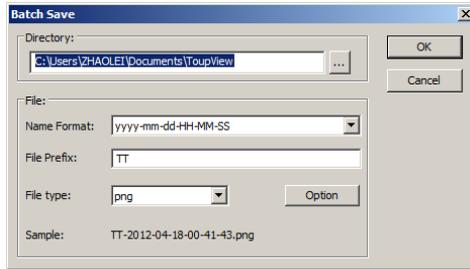

Directory: Enter the name of the drive and directory where your captured images will be saved. You may either type the path information, or use the **Browse** button to locate it from a standard **Browse Folder** dialog.

Name Format: The **year**, **month**, **date**, **hour**, **minute** and **second** or **nnnn(sequence)** are used to as part of the file name. If more files are saved with in a second, a **(xx)** suffix is attached to the end of **Name Format** to avoid the possible name conflict. For the **nnnn(sequence)** "**Name Format**", no suffix is needed.

File Prefix: Enter a file name prefix for **Batch Save** when generating files names for a series of images. This prefix will be combined with **Name Format** to form a final file name naming paradigm.

File Type: In this combobox, select the format in which you want the image to be saved (can be **BMP**, **JPG**, **PNG**, **TIF**). Click **Option** button to set the different parameters for encoding the file (For **BMP format**, the **Option** will be disabled. See **File>Save As...** menu about the details of the format encoding methods); If there are Measurement Objects over the image, the **File Type** will be chose as "**ToupView File Type (*.tft)**" by ToupView.

Sample: The final file name is shown at the right of the **Sample** label for reference.

4, If everything was finished, click **OK** button to begin the file batch save process or **Cancel** to cancel the **File>Batch Save...** command and return to the application.

Note: a) In the process of the **File>Batch Save...** command, the title on the image tab or image window will be modified with the file name formed in the **Batch Save** dialog. b) The **File>Batch Save...** will perform no saving operation if the file is not

modified.

6.9 Paste as New File

The **File>Paste as New File** menu will be enabled when there is valid image data on the clipboard first (see the **Edit>Copy** menu). If there is no image data on the clipboard, the **File>Paste as New File** menu will be disabled.

Choose the **File>Paste as New File** command to place the contents of the clipboard image into a new image window, which becomes the active image.

The new image type will be the same as that of the original image. ToupView will accept image data from other applications via the clipboard as long as it is in **Windows Bitmap (DIB)** format.

Note: ToupView will assign a digit to the **Paste as New File** command created image window title bar.

6.10 Print Setup●●●

Choose the **File>Print Setup●●●** this command to access the setup panel for the printer that has selected. ToupView will present the standard setup panel for the particular printer (this is the same panel one would receive if one were setting up the printer from the **Windows Control Panel**). Change printer's setup to satisfy the requirements, click **OK** button to return.

6.11 Print Preview●●● Ctrl+Shift+P

Choose the **Print Preview** command to see the real-time effect of the printer without actually printing it out.


6.12 Print●●● Ctrl+P

Choose the **File>Print●●●** command to print one or more copies of the current image to the selected output device. The ToupView **File>Print●●●** command lets one take full advantage of the printer's capabilities. If the printer has built-in half-toning or color dithering capabilities, use them or instruct ToupView to perform these processes before sending the image to the device.

The **File>Print●●●** command also has facilities that let one adjust the size and position of the image on the printed page.

6.13 Recent Files

ToupView maintains 4 (default) most recently opened document files under the **Recent Files** menu. Choosing one of these submenus will reopen that file immediately.

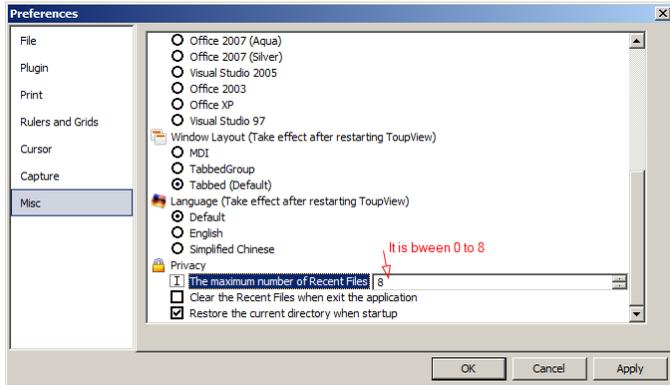

1. The maximum number of Recent Files can be modified by choosing the **Options>Preferences...** command and clicking the **Misc** page. Here, clicking the **4** (default) edit box will allow you to enter the number of the **Recent Files** submenu that you want. The value ranges from **0** to **8**, the default is **4**;
2. One can also check the **Clear the Recent Files when exit the application** to clear the **Recent Files** after exit ToupView application.

6.14 Exit

Choosing the **File>Exit** command will close video, all of the image windows and **Browse** window and remove their windows from the screen. After all of the windows are closed, ToupView will end itself.

Note: If an image has been modified before attempting to **Exit** it, ToupView will issue a warning to ask if user want to save the image or not first.


7 Edit

7.1 Undo/Redo **Ctrl+Z**

If the **Image** or **Processing** menu's submenu command is exerted on the image, the **Edit>Undo** menu will be enabled.

Most of the operations in ToupView can be undone. Alternatively, one can restore all or part of an image to its last saved version. The basic **Undo** demo process is:

1. Choose the **File>Open Image...** command to open an image;
2. Choosing the **Image>Adjust>Auto Level** command, then the **Edit>Undo** menu will be enabled;

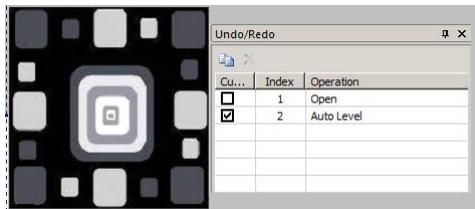

3. Choosing the **Edit>Undo** command will cancel the **Image>Auto Level** operation and return the **Operation** to its initial opened state.



If return to the **Edit** menu, one will find that the **Edit>Undo** menu now is changed to **Edit>Redo**. You can select one of these two operations to understand the **Undo** and **Redo** mechanisms:

1. Choosing the **Edit>Redo** command will return the image to the **Image>Auto Level Operation** and the **Edit>Undo** menu will be enabled again;
2. Choosing the **Image>Auto Contrast** command, then the **Edit>Redo** menu will change to **Edit>Undo**;

Note: ToupView supports only one step **Undo** and **Redo** operations.

7.2 Forward


The **Edit>Forward** menu will be enabled if it is not in the final step in the **Undo/Redo Sidebar**.

The **Edit>Backward** menu will be enabled if it is not in the first step in the **Undo/Redo Sidebar**.

The **Edit>Forward** command will forward the current displayed image to the next operation listed in the **Undo/Redo Sidebar** (If it is not in the final step).

Forward Demo

1. Open an image, choose the **Image>Adjust>Auto Level** and the **Image>Adjust>Invert** commands; check **Current** on **Index 1(Open Operation)**. The **Undo/Redo Sidebar** is shown below in Fig.1

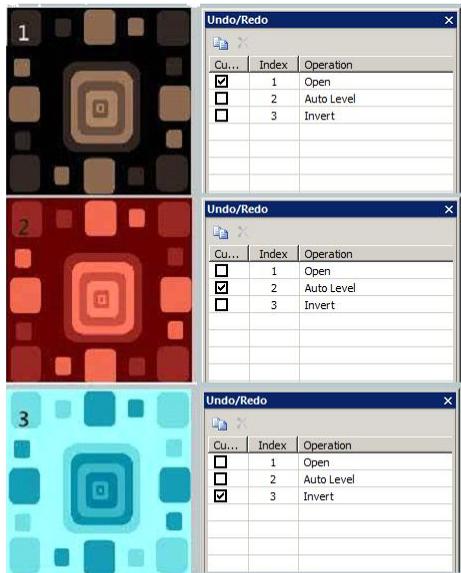

Now we continue the **Edit>Forward** menu demo. Since it is in **Index 1**, the **Edit>Backward** menu is disabled and the **Edit>Forward** menu is enabled.

2. Choosing the **Edit>Forward** command, the image and the **Index** will **Forward** to

Fig.2. Now the **Edit>Backward** menu is enabled;

3. Choose the **Edit>Forward** menu again to go to **Index 3** as shown in Fig 3. Now the **Edit>Backward** menu is still enabled, but because it is in the final operation, the **Edit>Forward** menu is disabled.

See **Edit>Backward** for further information.

7.3 Backward

The **Edit>Forward** menu will be enabled if it is not the final operation in the **Undo/Redo Sidebar**.


The **Edit>Backward** menu will be enabled if it is not the first operation in the **Undo/Redo Sidebar**.

The **Edit>Backward** command will move the current displayed image to the previous **Index** listed in the **Undo/Redo Sidebar** (If it is not in the "Open" status).

Backward Demo

1. Choose the **Open Image●●●** command to open an image;
2. Exert the **Image>Adjust>Auto Level** command on the image;
3. Choose the **Image>Adjust>Invert** command, the final image is shown in Fig 3. Since it is in the final operation state (Fig.3), the **Edit>Backward** menu will be enabled;
4. Choosing the **Edit>Backward** command, the image and the **Index** will **Backward** to **Index 2** as shown in Fig 2. Since it is in **Index 2**, the **Edit>Forward** menu will be enabled;
5. Choosing the **Edit>Backward** command again, both the image and the **Index** will **Backward** to **Index 1** as shown in Fig.1. Since it is in **Index 1**, the **Edit>Forward** menu will still be enabled, but the **Edit>Backward** menu will be disabled.

See **Edit>Forward** for further information.

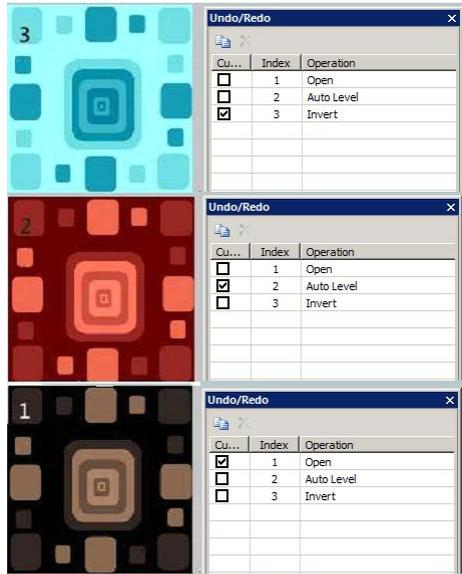

7.4 Cut Ctrl+X

The **Edit>Cut** menu will be enabled only when an **Object** or **Objects** on the **Layer** over the image is or are selected.

Check the **Measurements>Object Select**  or **Edit>Select All** menu about how to select **Layer** objects for the **Edit>Cut** menu function.

Choose the **Edit>Cut** command to copy the selected **Measurement Objects** to the clipboard and delete the selected **Objects** on the image. Any data already exist on the clipboard will be replaced.

The data copied to the clipboard can be pasted into the active window or into another opened video/image window on the layer above the image using the **Edit>Paste** command (when there is no **Layer** over the **Background**, create a new **Layer** with the **Layer>New...** command first).

Note: The **Edit>Cut** command does not support **Background Layer** (image) **Cut** operation.

7.5 Copy Ctrl+C

When the **Measurement Object(s)** is (are) highlighted, the **Edit>Copy** menu will be

enabled.

When an image ROI is selected and the **Current Layer** is the **Background Layer**, the **Edit>Copy** menu will be enabled.


Choose the **Edit>Copy** command to **Copy** the highlighted **Objects** (on **Measurement Layer**) or an image's selected area on the **Background Layer** to the clipboard.

Copy the selected area on the background layer to the clipboard.

1. Select the image area using the **Image Select**  button on the toolbar. The **Edit>Copy** menu will be enabled;
2. Choose the **Edit>Copy** command to copy the selected image area to the clipboard.

Copy object(s) on the measurement layer to the clipboard.

1. For the **Layer** operation, see the **View>Sidebar>Layer** menu in Sec.8 and the **Layer** menu in Sec.13 for details;
2. For the **Measurement** operation, see the **View>Sidebar>Measurement** menu in Sec.8 and the **Measurements** menu in Sec.14;
3. After the **Measurement** operation has been done, choose **Measurements>Object Select** command or check the **Object Select** button , the cursor will change to ;
4. Move the mouse until the cursor becomes , this means the cursor is right on the **Object**. Clicking it will highlight **Object** and select the **Object**;
5. Option 1: Continue to move the mouse until the cursor becomes  again, this means the cursor is right on another **Object** again. Clicking it with **SHIFT+left mouse button** and the second **Object** will be selected and highlighted;
6. Option 2: a) Move the cursor over the image, click down the left mouse button; b) Drag the mouse to draw a rectangle on the image. A dotted rectangle will appear around the selected area; c) Release the mouse and all of the **Measurement Objects** within the dotted rectangle will be highlighted and selected;
7. After the **Measurement Objects** are selected, the **Edit>Copy** menu will be enabled;

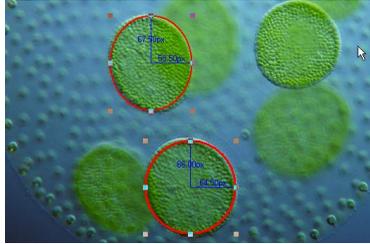

8. Choose the **Edit>Copy** to **Copy** the **Object(s)** to the clipboard. Then the **Edit>Paste** menu will be enabled. One can then **Paste** the objects onto the **Current Layer** or onto the other **Measurement Layer** in the same video/image window. If one switches to the **Background Layer**, the **Edit>Paste** menu will be disabled, but if one returns to the **Measurement Layer**, the **Edit>Paste** menu will be enabled again.

Note: a) The **Edit>Copy** command will not delete the **Measurement Objects** over the image. Any data existing on the clipboard will be replaced with the new data; b) The copied **Object(s)** can be pasted into the active window or into another opened window using the **Edit>Paste** command as long as the current window is not on the **Background Layer** (the **Edit>Paste** menu will be disabled if the **Background Layer** is active). See the **View>Sidebar>Layer** command in Sec.8 and the **Layer** menu in Sec.13 for details.

7.6 Paste **Ctrl+V**

If there is **Measurement Object** data in the clipboard and the **Current Layer** is not the **Background Layer**, the **Edit>Paste** menu will be enabled.

Choose the **Edit>Paste** command to **Paste Objects** from the clipboard onto the active image's **Measurement Layer**. One can also choose **Edit>Paste** command to transfer a layer's **Measurement Objects** from one image window's **Measurement Layer** to another image window's **Measurement Layer**.

Note: The **Edit>Paste** command does not support the image area **Paste** operation.


7.7 Image Select

The **Edit>Image Select** command can be used to mark **ROI** and **Copy** the selected **ROI** to the clipboard. This command is only used to select the **ROI** on the **Background Layer**.

Choosing the **Edit>Image Select** command will check this menu (or click  on the toolbar will keep it down) and the cursor will turn into “+”. After the area is selected and the **Current Layer** is the **Background Layer**, the **Edit>Copy** menu will be enabled and then the selected area can be copied to the clipboard for further application.

To select an area,

1. Choose **Edit>Image Select** command (Or click the **Image Select** button  on the toolbar)
2. Drag the mouse cursor across the image with the left button held down until the area is selected.
3. Release the button and the area will be marked. Handles will appear on the area that will allow altering the selection after it is marked.

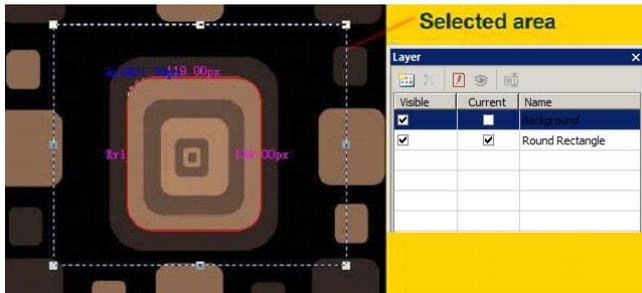

Note: Only when a ROI is selected and the **Current** item is checked on the **Background** in the **Layer Sidebar**, the **Edit>Copy** command can copy the selected image ROI to the clipboard.

7.8 Select All Ctrl+A

7.8.1 Select all on the background layer

When the video/image window is active and the **Background Layer** is checked, choosing the **Select All** command to select all pixels on the **Background layer** within the canvas (shortcut: **Ctrl+A**).

7.8.2 Select all objects over the background layer

When the video/image window is active and the **Background Layer** is not checked, choosing the **Edit>Select All** command will select all of the **Objects** on the **Current**

Layer (shortcut: **Ctrl+A**).

7.8.3 Select all files in the browse window

When the **Browse** window is active, choosing the **Edit>Select All** command will select all of the files in the **Browse** window under the specified directory in the **Folders Sidebar**. See **Edit>Invert Selection** for select all files method.

7.9 Select None **Ctrl+D**

Deselect any selected area on the image or the **Measurement Objects** on a Layer.

1. When the **Current Layer** is the **Background Layer** and an image area is selected, the **Select None** menu will be enabled. Choosing the **Edit>Select None** command will delete the dotted rectangle representing the selected area;
2. When the **Current Layer** is not the **Background Layer** and the **Measurement Objects** are selected, the **Edit>Select None** command will be enabled. Choosing the **Edit>Select None** command will deselect all of the selected **Measurement Objects**.

Note: See **Edit>Image Select**, **Edit>Select All** and **Measurements>Object Select**  to understand how to perform select operations.

7.10 Inverse Selection

This command is for **Browse** window only.

When organizing files, one can invert a selection in order to select all of the other files that were not previously selected with the **Edit>Invert Selection** command.

When no files were selected, choosing the **Edit>Invert Selection** command will select all of files under the **Folders Sidebar** directory which is just as the **Edit>Select All** command.

7.11 Delete File **Delete**

The **Edit>Delete File** menu will be enabled if the file(s) is (are) highlighted in the **Browse** window. This command is for **Browse** window only.

You can **Delete** or remove one or more files from the **Browse** window. The steps are as follows:

1. Select one or more files by a) Clicking the displayed file icons, a single file will be

highlighted; b) Clicking the file one by one with **CTRL + left mouse button**, all of the clicked files will be highlighted; c) Clicking the displayed file icons, the first clicked file will be highlighted, clicking the end file with **Shift + left mouse button**, all of the files among the first and last will be highlighted. d) Dragging the mouse to draw a dotted line rectangle across the files you wish to delete, all of the files in the rectangle will be highlighted;

2. a) Press the **Delete** key to delete the selected files; b) Click you right mouse button to bring up a context menu, choose the **Delete** command to delete the highlighted files. A **Confirm File Delete** dialog will bring up.

In the **Confirm File Delete** dialog, click **Yes** to move the files to the desktop recycle bin, or **No** to cancel the **Delete** operation.

8 View

8.1 Browse Ctrl+B

8.1.1 Open the browse window


1. Choose the **View>Browse** menu from the **View** menu or click the **Browse** toolbar button  to browse images under the specified directory in the **Folders Sidebar**;
2. Click the **Folders Sidebar** to activate it and double-clicking the listed directory in the **Folders Sidebar** will create the **Browse** window.

After creating the **Browse** window, ToupView will display a **Browse** window that looks like windows explorer. The child window on the left part of the **Browse** window called **Folders Sidebar** is used to locate the directory on the hard disk. Images in the current directory are displayed in **Large Icons** or **Small Icons** mode on the right side of the **Browse** window.

Their order can be set in **Forward** or **Reverse** format according to **Sort by Name, Type, Size, Width** or **Height** et al.

8.1.2 Browse window right mouse button context menu

Clicking the right mouse button on the listed **Icon** in the **Browse** window will bring up a right mouse button context menu as shown below:

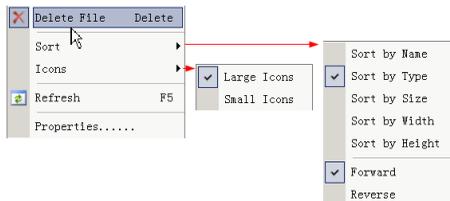

These context menu functions are described in

View>Delete File

View>Sort>Sort by Names

View>Sort>Sort by Type

View>Sort>Sort by Size

View>Sort>Sort by Width

View>Sort>Sort by Height

View>Sort>Forward

View>Sort>Reverse

View>Icon>Large Icons

View>Icon>Small Icons

View>Refresh for details

Note: The **Folders Sidebar** and the **Browse** can be used to perform tasks such as creating new folders, renaming, moving, and deleting files. Individual file information and import data from digital cameras can also be displayed. Double-clicking the left mouse button on the icon will open the image as an active image in full size. See **Image window UI** in Sec.3 for more details.

8.2 Sort>Sort by Names

This command is for the **Browse** window only.

Sort the image files in order of names in the **Browse** window.

8.3 Sort>Sort by Type

This command is for the **Browse** window only.

Sort the image files in order of type in the **Browse** window.

8.4 Sort>Sort by Size

This command is for the **Browse** window only.

Sort the image files in order of size in the **Browse** window.

8.5 Sort>Sort by Width

This command is for the **Browse** window only.

Sort the image files in order of image width in the **Browse** window.

8.6 Sort>Sort by Height

This command is for the **Browse** window only.

Sort the image files in order of image height in the **Browse** window.

8.7 Sort>Ascending

This command is for the **Browse** window only.

Sort the image files in order of the **Ascending** mode (i.e. 1, 2, 3 and 4) in the **Browse** window.

8.8 Sort>Descending

This command is for the **Browse** window only.

Sort the image files in order of the **Descending** mode (i.e. 4, 3, 2 and 1) in the **Browse** window.

The **Sort** settings are saved until they are changed. For example, if you sort images in the **Browse** window according to the **Type**, the images will remain sorted according to **Type** until the **Sort** settings are changed.

8.9 Icon>Large Icons

This command is for the **Browse** window only.

The **Thumbnail** mode displays small preview images; you can select different view modes in the **Browse** window.

Choosing the **Icons>Large Icons** will display the image files in **Large Icon** mode in the **Browse** window.

8.10 Icon>Small Icons

This command is for the **Browse** window only.

The **Thumbnail** mode displays small preview images; you can select different view modes in the **Browse** window.

Choosing the **Icons>Small Icons** will display the image files in **Small Icon** mode in the **Browse** window.

8.11 Refresh F5


This command is for the **Browse** window only.

If the files under the **Folders Sidebar**'s directory are altered outside of ToupView, after switch back to ToupView, one can **Refresh** the image files in the current directory to update the **Thumbnails** with **View>Refresh** command.

8.12 Measurement Sheet

Help Manual

When choosing the **View>Measurement Sheet** command, the **Measurement Sheet** will be active. The **Measurement Sheet** shows the **Object's** possible features, such as **Name, Center, Radius, Area, Perimeter, Angle, Start Point, and End Point** overlaid on the extra layer.

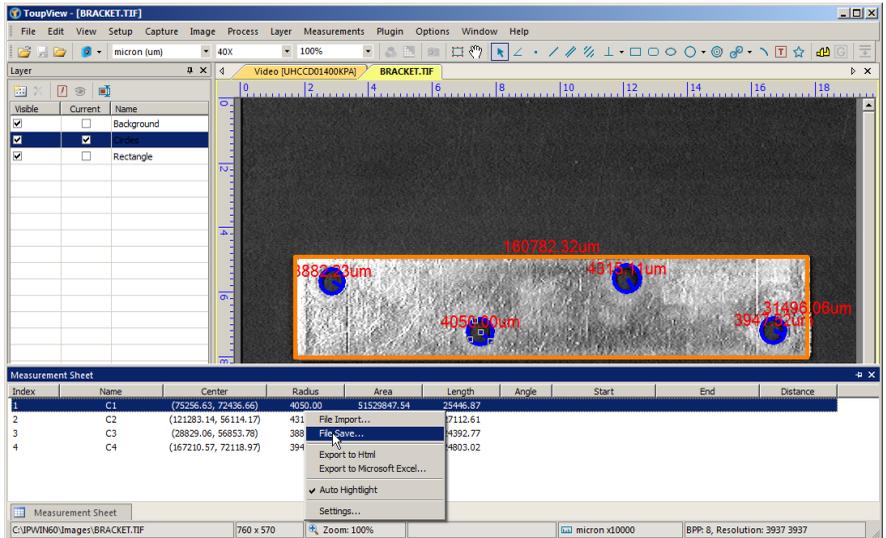

Clicking the right mouse button on the **Measurement Sheet** and the right mouse context menu will bring up on the **Measurement Sheet** window as shown above. They are explained below:

8.12.1 File Import

Choose this command to load a **Measurement files (*.measurement)** and overlay the **Measurement Objects** over the current video/image background.

8.12.2 File Save

Choose this command to save the **Measurement Objects** over the current video/image on the **Current Layer** to a **Measurement files (*.measurement)**.

8.12.3 Export>To Html File

Export all the **Layer Measurement Objects** to the ***.html** file in tabbed format.

Help Manual

Layer Name	Index	Name	Center	Radius	Area	Length	Angle	Start	End	Distance
Layer 1	1	L1	(837.00, 627.00)			712.12	58.32	(650.00, 930.00)	(1024.00, 324.00)	
Layer 2	1	R1	(1268.00, 400.00)		259072.00	2176.00		(1636.00, 576.00)	(900.00, 224.00)	
	2	R2	(1755.00, 1110.00)		186416.00	1740.00		(1946.00, 1354.00)	(1564.00, 866.00)	

Length Unit:P, Angle Unit:Pi

Note: This menu will be enabled only when there are **Measurement Objects** over the **Background Layer** (video/image).

8.12.4 Export to Excel

If this command is chosen, the **Current Layer Measurement Objects** over the video/image will be exported to **Microsoft Excel**. The **Objects'** parameters in the **Measurement Sheet** will also be arranged as a table.


A	B	C	D	E	F	G	H	I
Name	Center	Radius	Area	Length	Angle	Start	End	Distance
R1	(1268.00, 400.00)		259072	2176		(1636.00, 576.00)	(900.00, 224.00)	
R2	(1755.00, 1110.00)		186416	1740		(1946.00, 1354.00)	(1564.00, 866.00)	

Note: This menu will be enabled only when there are **Measurement Objects** on the **Current Layer** (video/image).

8.12.5 Auto Highlight

When this menu is checked, clicking the row in the **Measurement Sheet** will **Highlight** the corresponding **Measurement Object** over the **Background Layer**. Clicking the **Object** over the video/image will **Highlight** the corresponding row in the **Measurement Sheet**.

8.12.6 Settings●●●

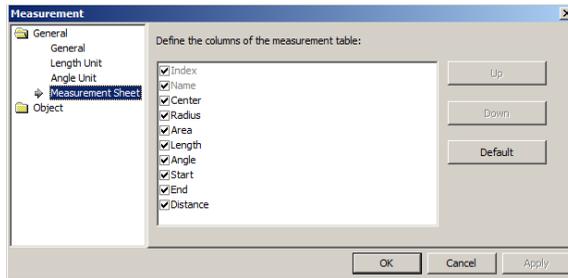

1. To modify the **Measurement Sheet's** item order, click the item to highlight it, the **Up** or **Down** button will be enabled (If the item is in the third position, the **Up** button will

be disabled, if the item is in the last row, the **Down** button will be disabled). Click the **Up** or **Down** button to modify the item order;


Note: The item **Index** and **Name** are always in grayed states, this means these two items cannot be modified (Both their hide/show states and their positions).

2. Checking/Unchecking the item will show/hide the item in the **Measurement Sheet**;
3. Clicking **Default** will return to the ToupView's default settings;
4. Item 1(**Index**) and item 2(**Name**) cannot be changed.

8.13 Sidebar

There are 5 **Sidebar** in the ToupView frame window in tabbed format. They are **Camera Sidebar**, **Folders Sidebar**, **Undo/Redo Sidebar**, **Layer Sidebar** and **Measurement Sidebar**.

8.13.1 Sidebar overview

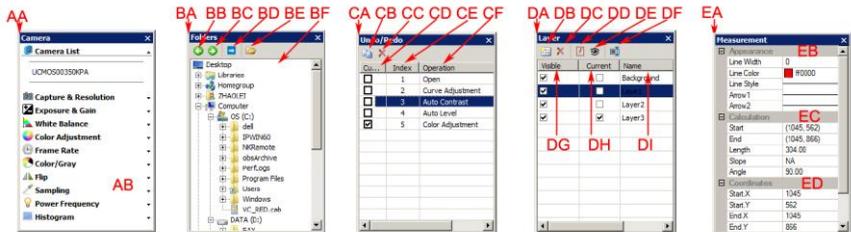

AA: **Camera Sidebar**; AB: **Camera Sidebar** groups for the control of the started up camera.

BA: **Folders Sidebar**; BB: **Back** to the previous folder;

BC: **Forward** to the next folder;

BD: **Browsing** the pictures under the directory of ToupView's directory;

BE: Auto hide button to show/hide the current selected directory, this will open the **Browse** window if it is not opened (Double-clicking on the selected directory will perform the same functions.);

BF: **Folders** to locate the **Browse** window's file directory.

CA: **Undo/Redo Sidebar**;

CB: **Operation Copy**: Copy the highlighted operation in the **Undo/Redo Sidebar** to a new image window; User can also drag the selected operation in the **Undo/Redo Sidebar** to the window area to create a new image window;

CC: **Remove** the highlighted operations from the **Undo/Redo** list (This button will be enabled only when the operation(s) is (are) selected, the opened image(the first operation) cannot be removed);

CD: Indicating the **Current** operation displayed in the image window;

CE: **Operation Index**; CF: **Operation** name.

DA: **Layer Sidebar**; DB: Make a **New** layer; DC: **Remove** a layer;

DD: Set as the **Current** layer; DE: **Show/Hide** a layer; DF: **Rename** a layer;

DG: Visibility control of the layer items;

DH: The **Current** active layer for operations;

DI: The layer **Name**. The image layer is always named as "**Background**".

EA: **Measurement Sidebar**;

EB: The **Appearance** of the highlighted **Object** on the **Current** layer; you can edit the **Appearance** by clicking its item and editing it;

EC: The **Calculation** of the highlighted item on the **Current** layer;

ED: The **Coordinate** of the highlighted item on the **Current** layer; you can edit the **Coordinate** by clicking the item.

8.13.2 Sidebar>Camera

The **Camera Sidebar** is mainly used for the control of Toupcam camera, it includes 11 groups. Every group can be expanded by clicking the group name or the arrow button at the right of the group name.

Checking the **Sidebar>Camera** menu will activate/show it on the frame.

Please check **Camera Sidebar** in Sec.5 for details.

8.13.3 Sidebar>Folders

The **Folders Sidebar** is mainly used for the image **Browse** control.

Checking the **View>Sidebar>Folders** menu will activate/show the **Folders Sidebar**.


Clicking its tree can navigate the file directories.

Help Manual

Double-clicking the directory in the **Folders** will create the **Browse** window. If there are image files under the tree that ToupView supports to browse, the image files will be displayed in **Large** or **Small** icons mode.

Their order can be set in **Forward** or **Reverse** format according to **Sort** by **Name**, **Type**, **Size**, **Width** or **Height** et al.

Clicking with the right mouse button on the directory will bring up the right mouse context menu as shown below:

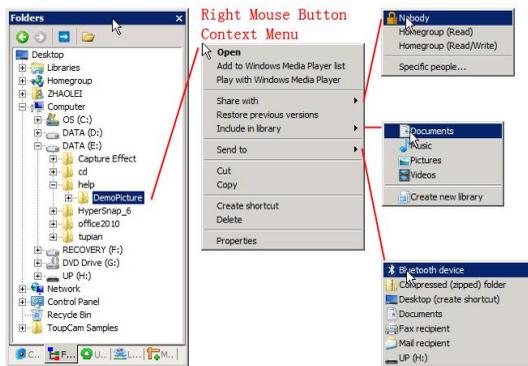

This is the basic window explorer menu and will not explain it at here in this manual.

8.13.4 Sidebar>Undo/Redo

The **Undo/Redo Sidebar** is used to **Undo/Redo** operations for the **Image** and **Process** menus' command.

Checking the **View>Sidebar>Undo/Redo** menu will activate/show the **Undo/Redo Sidebar**.

8.13.5 Sidebar>Layer

Layer Sidebar is used for the management of **Layer** operations. This operations including making a **New Layer**, **Removing** a **Layer** or **Renaming** a **Layer** and **Layer** visibility controlling et al.

Checking **View>Sidebar> Layer** will activate/show the **Layer Sidebar**.

8.13.6 Sidebar>Measurement

The **Measurement Sidebar** is used to check or edit the selected **Layer Objects** on the **Layer**.

Checking **View>Sidebar>Measurement** will activate/show the **Measurement Sidebar**.

8.14 Ruler and Grid


Ruler and Grid menu has 3 submenus, they are:

8.14.1 Grids>No Grids

Choosing this command will remove both **Manual Grids** and **Auto Grids** overlaid on the video/image window.

8.14.2 Grids>Manual Grids

Choosing this command will display two small **Right Arrow** and **Down Arrow** overlaid on the top of the **Vertical Ruler** and on the left of the **Horizontal Ruler** as shown below:

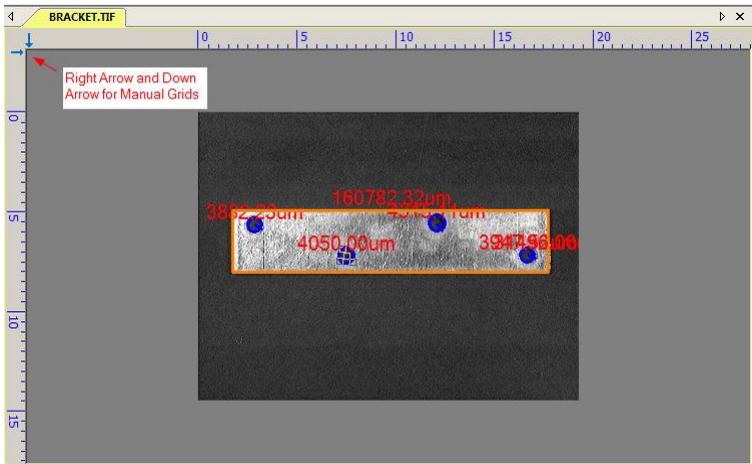


Move the mouse over the **Down Arrow** will show a horizontal drag icon. Drag the **Down Arrow** along the **Horizontal Ruler** to where ever you want. When it is dragging over the video/image, there will be a **Vertical** line appeared to let you judge where to release this line on the video/image. You can drag any lines to overlay them on the video/image.

Help Manual

The overlaid lines can be adjusted by dragging the specific **Down Arrow** again on **Horizontal Ruler**.

Move the mouse over the **Right Arrow** will show a vertical drag icon. Drag the **Right Arrow** along the **Horizontal Ruler** to where ever you want. When it is dragging over the video/image, there will be a **Horizontal** line appeared to let you judge where to release this line on the video/image. You can drag any lines to overlay them on the video/image.

The overlaid lines can be adjusted by dragging the specific **Right Arrow** again on **Vertical Ruler**.

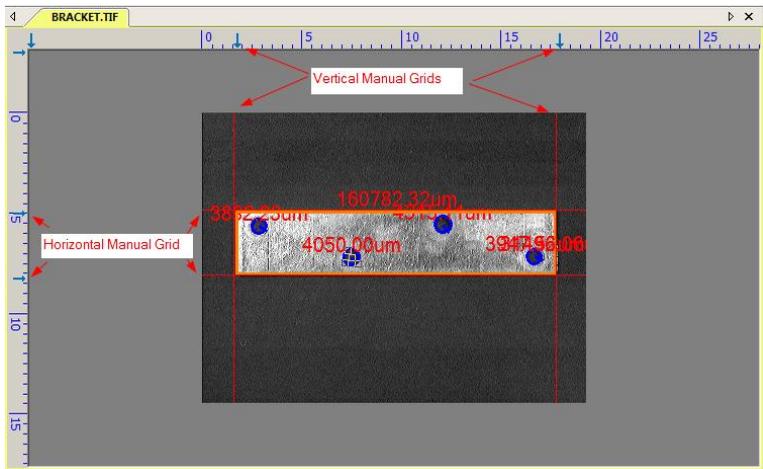

Choose the **Grids>No Grid** command to remove the **Manual Grids** overlaid on the video/image window. The menu will be checked.

Choose the **Grids>Manual Grids** command again (If it is unchecked) to display all the previous **Manual Grids** overlaid on the video/image window. The menu will be checked.

8.14.3 Grids>Auto Grids

Choose the **Grids>Auto Grids** command to overlay the grids on the video/image automatically.

The **Auto Grids** can be set in the **Ruler and Grids>Setting●●●** menu.

8.14.4 Grids>Remove All Grids


Remove (Delete) all of the **Manual Grids** or **Auto Grids** overlaid on the image/video

window.

8.14.5 Settings...

Choosing the **View>Ruler and Grid>Setting...** command, or choosing the **Preference>...** command and clicking the **Rulers and Grids** page realize the same function.

Choosing the **Preference>...** command and clicking the **Rulers and Grids** page can set **Ruler Color**, **Cursor Color** of the **Horizontal Ruler** and **Vertical Ruler** for the video or image windows. The **Grid Style**, **Line Style**, and **Line Color** for the **Grid** overlaid on the video and image window can also be set.

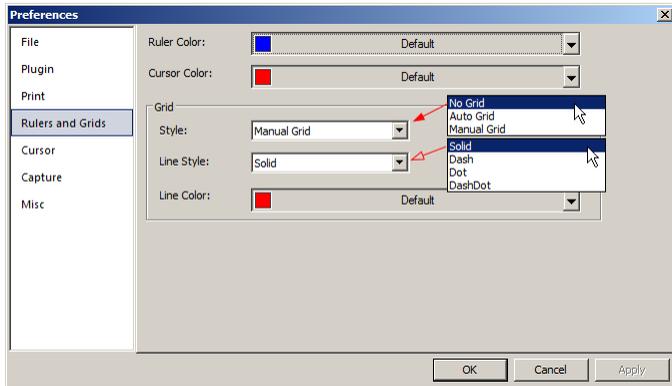

Ruler Color: The **Horizontal Ruler** color and **Vertical Ruler** color; Default is blue (0,0,255);

Cursor Color: The **Cursor** color; When you are moving your mouse over the video/image window, there will be **Horizontal Cursor** and **Vertical Cursor** move on the **Horizontal Ruler** and **Vertical Ruler** moving at the same time indicating the current **Cursor** position in both directions. Default is Red (255,0,0);

Grid Style: The **Grid Style** can be **No Grid**, **Auto Grid** or **Manual Grid**. Default is **No Grid**;

Grid Line Style: The **Line Style** for the grid can be **Solid**, **Dash**, **Dot** or **DashDot** et al. Default is **Solid**;

Grid Line Color: The color of the grid line. Default is Red (255,0,0).

8.15 Best Fit

Choose the **View>Best Fit** command to automatically resize the video/image to fit in the window.

Note: Choosing this command will enable **View>Actual Size** menu.

8.16 Actual Size

Note: This option will be disabled if the image is currently viewed at 100%. At any other zoom ratio, **View>Actual Size** will be enabled.

Choose the **View>Actual Size** command to set the active image to its **Actual Size** (e.g. 100%).

8.17 Track

Note: If the video/image size is smaller than the window size. The track operation will be disabled.

If the image's actual size is larger than the video/image window, check this command to position the video/image that does not fit entirely within in the video/image window. Its function is similar to the scroll bars. It is an alternative to using the arrows on the scroll bars for positioning the video/image within the window.

Checking the **View>Track** menu will change the cursor to  and the button on the toolbar will be checked.





Then keep down the mouse button to drag the region of interest on the video/image to any location in the video/image window.

8.18 Properties

If an image file listed in the **Browse** window is highlighted. Choosing the **View>Properties** command or clicking the selected file in the **Browse** window with the right mouse button will bring up a **Properties** dialog as shown below:

The file **Properties** dialog including 4 pages. They are **General**, **Security**, **Details** and **Previous Versions** pages. These **Property pages** may depend on the operating system and we will not discuss it in this help manual.

Help Manual

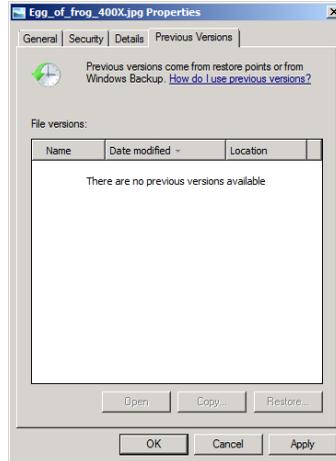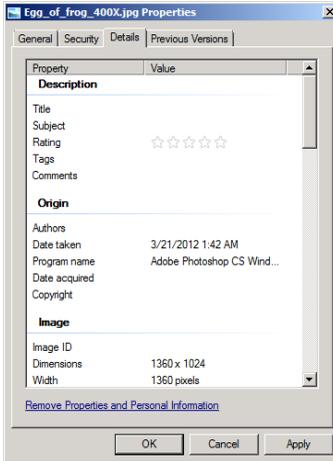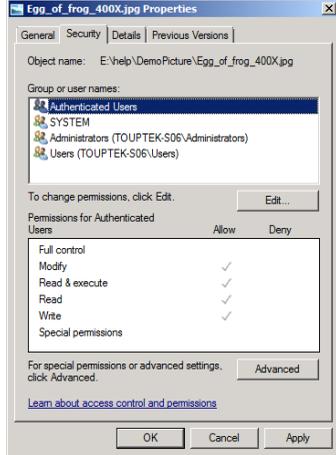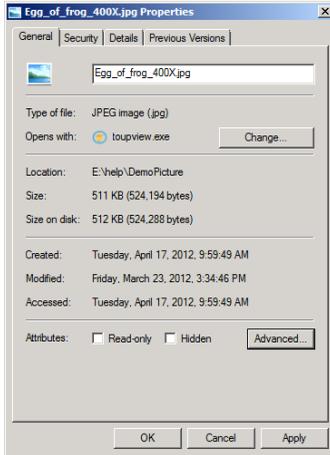

9 Setup

9.1 Start/Pause Pause

If the window is running, one can choose the **Setup>Start/Pause** command to pause the video and choose the **Setup>Start/Pause** command again to start the video.


9.2 Full Screen ESC

Choosing the **Setup>Full Screen** command will display the video window in full screen style. Choosing the **Setup>Full Screen** command again will return to the default video window.

Note: The window should be in focus to make the **Setup>Full Screen** or **ESC** command enabled.

9.3 View Properties●●●

The **Setup>View Property●●●** will help you to understand the camera statistical properties. Choose the **Setup>View Property●●●** command to invoke the **View Property** dialog as below:

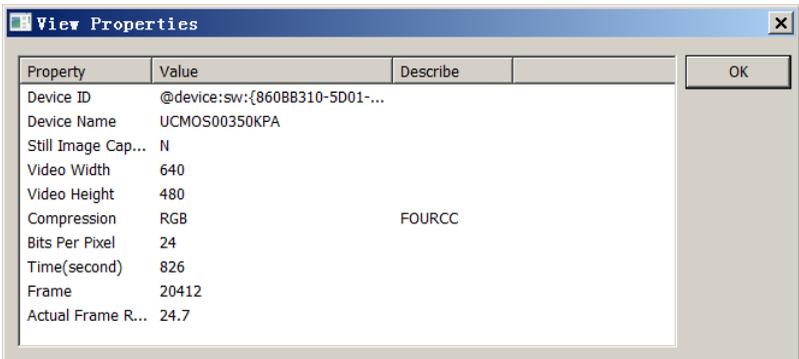

The items in the dialog are described in the following table:

Item	Description
Device ID	Unique ID to identify the camera device.
Device Name	Human readable string to identify the name of the camera device.

Help Manual


Still Image Capture	Whether or not the camera supports Still Image Capture . Still Image Capture is used for high resolution camera to capture an image with a different resolution from the video. This feature is mainly used to capture high resolution image under low resolution video to compromise the frame speed and the image resolution.
Display Width	The video window width.
Display Height	The video window height.
Video Width	The actual video window video width.
Video Height	The actual video window video height.
Compression	The compression format of the video stream.
Bits Per Pixel	Indicate how many bits are used to store on pixel.
Time (second)	Seconds elapsed since the video has been started.
Frame	Frames acquired since the video has been started.
Actual Frame Rate	Actual frame Rate of the video stream.

Note: The **Actual Frame Rate** is listed for reference. It varies depending on the computer's configuration. Different hardware configurations may have different **Actual Frame Rates**.

9.4 Video Overlay●●●

9.4.1Video Overlay: Overlay

Choose the **Setup>Video Overlay●●●** command and click the **Overlay** page to overlay **Scale**, **Magnification**, **Date** and **Clarity Factor** on the video window.

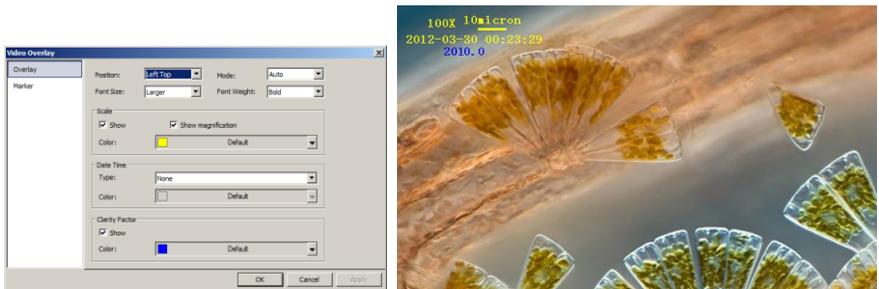

This command will invoke **Overlay** page as above. The **Position**, **Font Size**, **Font Weight** of the **Scale**, **Magnification** and **Date**, and **Clarity Factor** can be defined


together. Their **Colors** can be defined separately.

Clicking the **OK** button and the **Scale**, **Magnification** and **Date**, and **Clarity Factor** will be overlaid on the **Video** window. The **Clarity Factor** can tell if the sample is in good focused state. The larger the **Clarity Factor**, the better the sample focused.


Note: To enable the **Scale** bar, the **Magnification** must be defined and chosen first. The **Unit** can be any unit except **Pixel**. There are two methods to set the **Unit**, they are:
a) Choosing the **Unit** in the **Unit** dropdown combobox (**micron (um)** **10X** **100%**) on the tool bar which is just on the left of the **Magnification** dropdown list box; b) Choosing the **Option>Measurement...** command, a dialog called **Measurement** will bring up, click the **Length Unit** page and check the **Unit** in the **Current** to set the **Unit**.

9.4.2Video Overlay: Marker...

Choose the **Video Overlay: Marker...** command to overlay **Video Marker** on the video window. The **Video Marker** type can be **Cross**, **Rectangle**, **Circle**, **Cross+Rectangle**, or **Cross+Circle**. The **Video Marker** is shown as below:

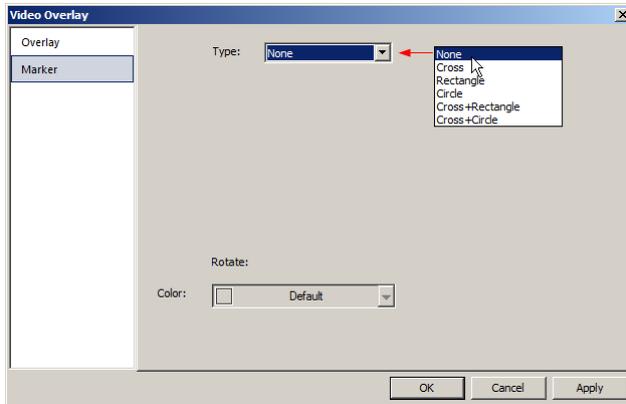


Choosing "**Cross+Retangle**" in the **Type** combobox and **Video Overlay: Marker** dialog will become:

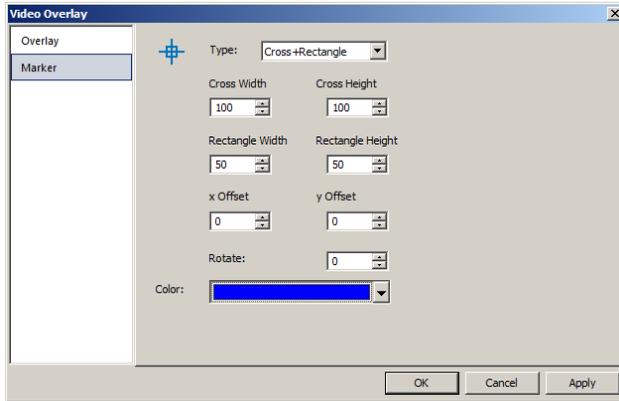

Edit the **Cross Width** and **Cross Height**, **Rectangle Width** and **Rectangle Height**, **x Offset** and **y Offset**, in their specific fields. Click **Color** to define the **Video Marker** color.

Click **OK** to end the **Video Marker** dialog and a **Cross+Rectangle Marker** will be overlaid on the video window as shown below:

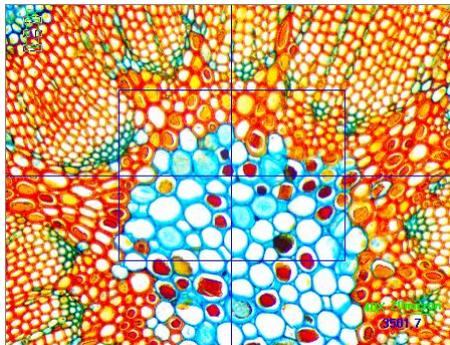

Click **Cancel** to cancel the **Video>Overlay: Marker** setup and return to the application area, or **Apply** to overlay the **Marker** on the video and keep the **Video Overlay** dialog there for further modification.

9.5 Video Watermark●●●

Fig.1 shows a **micro ruler**. The dark lines can be extracted as **Video Watermark** and overlaid on the video window. The steps are:

1. Choose the **Capture>Capture Image** command or click  to capture the **micro ruler** image as shown in Fig.1;
2. Choose the **Process>Binary** command to binarize the image as shown in Fig.2;
3. Choose the **Image>Adjust>Invert** command to invert the image and choose the **Image>Color Quantize** command to convert the image to 24 bits as in Fig.3. Choose the **File>Save As** command to save the image in **24 bit BMP** format;

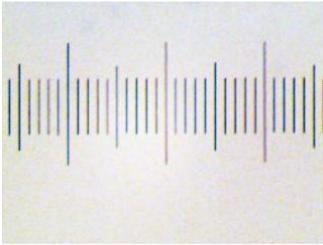


Fig.1 Captured **Micro Ruler**

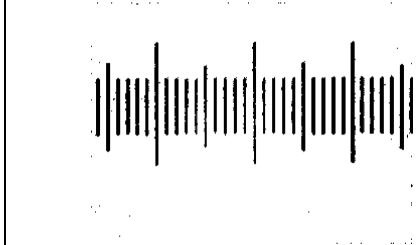


Fig.2 **Micro ruler** after being binarized

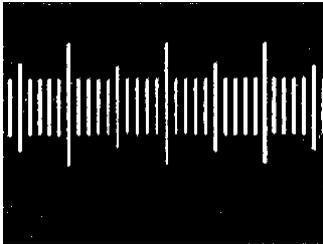


Fig.3 Inverted 24 bits image

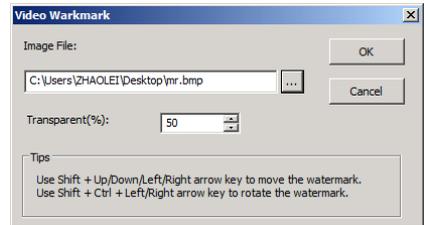


Fig.4 **Video Watermark** setup dialog

4. Choosing the **Setup>Video Watermark** command and a dialog called **Video Watermark** is brought up as shown in Fig.4. Click the  button to locate the image saved in step 3. Use the defaults **Transparent (%)**(50). If everything is ok, click **OK** button. The final **Video Watermark** is overlaid on the video window as shown in Fig.5.

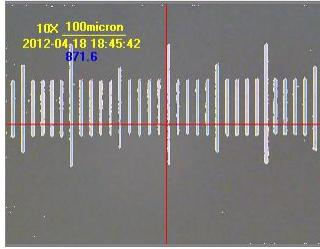



Fig.5 Video Window with Video Watermark overlaid

9.6 Move Watermark

9.6.1 Move to...

If there is **Watermark** overlaid on the video window, this menu will be enabled.

Choosing the **Setup>Watermark...** command will bring up a **Move** dialog. Where one can enter the **X:** and **Y:** offset value in their fields for the desired pixel move distances. The default **X** and **Y** offsets are 0;

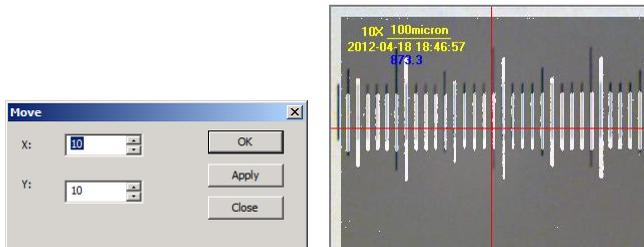

9.6.2 Move to zero


If the **Watermark** was moved, this menu will be enabled. Choosing this menu will move the **Video Watermark** to its original coordinates (0, 0).

9.7 Rotate Watermark

9.7.1 Rotate to...

If there is **Watermark** overlaid on the video, this menu will be enabled.

Choosing the **Setup>Rotate Watermark>Rotate to...** command will bring up a **Rotate** dialog, where one can define an **angle** to **Rotate** the **Video Watermark** a specified angle around the video center (0, 0) (default value).

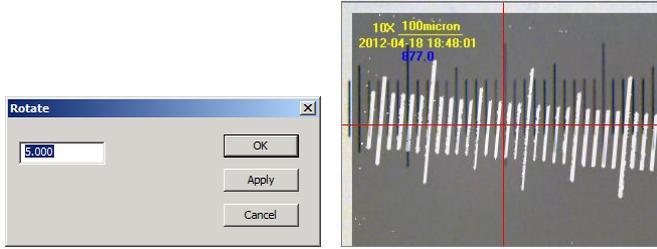


9.7.2 Rotate to zero

If the **Video Watermark** was rotated, the **Rotate to zero** menu will be enabled. Choosing this menu will rotate the **Video Watermark** to zero degree.

9.8 Gray Calibration●●●

This function can make the image brightness a desired value among various scenarios in a specified area, achieving the continuity requirement of the observation. The **Gray Calibration** steps are summarized as follows:

1. Click the **Exposure & Gain** group name on the **Camera Sidebar** to expand the **Exposure & Gain** group, uncheck the **Auto Exposure** box (If it is checked).

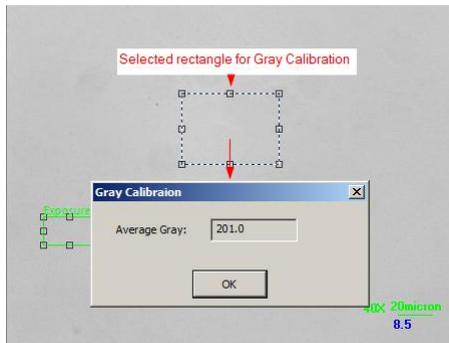

2. Choose the **Edit>Image Select** command or click the **Image Select** button  on the toolbar to select a reference region, and choose the **Setup>Gray Calibration●●●** command. A dialog called **Gray Calibration** will be brought up to display the current **ROI Average Gray**. Now the brightness of the microscope can be adjusted until the **Average Gray** reaches to the desired value. Click "OK" to finish the calibration and return to the application area. The desired value here is around is 200.

9.9 Manual Fusion●●●



Make sure that the **ToupView** package and **UCMOS** or **UHCCD** camera are correctly installed. Turn on the microscope's light.

1. Run ToupView and start the camera;
2. Choosing the **Setup>Manual Fusion●●●** command or clicking  on the toolbar will bring up a dialog called "Manual Fusion";
3. Use the microscope **coarse** or **fine** focus knobs to move the sample stage up and down, in order to find the positions where the clearest regions of the whole sample can be seen on the video window;
4. Click the **Capture** button to capture an image into the image list which will be used for the fusion operation.



5. Unless there is 1 captured image, the **Clear** buttons will not be enabled. Unless there are more than two captured images, the **Fusion** buttons will not be enabled. If the captured images are not satisfactory, click the **Clear** button to clear the captured image(s), and capture new image(s).
6. If more than two images are captured, click the **Fusion** button to do the image fusion and the **Manual Fusion** dialog will be closed automatically. If the **Fusion** button is clicked, please wait for some time to get the final fusion result. The fused image will be displayed in a new active image window and its title bar will be associated with a digit assigned by ToupView. Clicking **Close** will end the **Manual Fusion** and return to the application area.

Note: a). Use the **coarse** and **fine** focus knobs to move the sample stage up and down;
b). Clicking the **Capture** button will capture the current image into the image fusion list;
c). Repeating steps 1-4 until there are enough images;
d). Clicking the **Fusion**


button will start image fusion. Waiting for some time and a fantastic fusion result will be displayed in a new image window;

9.10 Video Source Property●●●

This menu will be displayed and enabled only when the camera that supports DirectShow interface is started.

Choose the **Setup>Video Source Property●●●** command to invoke the **Video Source Property** dialog:

The **Video Source Property** includes several categories such as **Color**, **Exposure**, **Extended**, **Misc** and **ROI** page et al. These pages are used for the control of your camera.

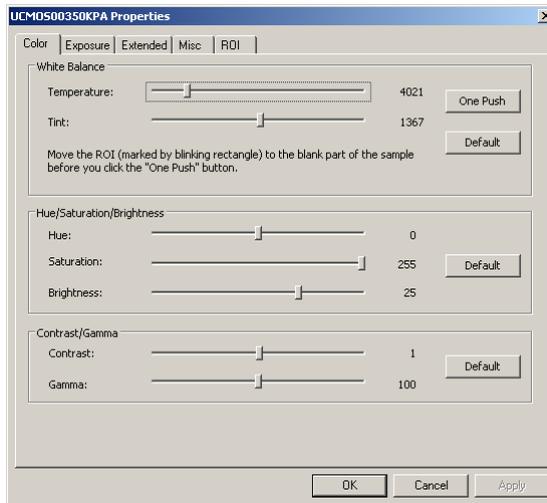


Note: Different camera may have different control contents. If you have any problems with these controls, try to contact your camera supplier to get the further technique help. ToupView has no responsibility to help you about these controls.

9.11 Video Stream Format●●●

This menu will be displayed and enabled only when the camera that supports DirectShow interface is started.

The ToupView **Video Stream Format** configuration dialog is shown below:

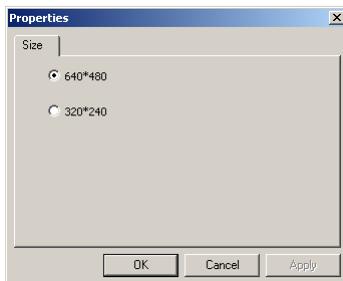

One can change the video format mode, frame rate, color space, compression options and so on. Here, we only have the **Video Size** to check. Selecting the desired one and clicking **OK**. This will give you a selected **Video Size** for the new video window.

Click **Cancel** to cancel the selection and return to video window without any changes or **Apply** to apply the current selection to the video with the **Properties** remain opened.

Note: Different devices may have different user interfaces and items. Contact the camera supplier for details.


9.12 Still Image Options●●●

Some cameras can capture a still image with its sizes different from the video stream. The camera may have a "pin" that acts as a hardware trigger, or it may support software triggering. A camera that supports still images capture will expose a still image pin to the designer.


To capture a still image, one should first set the size of the image to capture. The size is determined by the camera hardware. Here we show an example with **UCMOS08000KPA** camera.

Choosing the **Setup>Still Image Options●●●** command will show the following **Properties** dialog:

Help Manual

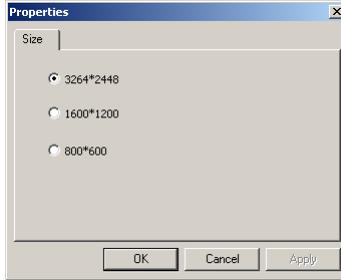

It has 3264*2448, 1600*1200, and 800*600 resolutions. Select the desired one and press **OK** to accept the selection, or **Cancel** to ignore the current selection. Click **Apply** to apply the selection.

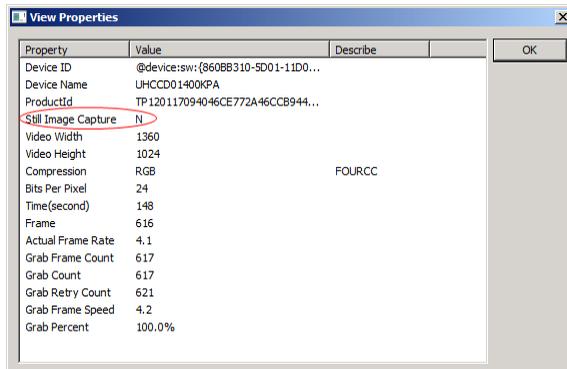

One can find if the device supports **Still Image Capture** or not by choosing the **Setup>View Property...** command.

If the **Still Image Capture's Value** is **Y** in the **View Property** dialog, this means that the camera supports the **Still Image Capture** operation.

10 Capture

10.1 Capture Image F8

During the video preview, you can always choose the **Capture>Image Capture** command to capture the video image.

After the image is captured, the captured image will be the current active window. The **Capture>Capture Image** menu will be disabled. If you wish to capture image again, click the video window title to activate the video window and the **Capture>Capture Image** menu will be enabled again.

Note: a) The "Snap" button  on the **Camera Sidebar** can continuously shoot the image even if the video window is not activated. Click this button on the **Camera Sidebar** to capture image quickly; b) Only when the video window is active, the **Capture>Capture Image** menu will be enabled; c) If the **Live** and **Snap** resolutions are different, ToupView need to switch the resolution from **Live** to **Snap** in the background to capture an image with **Snap** resolution. After the **Snap** is finished, ToupView will switch back to the **Live** resolution to continue the video stream process. Thus it will take more time to capture a still image.

10.2 Time-lapse (Auto Capture)●●●

This function can capture a sequence of pictures with the same time interval; you can precisely set the time interval (2 to 3600 seconds) and the total number of images. Choosing the "**Capture>Time-lapse (Auto Capture)●●●**" command will bring up **Time-lapse (Auto Capture)** dialog shown below:


Directory: The file **Directory** can be selected by clicking the **Browse** button.

File: The **File** name including **Name of Format**, **File Prefix**, **File type**. It can be a combination of prefix, time and type and is shown in the **Sample**.

Time Slot: **Time slot(Second,2-3600S)** is a time segment to capture an image.

Total Images: Checking **Total Images** will enable its edit box. You can enter the **Total Images** (1-9999) to be captured. ToupView will stop the **Time-lapse** capture process automatically when the **Total Images** are reached.

If **Total Images** is unchecked, ToupView will capture the images continuously until you choose the **Capture>Stop Time-lapse (Auto capture)** command again to stop the **Time-lapse** capture.

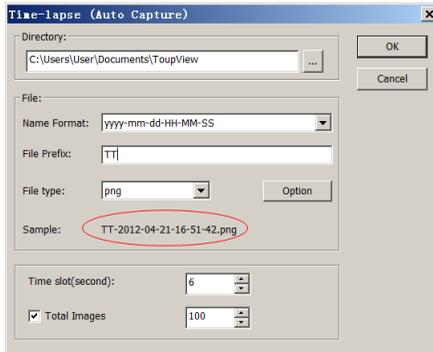

Click **OK** to begin the **Time-lapse** capture, or **Cancel** to cancel the **Start Time-Lapse(Auto Capture)...** command.


After the **Time-lapse** capture is started, the **Capture>Start Time-Lapse (Auto Capture)...** menu will be changed to the **Capture>Stop Time-lapse (Auto capture)** menu. Choosing this command will stop the **Time-lapse** capture.

There are a variety of image format available (they are **bmp**, **jpg**, **png** and **tif**) to save the captured image. For example, when choosing **jpg** format, you can set the parameters of "**Option**" to adjust its compression quality or encoding method. Please check the **File>Save As...** menu for details.


10.3 Start Record... F9

1. You can a) choose the **Capture>Start Record...** command; b) click "**Record**" button  on the **Camera Sidebar** ; c) use the shortcut key "**F9**" to start recording movies. The video format can be **wmv/asf**(recommended) or **avi**;


Help Manual

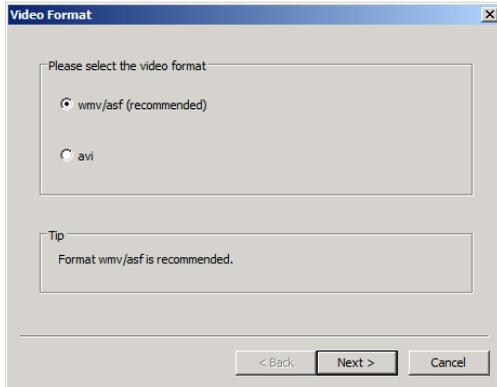


2. Clicking **Next** on the **Video Format** dialog will bring up **Video File** dialog as below; Enter the video file name under **1. Set the name for the captured video file** field and click the **Browse...** button under **2 Select the directory for the video file** item to locate the video file directory. Click **Back** to return to the **Video Format** dialog, or **Next** to the next step;

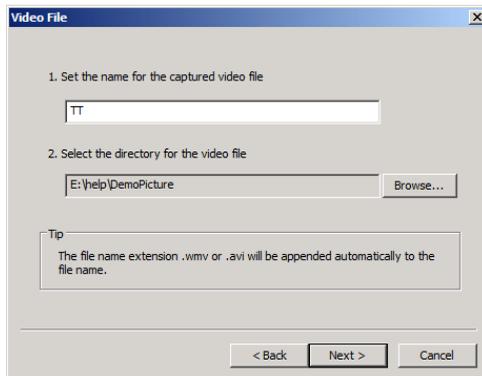


3. An **Encoder** dialog will be brought up. Here you can select the **Encoder** format, set the **Bitrat**(Kbps)(256-16384), **Quality** (1-100) and **Key Frames Spacing** (1-30). Click **Back** to return to the **Video Format** dialog, or **Next** to the next step;

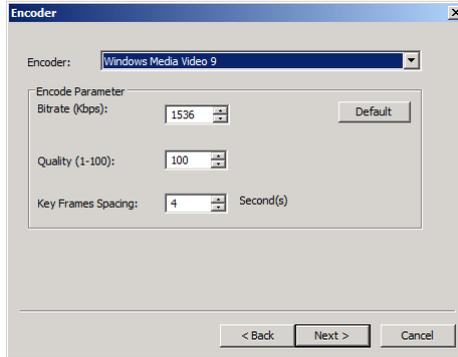

4. A dialog called **Display Information** will be brought up. Here you can enter **Title**, **Author**, **Copyright** and **Description** to their fields. Click **Back** to return to the **Encoder** dialog, or **Next** to the next step;

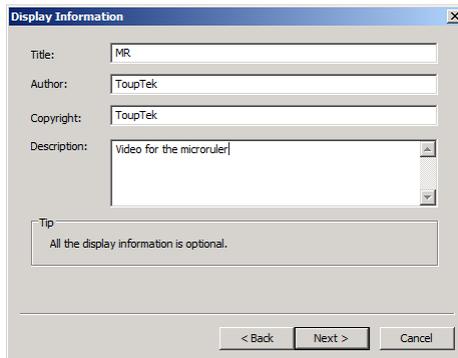


4. A dialog called **Start to Capture** will be brought up. Here you can check **Time Limit (Minutes 1-1440)** and enter recording time (If checked); Input **Time-lapse(1-100)**. There is a **Summary** text to display what you have been defined. Click **Back** to return to the **Encoder** dialog, or **Finish** to end the setup;

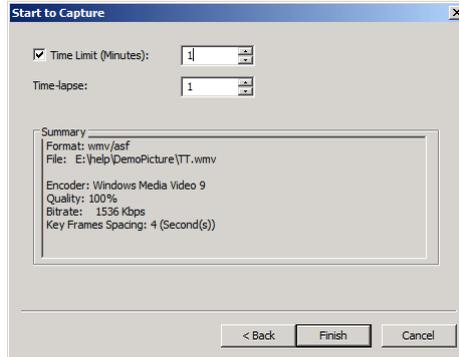

5. After the **Video** capture is started. The  button on the **Camera Sidebar** will become . Clicking  will stop the capture process, otherwise, it will stop until the **Time Limit** is reached. After the **Video** capture process is finished. The  on the **Camera Sidebar** will become  for the future **Record** process;

6. You can choose the **File>Open Video...** command to display the captured video file in the video window;

7. Clicking on the video window with your right mouse button will bring up a context menu. It includes the **View Properties...** menu to check the **Video Properties**, the **Start/Pause** menu to start or pause the video and the **Full Screen** menu to display the video in full screen mode. These commands can also be performed by choosing the **Setup>Start/Pause**, **Setup>View Properties** or **Setup>Full Screen** menu.

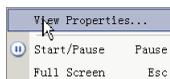


11 Image

11.1 Mode

11.1.1 Color Quantize●●●

The **Color Quantize●●●** command is widely used to change the image bit. ToupView supports the mutual changes among **24 bit**, **8 bit**, **4bit** or **1 bit** images.

When the **Colour Bits** dialog is opened, the default checked color bit is the current image's color bit. Check the desired bit and click **OK** to end the command. The image will be converted to the selected color bits in the image window.

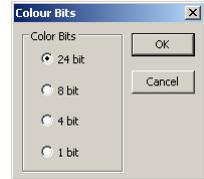


11.1.2 Gray Scale

Choose the **Gray Scale** command to convert a color image (true color image or index color image) to a gray scale image. If the original image is 24 bit, the new image is 8 bit. Otherwise the bit of the image will not be modified.

11.2 Adjust

11.2.1 Curve●●●

Choose the **Image>Adjust>Curve●●●** command to adjust the entire tonal range of an image. But instead of making adjustments using only three variables (highlight, shadow, midtone), one can adjust any point on the curve along a 0-255 scale while keeping up to any other values constant. One can also use **Curve** to make precise adjustments for individual color channels on an image. The **Curve** dialog is shown below:








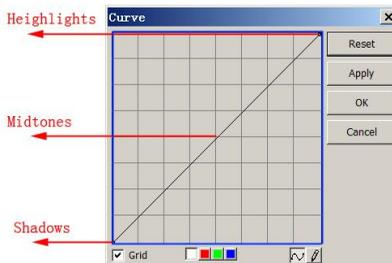


Horizontal	The horizontal axis of the graph represents the original intensity values of
-------------------	--

Help Manual

Axis	the pixels (Input levels).
Vertical Axis	The vertical axis represents the new color values (Output levels). In the default diagonal line, all of the pixels have identical Input and Output values.
Curve area	<p>Area (specified with the blue rectangle) to setup a map between input and output;</p> <ul style="list-style-type: none">a) When you open the Curve dialog you see, perhaps surprisingly, not a curve but a straight line. As a default, the input values are the same as the output, so there's no change;b) By clicking in the center of the line and dragging upwards, we make our first curve. Raising the curve increases the overall brightness of the scene.c) Conversely, clicking in the center and dragging downwards lowers the brightness, producing an image that's darker overall;d) As well as clicking in the center, we can also adjust the endpoints. Clicking the top right point and dragging down limits the brightest part of the image, reducing contrast.e) If we drag that top right point to the left rather than down, we produce the opposite effect—increasing the contrast of the image. This is a very useful and controllable quick fix;f) By dragging the top and bottom points towards the center, we create a stylized, posterized effect that turns any photograph into more of a graphic object.g) Click once in the center of the RGB curve to "pin" that midpoint; now drag just the top half of the curve to make this S shape, and the result is to increase the overall contrast.h) You can add up to any control points to the curve, locking those values. To remove a control point, 1)drag it off the graph or 2)click the right mouse button on that point. You cannot delete the endpoints of the curve.

Help Manual

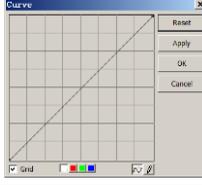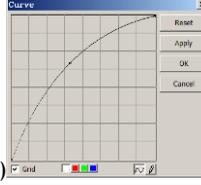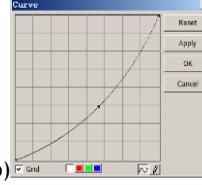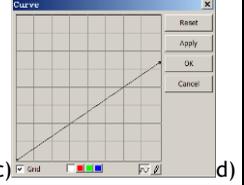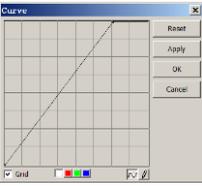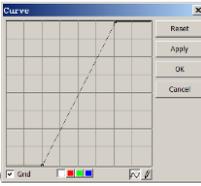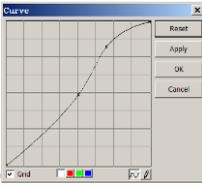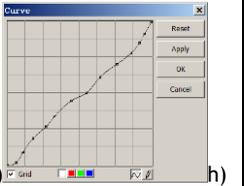
			
			
Grid	Check to overlay the grid on the Curve window;		
Channel 	To adjust the color balance of the image, check the channel (R , G or B) from the Channel button. Check the white button to select RGB channels at the same time, which is located on the left of the R (Red), G (Green) and B (Blue) buttons;		
Curve 	Click the Curve with your left mouse button, a spline curve (straight line) will setup the initial input-output relation. Click on the Curve will mark the Curve with a point. The point number can be any as long as it can be inserted on the Curve . Drag the marked point until the image looks satisfactory. The point can be deleted by move the mouse on it and click the right mouse button on it.		
Pencil 	Check the pencil button at the bottom of the dialog, and use your left mouse button to draw a new arbitrary Curve in the Curve area.		
Reset	Click the Reset button will reset the Curve to the initial straight line.		
Apply	Apply current Curve map on the image.		

11.2.2 Auto Level

The **Image>Adjustment>Auto Level** command moves the level's sliders automatically to set highlight and shadow. It defines the lightest and darkest pixels in each color channel as white and black and then redistributes the pixels' color values proportionately. Since **Auto Level** adjusts each color channel individually, it may remove or introduce color casts. The **Auto Level** command moves the level's sliders automatically to set highlight and shadow. It defines the lightest and darkest

pixels in each color channel as white and black and then redistributes the pixels' color values proportionately. Since **Auto Level** adjusts each color channel individually, it may remove or introduce color casts.

By default, this feature clips the white and black pixels by **0.5%**--that is, it ignores **0.5%** of the lightest pixels and **0.5%** of the darkest pixels when identifying the lightest and darkest pixels on the image. Choose the **Options>Auto Correction...** command to modify this default setting. This ensures that white and black values are representative without being determined by extreme pixel values. The **Auto>Correction...** dialog is shown below:

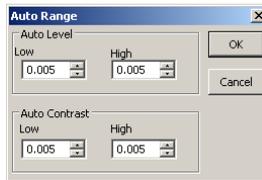

The **Image>Adjustment>Auto Level** command gives good results when an image with an average distribution of pixel values needs a simple contrast adjustment or when an image has an overall color cast. However, adjusting the **Curves** manually is more precise.


See **Image>Adjustment>Auto Contrast** for another auto adjust command.

11.2.3 Auto Contrast

The **Image>Adjustment>Auto Contrast** command automatically adjusts the overall contrast and mixture of colors in an RGB image. Since it does not adjust channels individually, **Auto Contrast** does not introduce or remove color casts. It maps the lightest and darkest pixels in the image to white and black, which makes highlights appear lighter and shadows appear darker.

When identifying the lightest and darkest pixels on an image, **Auto Contrast** clips the white and black pixels by **0.5%**--that is, it ignores the first **0.5%** of either extreme.

Choose the **Options>Auto Correction...** menu to modify this default setting. This ensures that white and black values are representative without being determined by extreme pixel values. The **Auto Correction...** dialog is shown below:

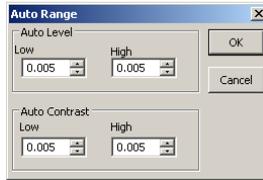

The **Auto Contrast** command can improve the appearance of many photographic or continuous-tone images. It does not improve flat-color images.


See **Image>Adjustment>Auto Level** for another auto operation.

11.2.4 Histogram Equalization

Histogram Equalization is a kind of histogram process. The histogram can reflect the statistical information for the **R, G, and B** of the pixels of the original image. The algorithm calculates each separately, equalizes the **R, G, and B** of the points linearly, and reassigns them

11.2.5 Brightness/Contrast●●●

The **Image>Adjustment>Brightness/Contrast●●●** command offers simple adjustments to the tonal range of an image. This command makes the same adjustment to every pixel in the image. The **Brightness/Contrast** command does not work with individual channels and is not recommended for high-end output because it can result in the loss of details about the image.

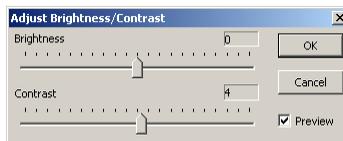

Preview: Check this button to display real-time effects when drags the slider bar. Default is checked

Brightness: Dragging the slider bar to the left decreases the level and dragging it to the right increases the level. The numbers on the right of the slider bar displays the **Brightness** value. Values can range from **-100** to **+100**. Default is **0**;

Contrast: Dragging the slider bar to the left decreases the level and dragging it to the right increases the level. The numbers on the right of the slider bar displays the

Contrast value. Values can range from **-100** to **+100**. Default is 0;

11.2.6 Color●●●


Choose the **Image>Adjustment>Color●●●** command to modify the overall mixture of the colors in an image.

There are four color modules:

RGB	<p>ToupView uses the RGB model. It assigns an intensity value to each pixel ranging from 0 (black) to 255 (white) for each of the RGB components in a color image.</p> <p>For example, a bright red color might have an R value of 246, a G value of 20, and a B value of 50. When the values of all three components are equal, the result is a shade of neutral gray. When the value of all components is 255, the result is pure white; when the value is 0, pure black.</p> <p>RGB images use three channels to reproduce up to 16.7 million colors on-screen. In addition to being the default mode for new ToupView images, the RGB mode is used by computer monitors to display colors. This means that when working in color modes other than RGB, such as CMYK, ToupView uses RGB mode for display on-screen.</p> <p>Although RGB is a standard color mode, the exact range of colors represented can vary, depending on the application or display device.</p>
CMYK	<p>The CMYK mode is based on the light-absorbing quality of ink printed on papers. As white light strikes translucent inks, certain visible wavelengths are absorbed while others are reflected back to the eyes.</p> <p>In theory, pure cyan (C), magenta (M), and yellow (Y) pigments should combine to absorb all light and produce black. For this reason these colors are called subtractive colors. Because all printing inks contain some impurities, these three inks actually produce a muddy brown and must be combined with black (K) ink to produce a true black. (K is used instead of B to avoid confusion with blue.) Combining these inks to reproduce color is called four-color process printing.</p> <p>The subtractive (CMY) and additive (RGB) colors are complementary colors. Each pair of subtractive colors creates an additive color, and vice versa.</p>

Help Manual

HSI	<p>Based on the human perception of color, the HSI model describes three fundamental characteristics of colors:</p> <p>Hue is the color reflected from or transmitted through an object. It is measured as a location on the standard color wheel, expressed as a degree between 0° and 360°. In common use, Hue is identified by the name of the color such as red, orange, or green.</p> <p>Saturation, sometimes called chroma, is the strength or purity of the color. Saturation represents the amount of gray in proportion to the hue, measured as a percentage from 0% (gray) to 100% (fully saturated). On the standard color wheel, Saturation increases from the center to the edge.</p> <p>Intensity is the relative lightness or darkness of the color, usually measured as a percentage from 0% (black) to 100% (white).</p>
HLS	<p>The HLS model is very similar to the HLS color model. The main difference between them is the calculation used to produce the brightness value. In the HLS model, a pixel's brightness (L) is derived from its three (R, G and B) color values. In the HLS model, a pixel's brightness (L) is determined by the minimum and maximum values of its three color values.</p>

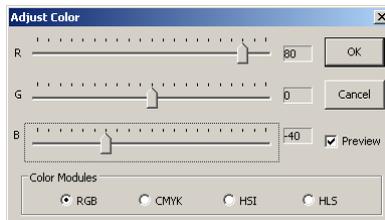

Preview: Check this button to display the real-time effect when the slider bar's position is modified.

The values beside the slider bar show the color changes in various color channels.

For **RGB** channel values, they can range from **-100** to **+100**. Default is 0;

For **CMYK** channel values, they can range from **-100** to **+100**. Default is 0;

For **HSI** channel values, the **H** value can range from **-180** to **180**, the **S** value can range from **-275** to **275**, and the **I** value can range from **-442** to **442**. Default is 0;


For **HLS** channel values, the **H** value can range from **-180** to **180**, the **L** value can

range from **-100** to **100**, and the **S** value can range from **-100** to **100**. Default is **0**;

11.2.7 HMS...


Choose the **Image>Adjustment>HMS...** command to adjust the **HL** (Highlight), **M** (Midtone), and **S**(Shadow) parts of the image. Each part's value ranges from **-100** to **100**. This command is only available for 24 bits true color image.

Preview: Check this button to display the real-time effect when one changes the slider bar's position. Default is checked.

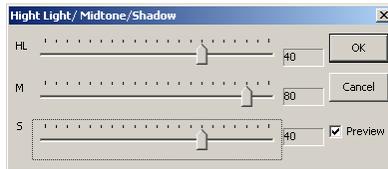

11.2.8 Gamma...

The **Image>Adjustment>Gamma** measures the brightness of midtone values produced by a device (often a monitor). A higher gamma value yields an overall darker image.

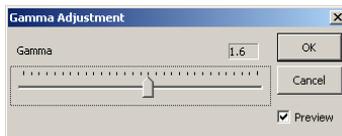

Preview: Check this button to display the real-time effects when one changes the slider bar's position. Default is checked;

Gamma: Dragging the slider bar to the left decreases the level, while moving it to the right increases the level. Values can range from **0** to **3.0**. Default is **1.0**.

11.2.9 Filter Color...

Choose the **Image>Adjustment>Filter Color** command to filter a special color channel from a color image. Check either **Red**, or **Green**, or **Blue** color to filter. For every pixel, if check **Red** color to filter, only information about the red channel will be discarded, and **Green** and **Blue** information will remain there.



See [Extract Color...](#) for another color operation.

11.2.10 [Extract Color...](#)

Choose the [Image>Adjustment>Extract Color...](#) command to extract a special color channel from a color image. Check either [Red](#) or [Green](#), or [Blue](#) color to extract.

For every pixel, if selecting [Red](#) color to extract, only information about the red channel will be kept, and [Green](#) and [Blue](#) information will be discarded.



See [Filter Color...](#) for another color operation.

11.2.11 [Invert](#)

Choose the [Image>Adjustment>Invert](#) command to reverse the pixel values of the active image without going through the lookup table.

11.3 [Rotate](#)

Choose the [Image>Rotate](#) command to rotate the entire image. One has the following submenus:

11.3.1 [90\(CW\)](#)

Rotate the image clockwise by a quarter-turn.

11.3.2 [180\(CW\)](#)

Rotate the image clockwise by 180 degrees.

11.3.3 [270\(CW\)](#)

Rotate the image clockwise by 270 degrees.

11.3.4 Arbitrary•••

Rotate the image by a specified angle. If choosing this option, enter an angle between 0 and 360 degrees in the angle text box, and check **CW** or **CCW** to rotate clockwise or counterclockwise. For the **Arbitrary•••** operation, it will invoke an **Image Rotate** dialog like below:



Degree: The degree that the image to be rotated.

CW: Rotates the image clockwise.

CCW: Rotates the image counterclockwise.

Quality: One can select one of the three methods for the image rotation among **Nearest Neighbor**, **Bilinear**, and **Bicubic**. The default is **Bilinear**.

11.3.5 Flip Horizontal

Reverses the image in the application area so that the top right corner of the original image is now the top left, and the top left corner of the original image is now the top right corner.

11.3.6 Flip Vertical

Reverses the image in the application area so that the top right corner of the original image is now the bottom right corner, and the top left corner of the original image is now the bottom left corner.


11.4 Crop **Shift+C**

Choose the **Image>Crop** command to remove the portions of an image that does not want so that the focus is on the part of the image that is left. This document instructs users on how to **Crop** an image in ToupView.

Crop Demo

1. Choose the **Open Image•••** command to open an image to **Crop**;

2. Choose the **Edit> Image Select** command or click the **Image Select** button  on the **Toolbar**, the cursor will change to a small cross.
3. Move the cursor over the image to the desired location, click the mouse button and hold it down to mark the first position.
4. Drag the mouse over the part of the image to be kept, a dotted rectangle appears around the selection and release the left mouse button.
5. **Optional 1:** To move the rectangle: a) Move the mouse over the selected area and when it becomes a move cursor, click and hold the left mouse button; b) Drag the selected area to the desired position.

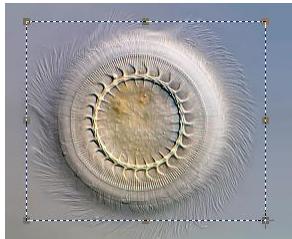

6. **Optional 2:** To change the size of the rectangle: a) Put the mouse cursor on one of the handles that appear on the edges of the selected area; b) Click and hold the mouse button; c) Drag the box to the desired size.

Note: Each of the handles that appear on the edges of the box sizes the box differently.

7. To **Crop** the image, choose the **Image>Crop** command or press **Shift+C** buttons.


Note: If the selected area is on the video, choosing the **Image>Crop** command will capture the selected area and display it on a new image window.

11.5 Image Scale●●●

Choose the **Image>Image Scale●●●** command to change the image to a specified size. This process actually changes spatial resolution by adding (replicating) or removing (decimating) pixels to achieve the specified dimensions.

Width and **Height:** When choosing the **Image Scale** command; the dialog displays

the dimensions of the original image in pixels. The **Width** and the **Height** can be set on the new image by adding or removing pixels. If **Constrain Proportions** is checked, the **Width** and **Height** will stay proportionate to each other. If **Constrain Proportions** is unchecked, the **Width** and the **Height** can set independently, but this will distort the image.

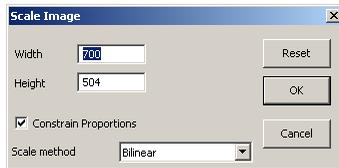

Reset: Reset the image **Width** and **Height** to the original ones.

Constrain Proportions: To maintain the current proportions of pixel **Width** and **Height**, check **Constrain Proportions**. This option automatically updates the **Width** as the **Height is modified**, and vice versa. Otherwise, uncheck the **Constrain Proportions** button.

Scale method: There are 3 options for the **Scale method**. They are: **Nearest Neighbor**, **Bilinear**, and **Bicubic**. The default is **Bilinear**.

11.6 Histogram●●●

A **Histogram** illustrates how pixels in an image are distributed by graphing the number of pixels at each color intensity level. The **Histogram** shows whether the image contains enough detail in the **Shadows** (shown in the left part of the **Histogram**), **Midtones** (shown in the middle), and **Highlights** (shown in the right part) in order to make a good correction.


The **Histogram** also gives a quick picture of the tonal range of the image, or the image key type. A low-key image has detail concentrated in the shadows, a high-key image has detail concentrated in the highlights, and an average-key image has detail concentrated in the **Midtones**. An image with a full tonal range has a number of pixels in all areas. Identifying the tonal range helps determine the appropriate tonal corrections.

Choose the **Image>Histogram●●●** command to open the **Histogram** dialog as shown

below.

Depending on the image's color mode, choose **R**, **G** and **B**, or **Luminosity** to view a composite **Histogram** of all the channels.

If the image is **RGB** true color, choose **Luminosity** to display a **Histogram** representing the luminance or intensity values of the composite channel.

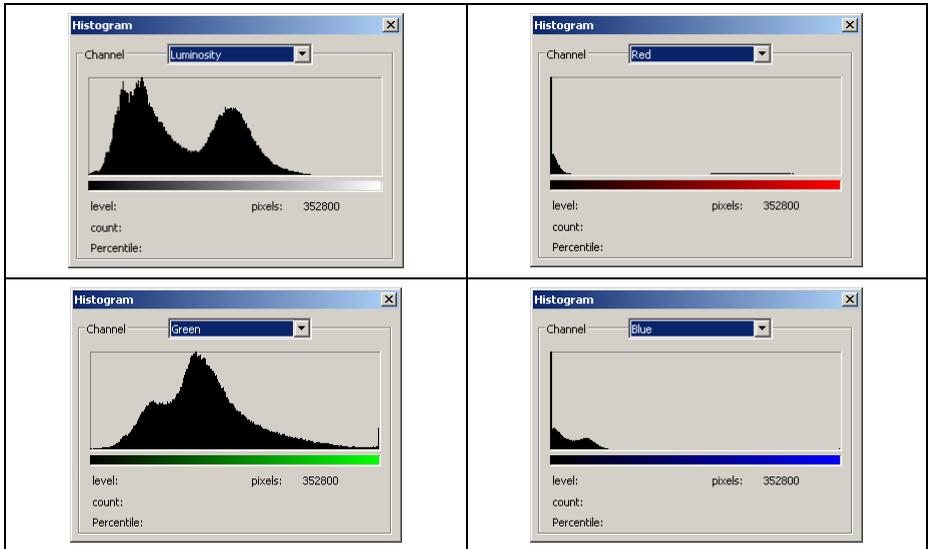

If the image is **RGB** true color, choose **R**, **G** and **B** to display a composite **Histogram** of the individual color channels in color.

Do one of the following:

To view information about a specific pixel value, place the mouse pointer in the **Histogram**.

To view information about a range of values, click down the left mouse button and drag it in the **Histogram** to highlight the range.

The dialog displays the following statistical information below the **Histogram**:

Pixels: Represents the total number of pixels used to calculate the **Histogram**.


Level: Displays the intensity level of the area underneath the pointer.

Count: Shows the total number of pixels corresponding to the intensity level underneath the pointer.

Percentile: Displays the cumulative number of pixels at or below the level underneath the pointer. This value is expressed as a percentage of all of the pixels in the image, from **0%** at the far left to **100%** at the far right.

11.7 Resolution●●●

Choose the **Image>Resolution●●●** command to set the image **Resolution** to calibrate the spatial scale. By default, ToupView expresses spatial measurements in terms of pixels. This **Resolution** command is used to change the terms in which ToupView reports such measurements. This command should be run first in order to measure objects in terms of units other than pixels.

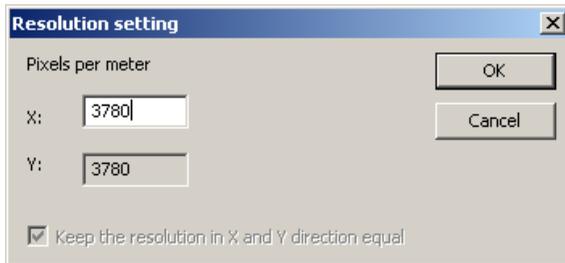

X: Horizontal **PPM (Pixels per meter)** of current **Resolution**.



Y: Vertical **PPM (Pixels per meter)** of current **Resolution**.

Note: The resolution in the **Y** direction need not to be filled. ToupView will always let it equal to the **X** direction.

After the new **Resolution** is set, all of the measurements will be calculated according to the new **Resolution**. (See more in the **Measurements and Options>Measurement●●●** menu).

11.8 Overlay Scale Bar●●●

The **Image>Overlay Scale Bar●●●** command will overlay a scale bar on an image. The scale bar **Position** can be **Top Left**, **Top Right**, **Bottom Left** or **Bottom Right**. The scale bar can be in **Auto** or **Fixed** mode. The **Font Size** and **Font Weight** can be chosen. Finally the scale bar **Color** and **Font Color** can be set

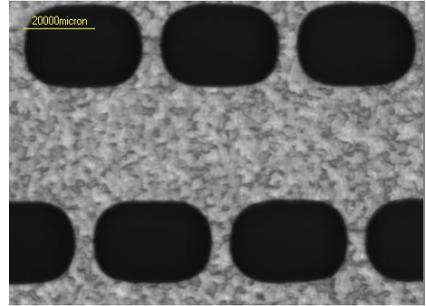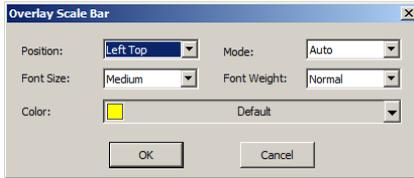


12 Process


12.1 Stitch

12.1.1 Video Stitch

When the video window is active, choosing the **Process>Stitch** command will invoke the video **Stitch** dialog as below:

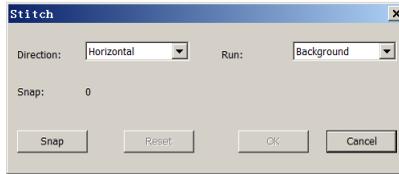

Direction	The image stitching direction, Horizontal or Vertical ; The default is Horizontal ;
Run	The stitching run method: Background or Foreground . The default is Background ;
Snap:	The captured images from the camera;
Snap	Click to capture image from camera;
Reset	When an image or images is or are captured, the Reset button will be enabled, click Reset will remove it or them from the memory; It will be disabled when there is no captured image;
OK	Click Ok to begin image stitching. This button will be enabled only when the snapped images are more than 2;
Cancel	Click to cancel the stitching process;

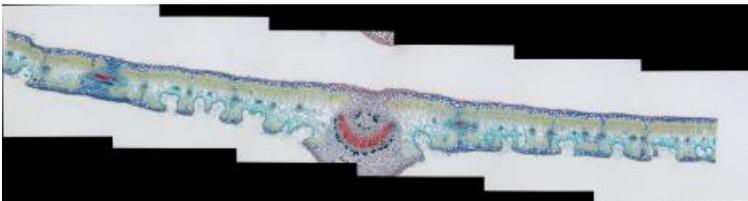


Stitching finished image

In the stitched image, there are some black areas. These are happened when there are translations between the images in both directions.

12.1.2 Image Stitch

Help Manual

When the image window is active, choosing the **Process>Stitch** command will invoke the image **Stitch** dialog

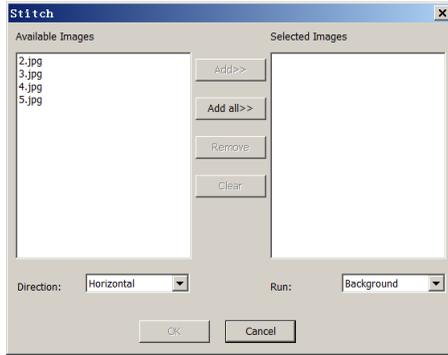


Available Images	The images opened by ToupView;
Selected Images	The images selected for stitching;
Add>>	When an image or images is or are selected and highlighted in the Available Images window, the Add>> will be enable. Click Add>> will add it or them into the Selected Images window;
Add All>>	Click Add All>> will add all of the Available Images into the Selected Images window;
Remove	When an image or images is or are selected and highlighted in the Selected Images window, the Remove button will be enabled, click Remove will remove it or them from the Selected Images window;
Clear	Used to clear all of the images in the Selected Images window. When no image in the Selected Images window, it will be disabled;
Direction	The image stitching direction, Horizontal or Vertical ; The default is Horizontal ;
Run	The stitching run method: Background or Foreground method. The default is Background ;
OK	Click Ok to begin image stitching. Only the Selected Images are more than 2, this button will be enable;


Help Manual

Cancel

Click to cancel the stitching process;

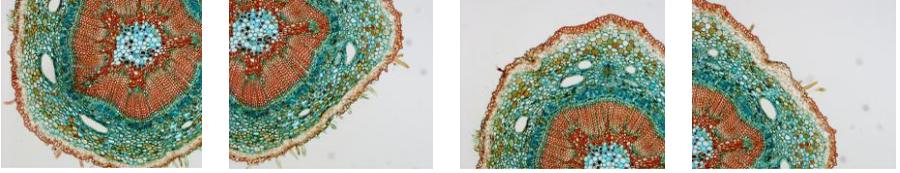

4 images for stitching

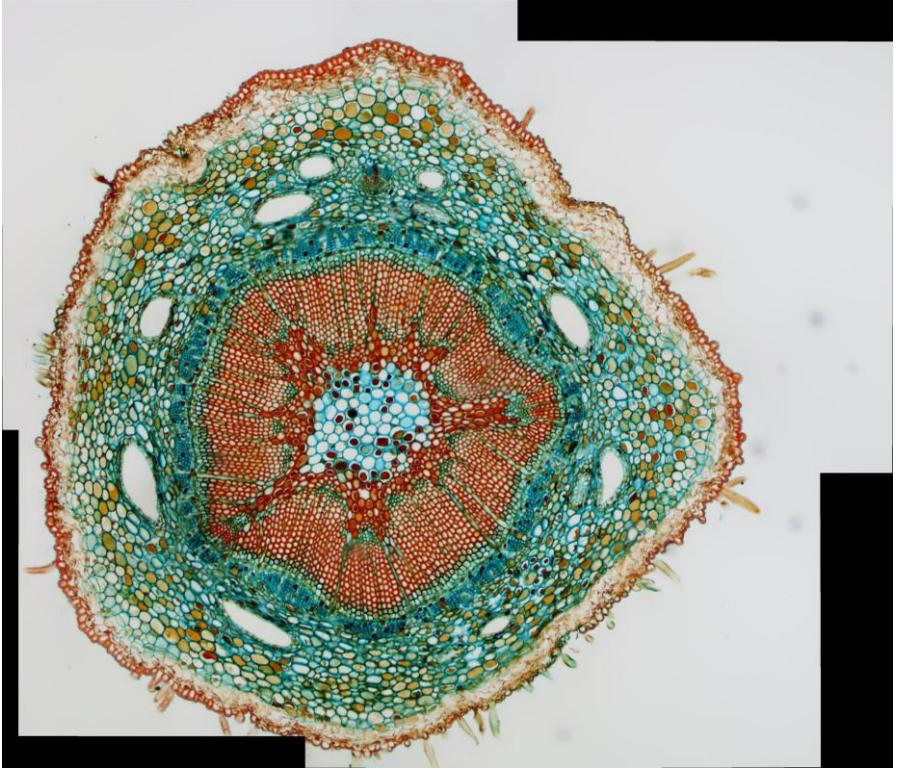

Stitching finished image

In the stitched image, there are some black areas; these are happened when there are translations between the images in both directions.


12.2 Filter●●● Shift+F

Choose the **Process>Filter** command to apply one of ToupView's numerous filter on the active image. If you are not familiar with the process and effects of filtering,

some discussions about spatial filtering should be reviewed. ToupView provides an extensive set of convolution and no convolution (morphological) filters. you can also create custom filter kernels and apply them with the **Filter** commands.

Choosing the **Process>Filter** command will open the **Filter** dialog. Each group of filters has its own property page or tab, where the **Filter** type and size can be selected. The filtered results are almost always written to the active image. The **Edit>Undo** command can be used to remove **Filter** operations that have been applied.

12.2.1 Image Enhance

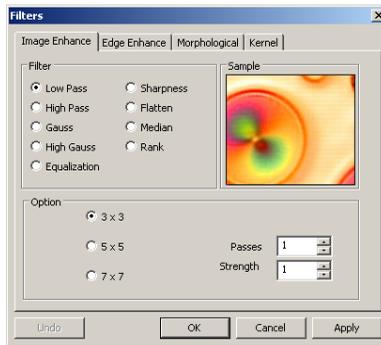

Low Pass	Check this filter to soften an image by eliminating high-frequency information (this has the effect of blurring sharp edges). The Low Pass filter replaces the center pixel with the mean value in its neighborhood. The Low Pass filter can also be used to remove noise.
High Pass	Check this filter to enhance high-frequency information. The High Pass filter replaces the center pixel with a convolved value that significantly increases its contrast from its neighbors. The High Pass filter leaves only elements of high contrast.
Gauss	Check this filter to soften an image by eliminating high-frequency information using a Gauss function. This has the effect of blurring sharp edges. The operation of the Gauss filter is similar to the Low Pass filter, but it degrades the image less than the Low Pass filter.
High Gauss	Check this filter to enhance fine details. This operation is similar to the unsharp masking technique (see the Sharpen filter), but it introduces less noise into the image. It uses a Gaussian curve type of kernel. Available in

Help Manual

	7x7 and 9x9 kernel sizes.
Equalization	This filter is used to enhance the contrast based on the histogram of the local neighborhood (See Option below).
Sharpness	Check this filter to enhance fine details, or refocus an image that is blurred. The sharpen filter sharpens the image using the unsharp masking technique.
Median	Check this filter to remove impulse noise from an image. The Median filter replaces the center pixel with the Median value in its neighborhood. It will also blur the image.
Rank	Check this filter to remove impulse noise from an image. The pixels in the kernel are ranked by order of intensity, and the pixel in that range at the rank percentage is chosen for comparison. For example, in a 5x5 kernel, there are 25 pixels. A rank percentage of 95% would choose second-brightest pixel for comparison. If the difference between the selected pixel and the center pixel is greater than the threshold value, the Rank filter replaces the value of the center pixel with the value of the selected pixel.

Option

1. If one of the **Enhancement** filters is checked, the following options will be displayed:

3 x 3	Check 3 x 3 kernel will produce a more subtle filtering effect.
5 x 5	Check 5 x 5 kernel will produce a moderate filtering effect.
7 x 7	Check 7 x 7 kernel will produces a more extreme filtering effect.
Passes	Set the filter applied times on the image. When a filter is applied multiple times, its effect is amplified by each pass. An image that has been softened by one pass of the Low Pass filter will be softened further by a second pass.
Strength	Enter an applied value from 1-10 that reflects how much of the filtering effect on the image. A value of 10 specifies the full strength (100%) of the filtered result applied to each pixel. Values less than 10 cut the full weight of the filter. A value of 1 indicates that only 10% of the difference between the filtered

Help Manual

	pixel value and the original pixel value should be applied, a value of 2 indicates that 20% of the difference should be applied, and so forth.
Rank	<p>This value specifies which pixel in the sorted array will be used to replace the center pixel. Pixels in the array will be sorted in ascending order. The pixels are indexed from 0 to $\text{Kernel Size} \times \text{Kernel Size} - 1$. In the pixel index 0 corresponds to the lowest pixel value.</p> <p>The Rank will be specified in terms of a percentage of the indexes ($\text{Kernel Size} \times \text{Kernel Size} - 1$). A 50% Rank means the middle of the array. 0% rank means the lowest index (lowest gray value), and 100% rank means the highest index (highest gray value).</p>


2. If the **Equalization** filter is checked, the options will relate to the histogram equalization. **Local Histogram Equalization** modifies the contrast of an image based on the pixel values in a small window surrounding each pixel.

Window	Image pixels statistics (min, max, histogram, mean, standard deviation, etc.) will be calculated on a small Window of the image. These measurements are then used to derive the local contrast for that area of the image. In short, an area of $\text{Window} \times \text{Window}$ around each pixel is all that is considered when modifying the intensities in the image. Larger Window produces smoother results, while small Window track small details more closely.
Best Fit	Choose the Best Fit button to optimize the values for the particular image. The results are achieved by stretching the local histogram to maximize the contrast between the brightest and darkest pixels in the local window region.
Linear	This option distributes the histogram linearly across the intensity scale. This function produces a high contrast image with the highest possible dynamic range.
Logarithmic	This option concentrates the histogram at the low end of the scale. This function produces a high contrast image with little dynamic image. It will tend to darken the image overall. It is useful for increasing the contrast in a very light image.
Exponential	This option concentrates the histogram at the high end of the scale. This

Help Manual

-tial	function produces a high contrast image with little dynamic image. It will tend to lighten the image overall. It is useful for increasing the contrast in a very dark image.
-------	--

12.2.2 Edge Enhance

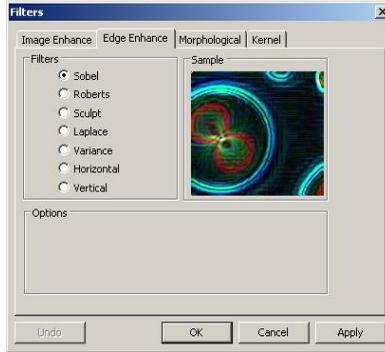

Sobel	Check this filter to enhance just the principal edges in an image. The Sobel applies a mathematical formula to a 3x3 neighborhood to locate and highlight its edges.
Roberts	Check this filter to enhance fine edges in an image. The Roberts filter is not a convolution filter. It applies a mathematical formula upon a 4 x 4 neighborhood to produce its effect. The upper left pixel in the neighborhood is the one that is replaced.
Sculpt	Check this filter to apply a sculpted effect on the image.
Horizontal	Check this filter to detect and emphasize horizontal edges.
Vertical	Check this filter to detect and emphasize vertical edges.

Options

1. If one of the **Edge** filters has been checked, the options will relate to kernel size and filtering strength. The following options will be displayed:


3 x 3	Check 3x3 kernels to produce a more subtle filtering effect.
5 x 5	Check 5x5 kernels to produce a moderate filtering effect.
7 x 7	Check 7x7 kernels to produce a more extreme filtering effect.
Passes	Enter the number of times that the filter will be applied to the image. When a filter is applied multiple times, its effect is amplified by each pass. An

Help Manual

	image that has been softened by one pass of the Image Enhancement Filter , will be softened further by a second pass.
Strength	Enter a value from 1-10 that reflects how much of the filtering effect to apply to the image. A value of 10 specifies that the full strength (100%) of the filtered result will be applied to each pixel. Values less than 10 cut the full weight of the filter - a value of 1 indicates that only 10% of the difference between the filtered pixel value and the original pixel value should be applied, a value of 2 indicates that 20% of the difference should be applied, and so forth.

2. If **Sobel** or **Roberts** is checked, no options are available.

12.2.3 Morphological

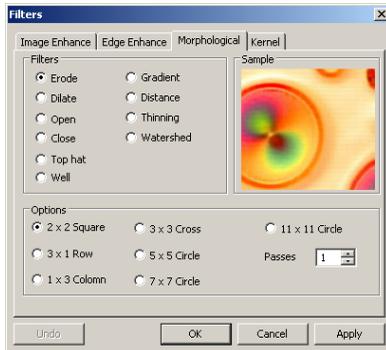

Erode	Check this morphological filter if one wants to modify the size of objects in the image. The Erode filter erodes the edges of bright objects and enlarges the edges of dark ones.
Dilate	Check this morphological filter if one wants to modify the size of objects in the image. The Dilation filter dilates bright objects and erodes dark ones.
Open	Check this morphological filter if one wants to modify the shape of objects in the image. Assuming the image contains bright objects on a dark field, the Open filter will smooth object contours, separate narrowly connected objects, and remove small dark holes.
Close	Check this morphological filter if one wants to modify the shape of the objects in the image. Assuming the image contains bright objects on a dark field; the Close filter will fill gaps and enlarge protrusions to connect

Help Manual

	objects that are close together.
Tophat	Check this filter to detect and emphasize points, or grains, that are brighter than the background. There are 3 kernel sizes for this processing. Click the radio button to change the kernel size to the value that most closely matches the size of the grains to detect.
Well	Check this filter to detect and emphasize points, or grains, that are darker than the background. There are 3 kernel sizes for this processing. Click the radio button to change the kernel size to the value that most closely matches the size of the grains to detect.
Gradient	Check this filter to enhance edges in an image.
Watershed	Check this filter to separate objects that are touching. The Watershed filter erodes objects until they disappear, then dilates them again, but will not allow them to touch. The Watershed filter will not operate upon True Color images. If one wants to separate objects in a True Color image, he must first convert it to Gray Scale (see Process Frame: Image>Gray Scale).
Thinning	Check this filter to reduce an image to its skeleton. When choosing this filter, one must set the threshold that determines whether a pixel is part of the subject or part of the background (see Options below). The Thinning filter will not operate upon True Color images. If one wants to thin a True Color image, he must first convert it to Gray Scale .
Distance	The distance filter is used to show the distances of pixels within blobs to the outer boundaries of those blobs. After applying the distance filter, the background will be black (i.e. pixels with value 0). Only the area within the blobs will have non-zero values (will be white). The values of each pixel within the blob will be a count of the shortest distance from that pixel to the edge of the blob. Thus, all pixels along the blob's border will have a value of 1 (since they are one pixel away from the edge of the blob); pixels that are a distance of 2 from the border will have the value 2, and so on. This creates a distance map of the image. The Distance filter will not operate upon True Color images. If one wants to use the Distance filter with a True Color image, he must first convert it to Gray Scale .

Options

1. If **Erode**, **Dilate**, **Open**, or **Close** filters is checked, the options will relate to the

Help Manual

kernel size and shape. The following options will be presented:

2 x 2 Square	Check to use the 2x2 square kernel configurations.
3 x 1 Row	Check to use the 3x1 row kernel configuration.
1 x 3 Column	Check to use the 1x3 column kernel configuration.
3 x 3 Cross	Check to use the 3x3 cross kernel configuration.
5 x 5 Circle	Check to use the 5x5 circular kernel configurations.
7 x 7 Circle	Check to use the 7x7 circular kernel configurations. This is a two-pass filter, accomplished using a 5 x 5 circle followed by a 3x3 cross.
11 x 11 Circle	Check to use the 11 x 11 circular kernel configurations. This is a three-pass filter, accomplished using a 5 x 5 circle followed by another 5 x 5 circle, followed by a 3 x 3 cross.
Passes	Set the number of times iterate the filter.

Note: The circular kernels are especially effective on round objects (cells, grains and so on) because their circular configuration preserves the circular shape of the objects better than square configurations.

2. If the [Tophat](#), [Well](#), or [Gradient](#) filter is selected, the options will relate to kernel size and shape. The following options will be presented:


3 x 3	Check to use the 3x3 square kernel configurations.
5 x 5	Check to use the 5x5 square kernel configurations.
7 x 7	Check to use the 7x7 square kernel configurations.

3. If [Watershed](#), [Thinning](#), or [Distance](#) filter is checked, the options will relate to the threshold. The following option will be presented:

Threshold	Enter a percentage value from 1-100 that specifies the intensity value to binarize the image. For example, a Threshold of 50% on a Gray Scale image would set all values ≤ 127 to 0 (black) and all values ≥ 128 to the maximum value for that image class (white).
-----------	---

12.2.4 Kernel

The [Kernel](#) page allows edit the kernel files for the morphological and convolution filters.

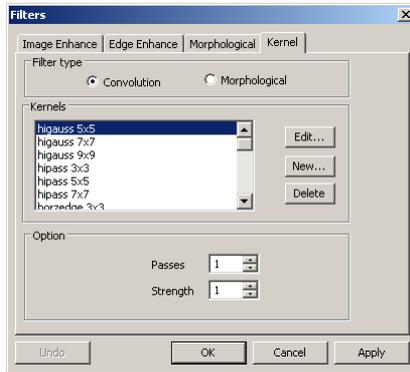

Note: The **HiPass**, **LowPass**, **Laplacian** and **Unsharp** kernel files are used by the **HiPass**, **LowPass**, **Laplacian**, and **Sharpen** options listed in the **Image Enhancement Filters** page dialog window (i.e. there is no difference between selecting one of these kernel files and selecting its **Option** button in the **Filter** window -- the two methods ultimately do the same thing). Because these kernel files are essential to the operation of these filtering options, they must not be deleted or renamed.

Filter type	Check to modify the kernel for a selected Filter type , either Convolution or Morphological filters.	
Edit...	Check to modify the selected filter kernel using the Edit Kernel dialog.	
	Name...	This combobox contains the name of the selected kernel file. If one wants to save the modified kernel file to the same file, leave it as it is. If one wants to save the file to a new location, enter the new filename here.
	Kernel Size	Click the spin buttons or enter the number to change the size of the kernel. Either direction may take into account one to nine pixels. As one modifies the Kernel Size , the shape of the kernel representation changes accordingly. In the center of the dialog, there are white boxes containing coefficients that will be multiplied with each pixel that will be taken into account by the filter kernel. One can change any coefficient by clicking on it and adjust it as desired.
	Fill	Click this button to fill every element of the kernel with a


Help Manual

		particular value. The Fill kernel dialog appears. One may enter a value between 0 and 10. Using the Fill button is useful for setting all coefficients to the same value. One may then change the coefficients that require a different value.
	Offset.	The pixel whose value is being modified is usually the center-most pixel. One may, however, designate any pixel. ToupView signals the pixel to be changed by putting a box around it. Choose the X and Y Offset spin buttons to apply
New...		Click to create a new filter kernel. The Edit Kernel dialog will appear. The functions of the dialog are the same way as the dialog for Edit... (described above), with the exception that the file name for the new kernel file must be provided.
Delete		Click to delete the selected filter kernel file.
Options		The choices in this group box will vary depending upon the kind of selected filter.

12.3 Range... Shift+R

The **Process>Range...** command allows set the intensity levels of the image to increase the contrast and enhance the display in low-light situations. Choose the **Range** command to open the **Range** dialog.

Two **vertical markers** show the **upper** and **lower** limits of the intensity levels. These markers can be moved with mouse through the drag and drop method. For a color image, the histogram will reflect the **red**, **green**, and **blue** values with corresponding colors lines.

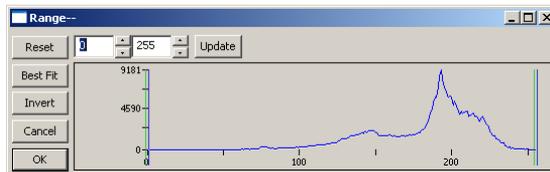

Two **Edit** controls indicate the values of the intensity levels. Choose the spin buttons to increase or decrease these values. All values between 0 and the lower limit will be black and all values between the upper limit and the upper end of the scale will be

Help Manual


white. These two edit control values corresponding to the two **vertical markers** show the **upper** and **lower** limits of the intensity levels. Defaults are 0 and 255 respectively.

Reset	The Reset button allows Reset the black and white levels to the high and low ends of the dynamic Range .
Best Fit	The Best Fit button automatically sets the intensity levels to the Best Fit . Best Fit instructs ToupView to optimize the brightness and contrast values for the particular image.
Invert	The Invert button reverses the color of the image.
Update	Update will refresh the display Range with the most current image information.

12.4 Segmentation••• Shift+S

Segmentation is a process through which certain colors (or gray levels) in an image can be visually identified when they are isolated from the image as a whole. Areas identified by **Segmentation** (classes) can either be removed from or kept in the image, while discarding the remainder of the image. Therefore, this process can be used for separating items or objects of interest from the "background noise" that normally occurs in most acquired images.


The process of identifying colors is the key to the operation of **Segmentation**. Due to the vast possibilities of differences in the images, and the color composition of the object(s) to be identified, ToupView provides **Histogram** based models for identifying the segmented area.

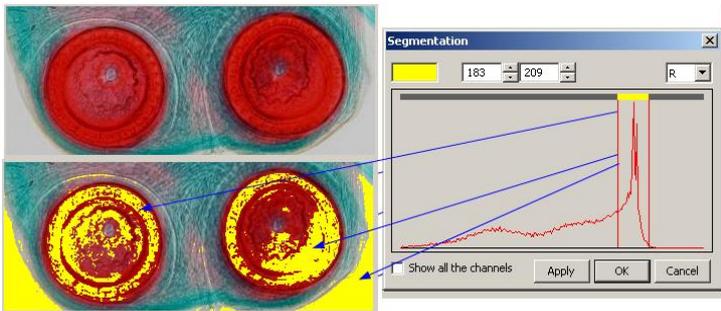

Choose the **Segmentation** command on a **True Color** image, select either **Red(R)**, **Green(G)**, or **Blue(B)** channel to do the operation.

12.5 Binary... Shift+B

Binary is a kind of gray level process. If the gray of the pixel is greater than the given threshold, the pixel's color will be changed into white. Otherwise, the pixel's color will be changed into black. Although the process may lose some information, it is an important step of other processes.

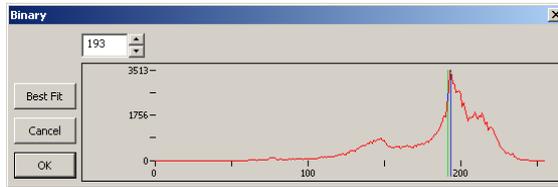

The curve on the **Binary** dialog shows the gray distribution of the image.


The **Line** in the dialog indicates the threshold value. Drag the line to change the threshold, or change the value in the **Line** position **edit** box (in the top left corner of the dialog) to move the **Line**.

The **Line** position **edit** box shows the current position value of the **Line**; Modify the value will move the **Line** at the same time;

Click the "**Best Fit**" button to apply the auto threshold process to the image. It uses an automatic threshold to make the image **Binary**.

12.6 Emboss... Shift+E

Emboss is a kind of artistic process. The process can make the image look like an embossed image. The **Preview** button allows previewing the image before creating it. The process supplies 3 kinds of convolutions including **Gradient**, **Different**, and **Prewitt**.

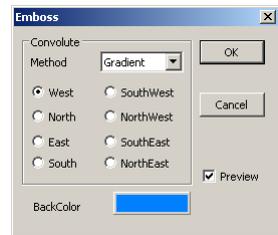

There are 8 directions (**West**, **North**, **East**, **South**, **Southwest**, **Northwest**, **Southeast**, **Noutheast**) in every kind of convolution. Users can get different effects with different convolution methods or directions;


Check **Preview** to preview the processing at real time;

Click on the  to change the **Back Color** of the **Emboss** processing;

12.7 Pseudo Color●●●

Note: Image must be in **Gray Scale**.


Choose the **Pseudo Color** command to "colorize" the active monochromatic image. This is used to highlight certain features in a gray scale image such as display all densities above a certain point in red, or, the imaging device recorded thermal information, all temperatures below a certain point can be revealed in blue color.

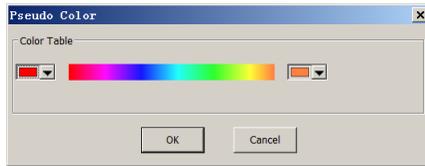

When **Pseudo Color** a monochromatic image, a special palette need to be build with which the monochromatic image is displayed. **Pseudo Coloring** an image does not modify the pixels' values in image bitmap (it does not convert image to true color or palette). It simply associates a **Pseudo Color** palette with the image that interprets the gray-level values in the image as color.

Pseudo Colored images are very similar in structure to palette class images, but they differ in a couple of important ways. First, the pixels' values in a **Pseudo Colored** image actually represent continuous-tone intensity information, whereas a palette image's pixels carry no intensity significance. Secondly, a palette image includes a palette table that is actually part of the image file.


The colors used to map the gray values can be selected. The buttons at each end of the color strips ( or ) will bring up the color dialoges separately, which allows select the start and end colors of the range.

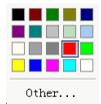

12.8 Surface Plot●●●

The **Process>Surface Plot●●●** (or 3-D Plot) tool creates a 3-D representation of the intensity of an image. When choosing the **Surface Plot** command, keep in mind that X=image width and Y=image height, and Z=pixel gray value.

In the viewpoint window, the **elevation** and **rotation** of the image can be adjusted by dragging the mouse on the image.

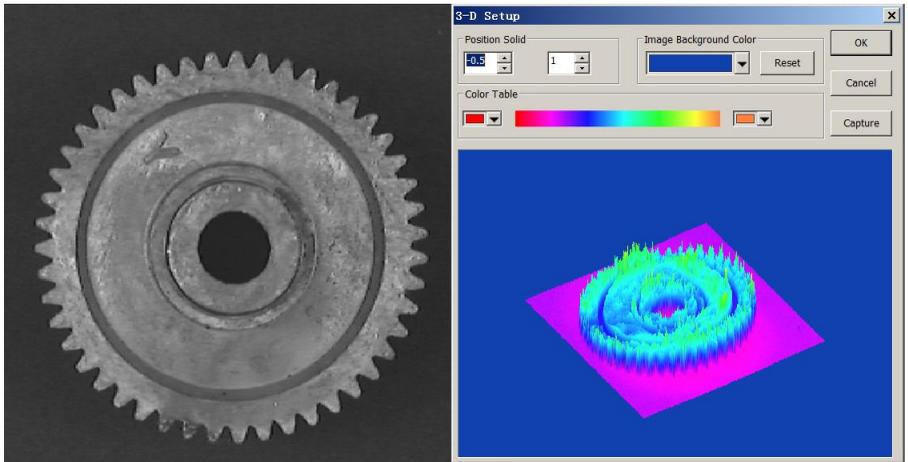

Position Solid: The left edit control indicates the relative position of the entire image in the viewpoint window, whose default value is 0.5. The right edit control indicates the relative height of the display of the Z scale, whose default value is 1.

Reset: Set the **Position Solid**'s two edit controls to their default values.

Image Background Color: Choose this command to display a color dialog where one can adjust the background color of the viewpoint window.

Capture: **Capture** the active image in the viewpoint window as a new image.


Color Table: Select the colors to map the gray values found in the surface plot. The button at each end of the **color table** brings up the color dialog, which allows select the start and end colors of the range. (Please refer to **Pseudo Color** for more information)

12.9 Line Profile●●●

Choose the **Process>Line Profile●●●** command to illustrate how pixels along a selected line are distributed by graphing the number of pixels at each color intensity level.

Choose the **Measurements>Arbitrary Line**, or **Measurements>Horizontal Line** or **Measurements> Vertical Line** command to draw a line on the image and choosing the **Process>Line Profile** will bring up a **Line Profile** dialog as below(The procedure can be

inversed):

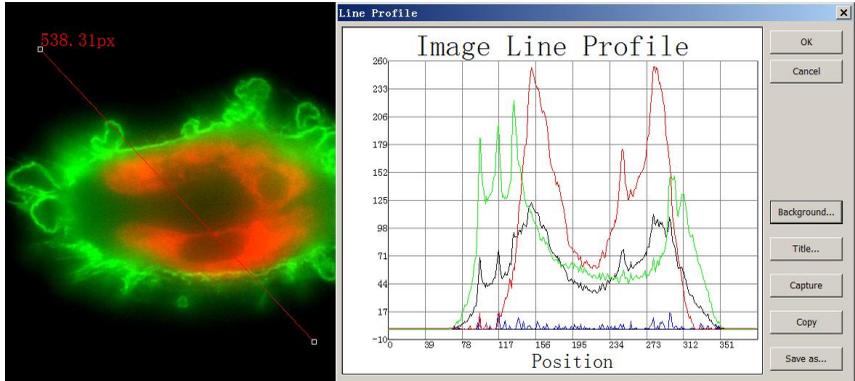

Line Profile: In a **Line Profile**, the X-axis represents the spatial scale, and the Y-axis represents the intensity values which range from 0 to 255. If the image is a gray scale image, the gray value will be profiled. If the image is a RGB color image, the R,G and B values will be profiled separately with the corresponding R, G and B colors

Background: Open the windows color dialog to set the background color of the profile window;

Title: Use this command to set a title on the image's **Line Profile**;

Capture: Capture the image in the **Line Profile** window as a new untitled image;

Copy: Copy the **Line Profile** window's content onto the clipboard;

Save as...: Save the **Line Profile** image in bmp format.

12.10 Diffuse... Shift+D

Diffuse is a kind of artistic process. It can diffuse the image. Users can adjust the parameter in the dialog to



Size: The **Diffuse** level, its range is between 1 and 29. Default is 1, means no diffusing.

Preview: Check it to display the real-time effect when drag the slider bar.

12.11 Granulate••• Shift+G

Granulate is a process that can make the image fuzzy. One can adjust the parameter in the dialog to control the degree of the fuzziness.



Size: The **Granulate** level, its range is between 1 and 20. Default is 1, means no granulating.

Preview: Check it to display the real-time effect when drag the slider bar.

12.12 Mosaic•••

Mosaic is a process that can combine the images opened into a new image. This will open the following dialog:

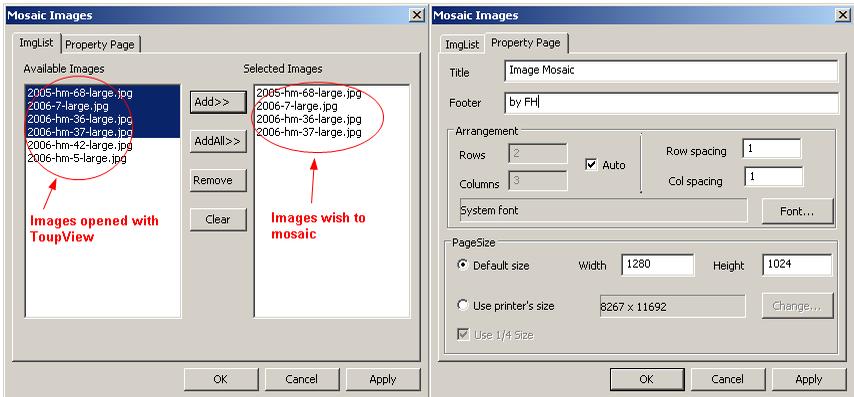


Image List page

Property page

ImgList page

Available Images	Images opened with Toupview;
Add>>	Add the selected images in the Available Images window to the Selected Images window. The Add>> button will be enabled when an image is highlighted;
Add All>>	Add all the opened images to the Selected Images window;
Remove	First select at least an image at the Selected Images window, then the

Help Manual


	Remove button will be enabled and click it to remove it from the Selected Images window. When at least an image is highlighted in the Selected Images window, the Remove button will be enabled;
Clear	Remove all the images from the Selected Images window. This button will be enabled when there is an image in the Selected Images window;

Property page

Title:	The title wanted
Footer	The footer wanted
Arrangement	The Mosaic images space Arrangement on the page. It can be arranged by a)specifying the image Row and Column numbers or b)determining the Row and Column numbers automatically(Check the Auto box; c) The spacing between the Row and Column , its unit is pixel; d) The Font for the arranged file names
PageSize	The page size for the Mosaic images. It can be a) Default , default is unchecked; The Width and Height can be changed. Defaults are 1024 and 1280 respectively. b) Use printer's size : This means the pixel resolutions for the Mosaic image can be calculated from the printer page's size and the printer resolution. For A4 and 300 DPI, the resolution should be c) Use ¼ size : If this box is checked, the size will be ¼ of Use printer's size

If everything is set, click **OK** to end the **Mosaic** operations and a new image window will be displayed and the final results should be


Color Composite Demo

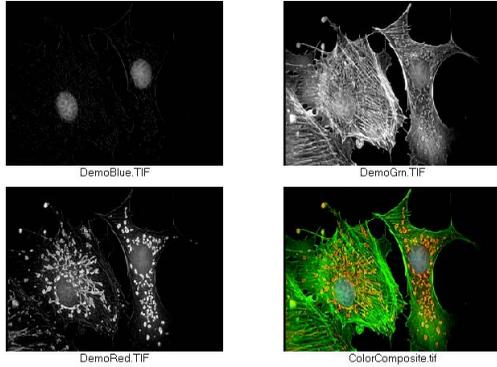

By ToupView Mosaic Processing

12.13 Fusion...

Similar to the dynamic multi-focus image fusion with the live video stream, static **Fusion** is a very useful tool to generate a clear image by combining a sequence of previously captured multi-focus images. Choosing the **Process>Fusion...** command, the following dialog will be brought up (assume 01.jpg 02.jpg ... 15.jpg are already opened in ToupView):

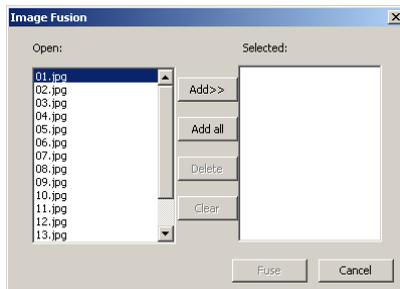

Clicking on the image file name in the **Open** combobox will highlight the image, and then the **Add>>** button is enabled (Click on the selected images will deselect them). Click the **Add>>** button, the highlighted images will be added into the **Selected** list box, which will be fused later.

Clicking the **Add all** button will add all images in the **Open** combobox into the **Selected** list box.

If images in the **Selected** combobox is highlighted, the **Delete** button will be enabled.

Help Manual

Click the **Delete** button, the highlighted images in the **Selected** combobox will be removed.

Clicking the **Clear** button will remove all the images in the **Selected** list box, including the unselected ones. The button will be disabled if there is no image in the **Selected** combobox

NOTE: Images used for fusing must be the same size; otherwise there will be a prompt when adding different sizes of images as below. The prompt dialog indicates which image is not the same size with the others.



When the desired images are all selected (If more than two image in the **Selected** List box, the **Fusion** button will enabled), the **Fusion** process can start. Clicking on the **Fuse** button will start the **Fusion** process, and the mouse cursor becomes an hourglass. When the fusion process finishes, the fused image will be generated in a new image window as below:





Clicking **Cancel** will cancel the **Fusion** processing and return to the application area.



12.14 Color Composite●●●

Help Manual

Use the **Color Composite** dialog box to create and configure color composites using black and white source images. You can access the **Color Composite** dialog box through the **Color Composite** menu item from the **Process** menu.

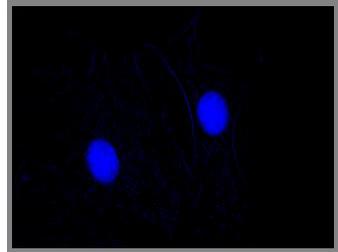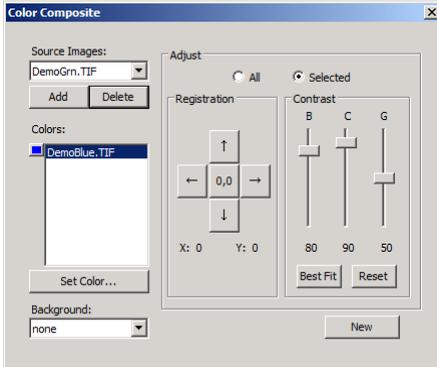

You can combine gray-scale images into a color composite. Any group of gray-scale images that are of the same size can be mixed in a color composite. Images of 8-, 12-, 16-bit integer or floating point format are combined into a 24-bit color composite. Each input channel will have individual LUT adjustments, as well as a registration offset to line it up with the rest of the images.

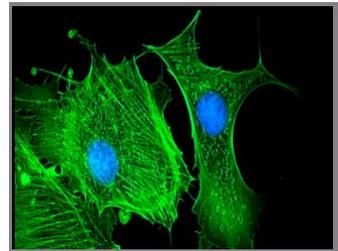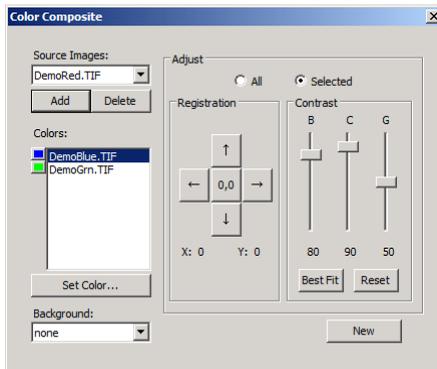

You can also combine individual channels from a single image. Each channel will be listed separately, as shown here:

Source Images: The **Source Images** combo box displays the images available for color mixing. The **Source Image** list is initially filled with a list of the gray scale


Help Manual

images that are currently open. The size of the color composite is determined by the first input image selected. Once this is entered, the list is filtered to contain only images of the same size as the initial selection.

Add: Click the **Add** button to add an image to the mix. You will be prompted for the color to be assigned to that image from the **Set Color** dialog.

Delete: Use this button to delete the selected image from the list.

Colors: The **Colors** combobox displays the images selected for color mixing.

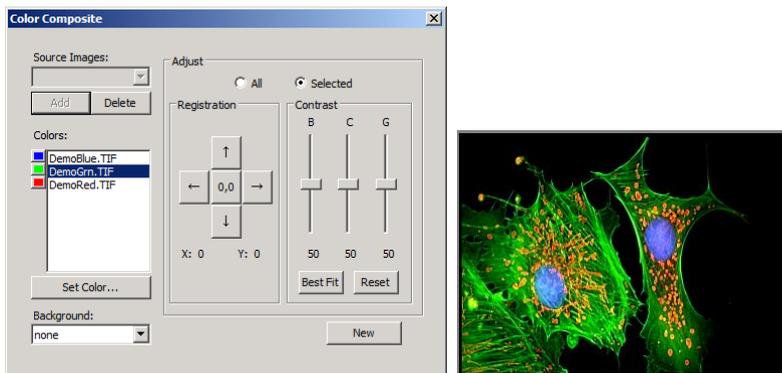


To change the color within the list box, double-click to select the image name (the **Select Color** dialog box appears) or highlight the name and click the **Set Color** button.

To delete colors/images from the **Colors** list box, first click on the color/image to highlight it then press the **Delete** button within the **Color Composite** dialog box.

Color Bar: The picture item box on the left side of the **Colors** list indicates the color associated with each image in the composite. The **Color Bar** will scroll with the contents of the **Colors** list box.

Set Color: This option determines the **Color** value used for the selected channel in the final image display. Clicking on the **Set Color** button brings up the **Select Color** dialog box.

Help Manual

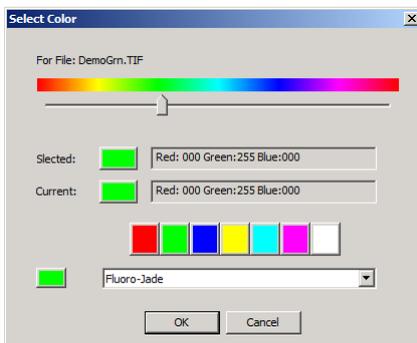

Directly select a color by:

- 1) Dragging the **Color Value Slider** to the desired color on the **Color Bar** (range of 0 to 359°): or
- 2) Clicking the appropriate **Color** button on the **Color Palette**. The **Color Bar** represents **Color** degrees in the standard color wheel representation, and the **Color Palette** contains a subset of possible **Colors** (pure red, green, blue, cyan, magenta, and white): or
- 3) Select a dye in the dye list, and click the **Select Dye** button.

Note: You can navigate through the **Select Color** dialog box by using the **<Tab>** and **<Shift + Tab>** keys to cycle through the items in the dialog box.

Intensity is determined by the image the **Color** is applied to, and **saturation** will always be 100%.

The only exception to this is the "white" **Color**, which has by definition a saturation of 0%. Selecting a value greater than or equal to 360 (dragging the **Color Value** slider over to the far right) is a choice of white, which results in a non-tinted (gray) image.

The elements of the **Select Color** dialog box are:

Color Bar: The **Color Bar** shows a rainbow displaying the **Colors** underneath the appropriate point of the **Color Value** slider.

Note:

If **Color Value Slider** is the selected (highlighted) selected control, then you can also control the slider with keyboard strokes:

Use the left/right arrow <-> or <-> keys to move the slider incrementally.

Press the <Page Up> and <Page Down> to move the slider in “chunk” sections along the **Color Bar**.

Use the <Home> and <End> keys to move the slider to the beginning and end of the **Color Bar**.

Selected: This color block displays the new **Color** selected by the **Color Value** slider bar or the **Color Palette**.

Selected Color: The **Selected Color** text field displays the corresponding numeric **Color** value of the **Selected** color block described above. The **Selected Color** value is expressed as an **RGB** color, with the red value first, followed by the green and blue values. If the color corresponds to a dye color, the dye name will be displayed. If the dye color corresponds to a color palette color, the name of the color will also be displayed.

Current: The **Current** color block displays the starting **Color** of the channel.

Color Palette: The **Color Palette** proves an alternate method of selecting a **Color**. Clicking on an item in the **Color Palette** will set the **Color Value** slider to the correct value and update the displays in both the **Selected** color block and **Selected Color** text field.

For example, clicking on the white button on the **Color Palette** will set the **Color Value** slider to the right, change the **Selected** color block to white, and cause the **Selected Color** text box to display “White”.

Adjust: The controls in this section are used to adjust the color composite contribution from each image, or for the color composite preview itself. You can adjust the input image that is currently selected in the **Colors** list by clicking the selected button. When you want to adjust the color composite preview, or you want no image is selected in the **Colors** list, use the **All** button.

Registration: **Registration** allows you to correct for mis-registrations of images caused by filter-induced optical shifts. Use the directional keys to shift the selected image in the x- and/or y- direction, with respect to the rest of the input images. One click moves that channel one pixel in the direction indicated. Blank areas are filled with black. The key labeled "0,0" will re-center the selected image.

The registration shift for the selected channel or the overall image is displayed in the static text box below the directional keys.

Contrast: Use the **BCG sliders** along with the **Best Fit** and **Reset** buttons to adjust the contrast of the selected image or the composite.

BCG sliders: The **Brightness**, **Contrast**, and **Gamma** sliders are the similar to the **Contrast Enhancement** feature in ToupView. You can also adjust each channel individually or collectively by using the **Selected** or **All** radio buttons (to the left of the **BCG** sliders in the **Adjust** group box).

Best Fit: Click this button to perform a **Best-Fit** contrast stretch on the selected input, automatically setting the **Brightness** and **Contrast** settings accordingly (**Gamma** will be reset to 50.). This method is identical to the **Best Fit** contrast adjustment under the standard **Contrast Enhancement** feature of ToupView. **Best Fit** is useful for first pass adjustments.

Reset: Pressing the **Reset** button automatically resets the **BCG** settings. If the **Adjust** radio button is set on **All**, the **BCG** settings for the overall color image will be reset.

Frame: Use the frame slider to adjust the frame of the selected image that is used for the preview, or the active frame of the composite. (Please see the discussion of **Color Composite** and **Image Sequences** that follows this section.)

Background: Select one of the available images to be the background color in this combo box. The rest of the image planes are subtracted from the background image — effectively punching "holes" into it and allowing the foreground colors to show without mixing with the background. This is most effective where the background covers large areas and the other images have smaller objects within the background objects.

Background allows the other inputs to be displayed “in front of” the background inputs, minimizing color mixing. For example, ToupView can display a red dot on a blue background without turning the dot to magenta.

Selecting **None** causes all images to be equally mixed.


New: Click on the **New** button to discard the current color composite image, and start over.

The **Source Images** combo box will reset to include all gray scale images (including floating point images). Selecting an image from the list will create a new composite the same size as the initial selection. As mentioned previously, the source images list will be filtered to contain only images of the same size as the initial selection.

13 Layer

13.1 About layer

ToupView **Layer** is like sheets of stacked acetate. You can see through transparent areas of a **Layer** to the **Layers** below. You add a **Layer** to position the image content on the **Layer**, like sliding a sheet of acetate in a stack. You can also hide/show a layer to make content invisible/visible.

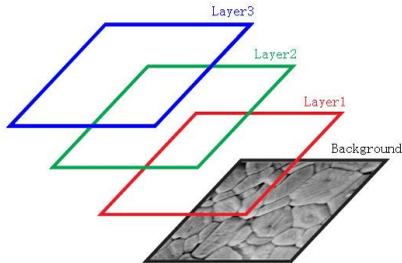

You use layers to perform tasks such as perform the measurement on the image without polluting the image and save it for the future adjustment for the further analysis, adding text to an image, or adding vector graphic shapes.

13.2 Organizing layers


A new image has a single **Layer** called **Background** layer. The number of additional layers you can add to an image is limited only by your computer's memory.

You work with layers in the **Layers Sidebar**. The **Layer Sidebar** help you organize and manage layers just like the **Layer** menu.

13.3 Layers for non-destructive measurement and label

Rather than edit image pixels directly, you can label the image with text and measure the image and leave the underlying pixels unchanged. Because of the non-destructively to images, you can later tweak or remove the label and the **Measurement Objects**.

13.4 Layer Sidebar

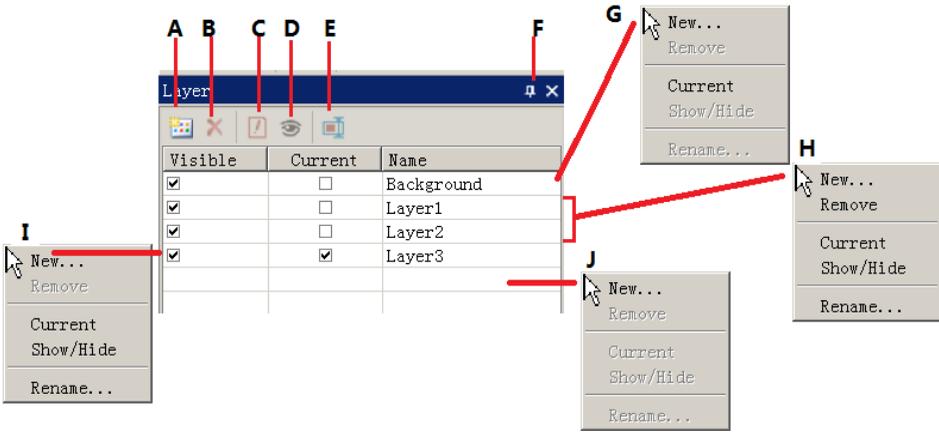

- A: Make a **New** layer;
- B: **Remove** a layer;
- C: Set the **Current** layer;
- D: **Show/Hide** a layer;
- E: **Rename** a layer;
- F: Auto hide the **Layer** Sidebar ;
- G: Right mouse button context menu for the **Background** layer;
- H: Right mouse button context menu for the **Current** not checked layer;
- I: Right mouse button context menu for the **Current** checked layer;
- J Right mouse button context menu in the blank area;

Note: a). Only one layer can be checked as **Current Layer**. If a layer's "Current" is checked, the **Visible** will be checked automatically; b). The **Current** and **Background** layers cannot be removed. Only the non-current and non- background layers can be removed; c). All the **Objects** on the **Current Layer** can be selected, edited or exported; d). The non-current **Layer** can be **Visible** (checked) or invisible (unchecked).

13.5 Layer menu and layer sidebar page context menu

The function of the **Layer** menu and **Layer Sidebar** right mouse button context menu are the same. They are:



13.6 New...

Setup a new layer. After the new layer is setup, it will be added to the end of the **Layer**

Sidebar's items and the **Visible** and **Current** will be checked.

13.7 Remove...

Remove the non-current **Layer**.

Note: the **Background** and **Current Layer** cannot be removed.

13.8 Current...

Set the selected layer as the **Current Layer**. User then can edit the **Object** on the **Current Layer** or added new **Object** on the **Current Layer**.

13.9 Show/Hide...

Set the non-current **Layer** visible or invisible. The **Current Layer** cannot be toggled between **Visible** and **invisible** state. It is always **Visible**.

13.10 Rename...

Rename a **Layer**.

13.11 Export to Image


Export all the layers' **Measurement Objects** to the image, this will pollute the image and cannot be recovered anymore.

13.12 Export to Microsoft Excel

Export the image plus all the **Objects** on the **Current Layer** to the excel file.

14 Measurements

The **Measurements** menu is mainly used for the image measurement application. With this menu items, you can measure the image with many geometrical shape at ease. ToupView use **Layer** technique to perform the measurement operations. This will never pollute the image pixels. The **Measurements** menus and its submenus are shown below.

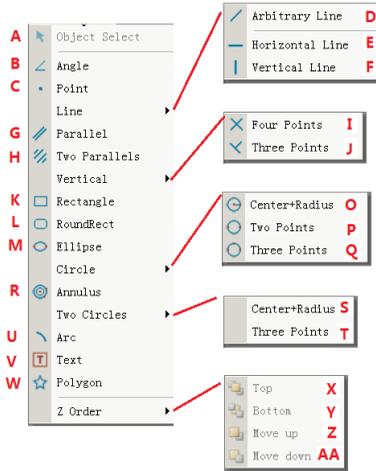

About the **Layer** technique, please check the **Layer** menu and **Layer Sidebar** in Sec.13.

About the **Measurements** setup, please check **Options>Measurement** and **Measurement Sidebar** in **View>Sidebar** menu in Sec.8.

About the **Measurements** menu's toolbar icons, please check ToupView's toolbar in Sec.4

14.1 Object Select




The **Measurements>Object Select** menu or the toolbar button will be enabled only when a new **Measurement** is made over the **Background Layer**.

After the **Measurement** is done on the specific layer, choose this menu to select the **Objects**.

1. The **Object** can be selected by clicking on it;
2. Select a group of **Objects** by including them in a rectangular area with **Object Select** command or by press down the **Shift** key and clicking the **Object** with left mouse button until all the desired **Objects** are selected.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding object parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.2 Angle

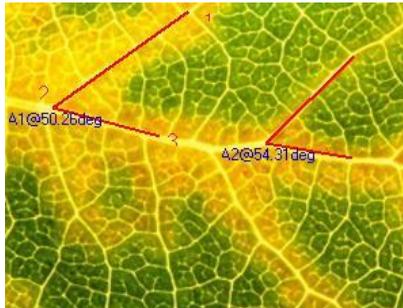

1. Move the mouse on the image to point 1, mark it by clicking the left mouse button;
2. Move the mouse to point 2, mark it by clicking the left mouse button;
3. Move the mouse to point 3, mark the final point by clicking the left mouse button. A label A1@50.26deg will be shown near point 2;

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the

selected **Object** type (The type could be the same or different).

14.3 Point



Move mouse to the point; click the left mouse to mark it. It will show the point **Label Pn** and its position x and y over the image.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).


14.4 Line

14.4.1 Line>Arbitrary Line



Draw **Arbitrary Line** between the two points on the specified layer.

1. Choose the **Measurements>Line> Arbitrary Line** command;
2. Move mouse to the 1st point; click the left mouse to mark it;
3. Move mouse to the 2nd point, click the left mouse to mark it again, a line with L1 and its length will be shown.

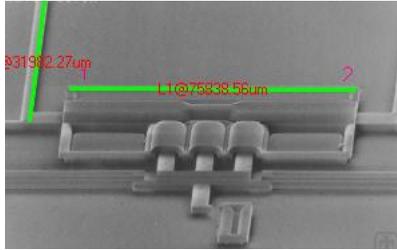

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.4.2 Line> Horizontal Line

Draw a **Horizontal Line** between the two points on the specified layer. When the first point is marked, the second point's y coordinate will be restricted to equal to the first point's y coordinate automatically.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.4.3 Line> Vertical Line

Draw a **Vertical Line** between the two points on the specified layer. When the first point is marked, the second point's x coordinate will be restricted to equal to the first point's x coordinate automatically.


Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated

automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.5 Parallel




1. Choose the **Measurements>Parallel** command;
2. Move mouse and click to mark the 1st point;
3. Move mouse and click to mark the 2nd point;
4. Move mouse and click to mark the 3rd point;
5. Move mouse again, one will find the 4th point is always restricted to parallel to the line connecting point 1 and point 2. Click to mark the 4th point. Two parallel lines will be labeled and numbered.

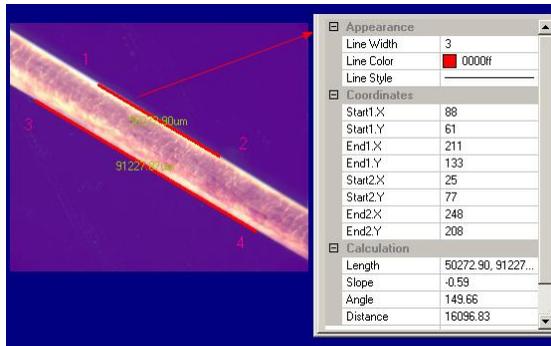

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.6 Vertical

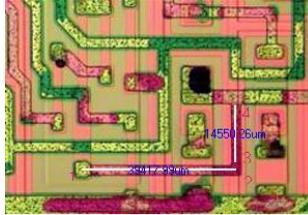

14.6.1 Vertical>Four Points.

1. Move mouse and click its left button to mark the 1st point.
2. Move mouse and click its left button to mark the 2nd point. A line (the 1st line) connect these two point be overlaid on the image
3. Move mouse and click its left button to mark the 3rd point.
4. Move mouse again, one will find the 4th point is always restricted to perpendicular to the line of point 1 and point 2. Click to mark the 4th point. Two vertical lines will be labeled and numbered.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).


14.6.2 Vertical>Three Points

1. Move mouse and click its left button to mark the 1st point;
2. Move mouse and click its left button to mark the 2nd point. A line (the 1st line) connecting these two points will be overlaid on the image;
3. Move mouse and click its left button to mark the 3rd point. The 2nd line will be overlaid on the image which is perpendicular to the 1st one.

Note: a) To check or modify the parameters of the selected **Object**, just select a

single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.7 Rectangle

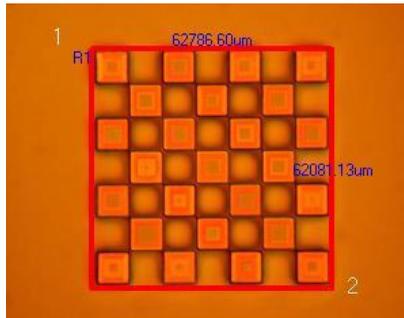

1. Move mouse to the 1st point; click its left mouse button to mark it;
2. Move mouse to the 2ed point; click its left mouse button to mark it. A **Rectangle** will be overlaid on the image according to these two diagonal points.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.8 RoundRect



1. Use **Arc** object to determine the round size. Its size is $21.39 \times 2 = 42.8$;
2. Choose the **Options>Measurement•••** command, click **Round Rectangle** to set the **Round x** and **y** to 43 and 43 respectively;

3. Move mouse to the 1st point; click its left mouse button to mark it;
4. Move mouse to the 2nd point, click its left mouse button to mark it, a round rectangle with Rr1 and its width and height the in the two dimensions will be displayed.




Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.9 Ellipse



1. Choose the **Measurements>Ellipse** command;
2. Move mouse to mark the 1st point;
3. Move mouse to mark the 2nd point;
4. If it is not superposing with the shape on the image, choose the **Measurements>Object Select** command to adjust the position to adjust the **Ellipse** in order to coincide with the image shape.

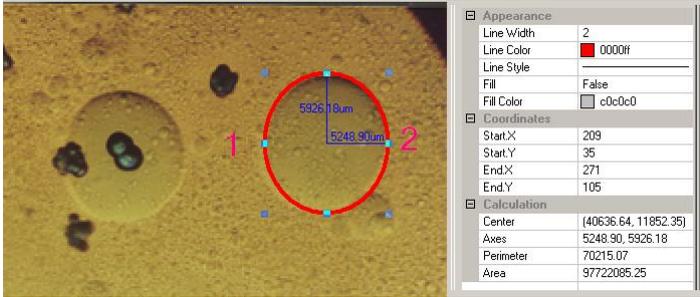

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.10 Circle

14.10.1 Circle>Center+Radius

Choose the **Measurements>Circle>Center+Radius** command to draw a circle with **Center+Radius** method on the specified **Layer**. Its radius is [C1@33583.88um](#).



Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.10.2 Circle>Two Points

Choose the **Measurements>Circle>Two Points** command to draw a circle with **Two Points** method on the specified layer.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance, Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).


14.10.3 Circle>Three Points

Choose the **Measurements>Circle>Three Points** command to draw a circle with **Three Points** method on the specified layer.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance, Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.11 Annulus

1. Find the **Annulus** center 1 and click mouse button to mark the **Annulus** center;
2. Move mouse to let the first circle superpose with the image circle, clicking the mouse left button;
3. Move mouse to let the second circle superpose the image circle, clicking the left mouse button. Two radiuses with number and the units will be displayed on the **Annulus**;

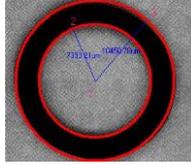

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.12 Two Circles

14.12.1 Two Circle>Center+Radius

Draw two circles with **Center+Radius** method on the specified layer. After the two circles are drawn, a line connecting these two circle centers will be drawn.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).


14.12.2 Two Circle>Three Points

Draw two circles with **Three Points** method on the specified layer. After the two circles are drawn, a line connecting these two circle centers will be drawn.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single

selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.13 Arc


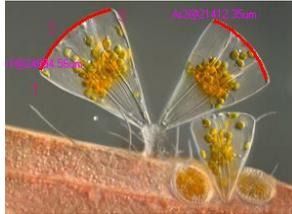


1. Move mouse to mark the 1st point;
2. Move mouse to mark the 2nd point;
3. Move mouse to mark the 3rd point, then an **Arc** with **Label** and its **Length** will be drawn connecting these three points.

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.14 Text


1. Choose the **Measurements>Text** command;
2. Move mouse to mark the 1st point;
3. Move mouse to mark the 2nd point, a rectangle with dash line restrict the **Text** window size. After the mouse button is released, a dialog called **Text** will bring up for you to enter the **Text** and define the **Text**, **Frame** styles and adjust the **Text** positions.
4. Enter the text and click the right mouse button to end the **Text** object.

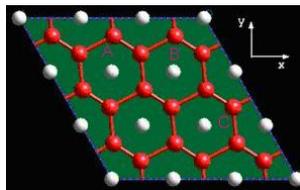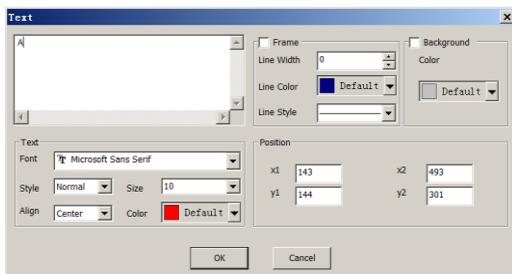

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.15 Polygon

1. Choose the **Measurements>Polygon** command;
2. Move mouse to mark the 1st point;
3. Move mouse to mark the 2nd point.
4. ●●●;
5. Move mouse to mark the nth point;
6. Repeat step1 to n to draw the other Polygon
7. Click the right mouse button to end the **Polygon** command

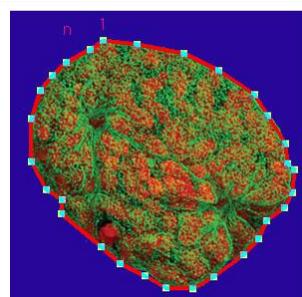

Note: a) To check or modify the parameters of the selected **Object**, just select a single **Measurement Object** and the **Measurement Sidebar** will be activated automatically. Here you can edit the corresponding **Object** parameters to modify its properties; b) The **Appearance**, **Coordinates** on **Measurement Sidebar** will be enabled only when a single **Object** is selected. The **Calculation** will be effective for a single selected **Object** or multiple selected **Objects**. Its calculated items will depend on the selected **Object** type (The type could be the same or different).

14.16 Z Order

Help Manual

Top	Modify the selected Object 's relative position to the uppermost place of all Objects on the Current layer
Bottom	Modify the selected Object 's relative position to the lowest place of all Objects on the Current layer.
Move Up	Modify the selected Object 's relative position to the higher place of the two on the Current layer.
Move Down	Modify the selected Object 's relative position to the lower place of the two on the Current layer


15 Options

15.1 Preferences●●●

There are 7 pages for the Preference dialog. They are File, Plugin, Print, Ruler and Grids, Cursor, Capture and Misc pages.


15.1.1 File

One can check a file Extension for the specified file Format and its Abbr.(abbreviation), to determine whether it will be displayed in the Browse window or not (The checked Format can be displayed in the Browse window).

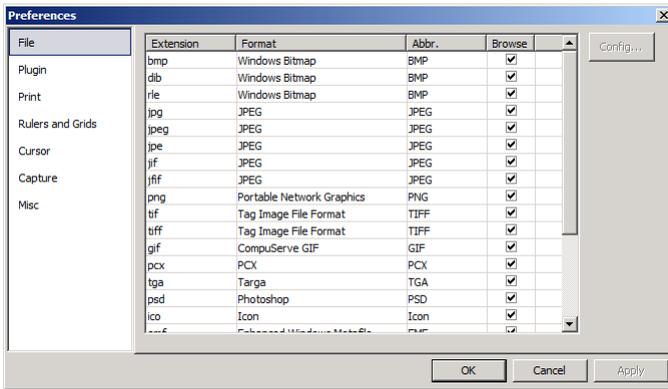


15.1.2 Plugin

The ToupView installed Plugins will be appended to the Plugin menu with the name specified in the plugin.XXX.cfg file under the installation directory. The current version includes 5 submenus as shown below:

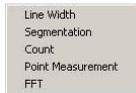

ToupView currently only provide Line Width, Segmentation, Count, Dimensioning, and FFT Plugins.


Every Plugin submenu can be Configured by choosing the Options>Preference●●● command and clicking Plugin page shown below:

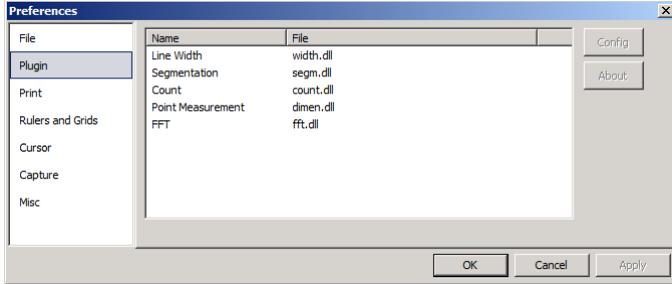

Click the item with the left mouse button to highlight it, the **Config** button will be enabled (This function is supported by the supplier, if this function is not provided; the **Config** button will be disabled). The same is true for the **About** button.

15.1.3 Print

One can set the **Page Header** and **Page Footer** in following format:

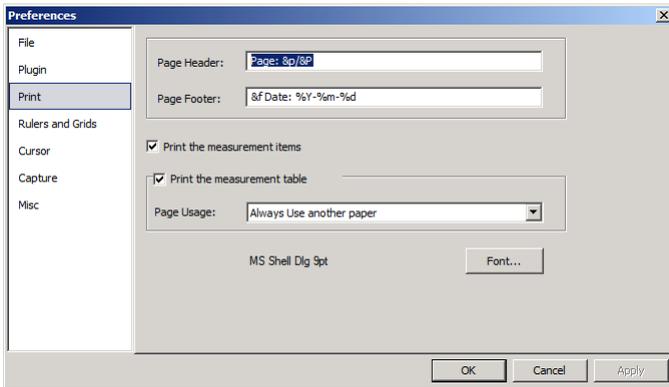

&f File Name

&p Current Page

&P Total Pages

%A Full weekday name

%b Abbreviated month name

%B Full month name

%c Date and time representation appropriate for locale

%d Day of month as decimal number (01 - 31)

%H Hour in 24-hour format (00 - 23)

Help Manual

- %I** Hour in 12-hour format (01 - 12)
- %j** Day of year as decimal number (001 - 366)
- %m** Month as decimal number (01 - 12)
- %M** Minute as decimal number (00 - 59)
- %p** Current local times A.M./P.M. indicator for 12-hour clock
- %S** Second as decimal number (00 - 59)
- %U** Week of year as decimal number, with Sunday as first day of week(00-53)
- %w** Weekday as decimal number (0 - 6; Sunday is 0)
- %W** Week of year as decimal number, with Monday as first day of week(00-53)
- %x** Date representation for current locale
- %X** Time representation for current locale
- %y** Year without century, as decimal number (00 - 99)
- %Y** Year with century, as decimal number
- %z %Z** Time-zone name or abbreviation; no characters if time zone is unknown
- %%** Percent sign

For example:

Choosing **&f Date:%Y-%m-%d** will print the file name and Date with yyyy-mm-dd format on the page header.

Choosing **Page:&p/&P** will print Page: 01/11 on the page footer(Suppose the current page is page 1 and the total pages is 11).

Print the measurement item: If there are measurement item overlaid on the image, check it will print these objects on the image. **Measurement Sheet** on the page;

Print the measurement table: If there are measurement item overlaid on the image, check it will print these objects' **Measurement Sheet** on the page with;

Always use another paper: Print the **Measurement Sheet** data on a separate page;

Auto: Print the **Measurement Sheet** data just behind the image (if the image and data can be arrange in a single page) of on a separate page (if not);


Do not use another paper: Print the **Measurement Sheet** data just behind the image page;

Font: Choose the **Font** for the data.

15.1.4 Rulers and Grids

Choosing the **View>Ruler and Grid>Setting...** command, or choosing the **Preference>...** command and clicking the **Rulers and Grids** page realize the same function.

Choosing the **Preference>...** command and clicking the **Rulers and Grids** page can set **Ruler Color**, **Cursor Color** of the **Horizontal Ruler** and **Vertical Ruler** for the video or image windows. The **Grid Style**, **Line Style**, and **Line Color** for the **Grid** overlaid on the video and image window can also be set.

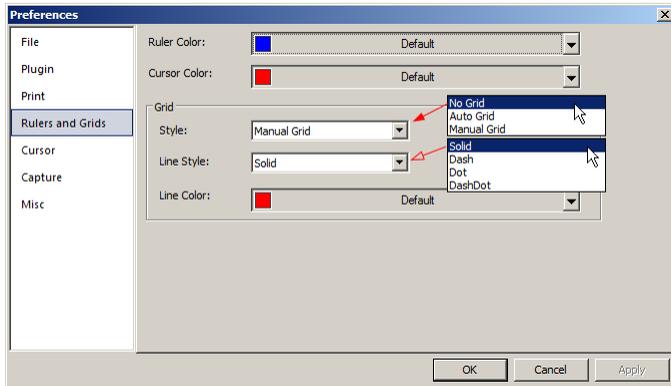

Ruler Color: The **Horizontal Ruler** color and **Vertical Ruler** color; Default is blue (0,0,255);

Cursor Color: The **Cursor** color; When you are moving your mouse over the video/image window, there will be **Horizontal Cursor** and **Vertical Cursor** move on the **Horizontal Ruler** and **Vertical Ruler** moving at the same time indicating the current **Cursor** position in both directions. Default is Red (255,0,0);


Grid Style: The **Grid Style** can be **No Grid**, **Auto Grid** or **Manual Grid**. Default is **No Grid**;

Grid Line Style: The **Line Style** for the grid can be **Solid**, **Dash**, **Dot** or **DashDot** et al. Default is **Solid**;

Grid Line Color: The color of the grid line. Default is Red (255,0,0).

15.1.5 Cursor

This command will set the mouse cursor for the video and image window operations.

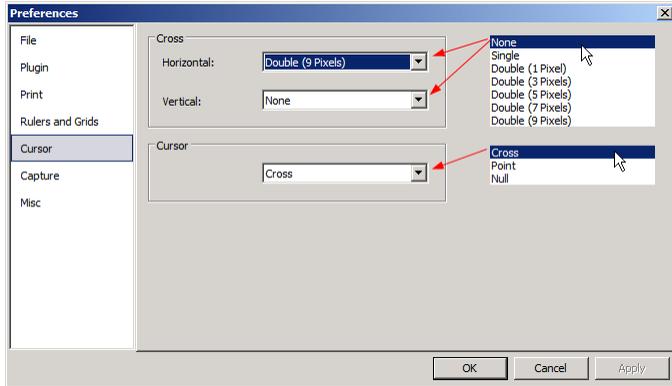

Select the **Horizontal** cross in:

None (window default), Single (single line), Double(1 Pixel), Double(3 Pixels), Double(5 Pixels), Double(7 Pixels), and Double(9 Pixels) formats.


Select the **Vertical** cursor in:

None (window default), Single (single line), Double(1 Pixel), Double(3 Pixels), Double(5 Pixels), Double(7 Pixels), and Double(9 Pixels) formats.

Single means single line. Double means two parallel lines. 1 Pixel means the line space between the two lines is 1 pixel in distance. The other sizes also have this meaning.

Select the cursor shape from **Cross**, **Point**, and **Null**.


In the figure below, a cursor with **Double horizontal and vertical lines with 9 Pixels cross** (for the alignment application, the longer one) and the **Cursor** as a **Cross** was defined

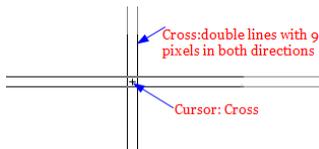

15.1.6 Capture

If one wish to capture an image with many **Measurement** objects overlaid on the video window, the following settings should be defined first.

1. Choosing the **Options>Preferences** command, a dialog called **Preferences** will be brought up as shown below;

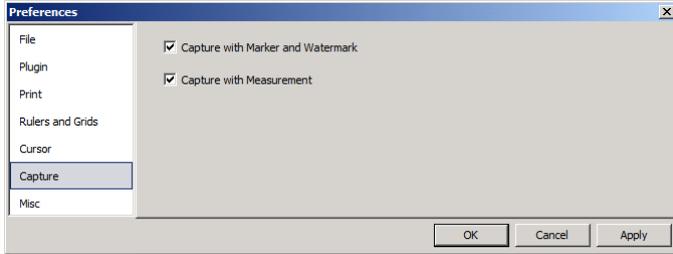

2. Clicking the **Capture** page and then checking **Capture with Marker and Watermark** or **Capture with Measurement**;

3. Click **Ok** to end the setup, or **Apply** to apply the setup, or **Cancel** to cancel the setup.

After the above steps are finished, if there are **Marker**, **Watermark** or **Measurement** objects overlaid above the video, choosing the **Capture>Capture Image** or clicking



on the **Camera Sidebar** will capture the image with **Marker**, **Watermark** or **Measurement** objects overlaid on the image.


15.1.7 Misc

The **Misc** page is mainly used for the control of the ToupView UI. It mainly includes:

1. ToupView' special file format warning information;
2. Sidebars;
3. UI Style;
4. ToupView Windows Layout;
4. Language;

5. Privacy;

settings et al.

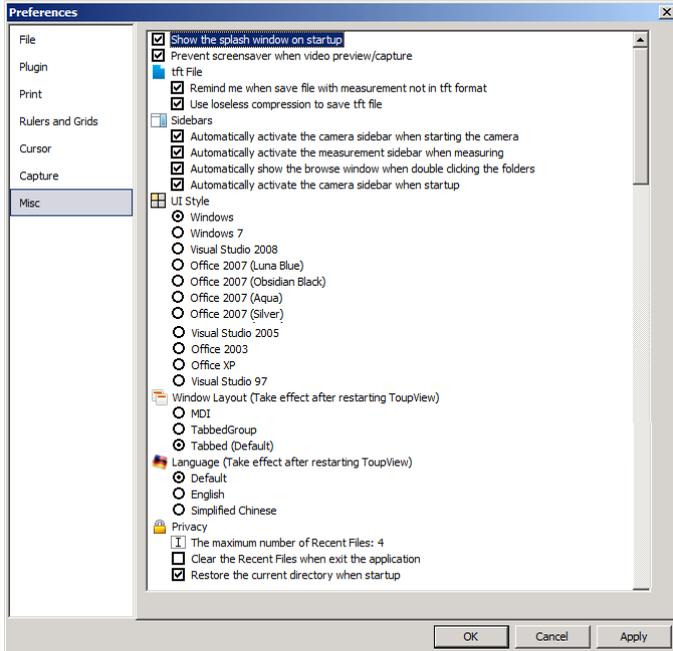


15.2 Measurements...

This sheet has many pages. They are:

15.2.1 General>General

The **General** page is shown below:

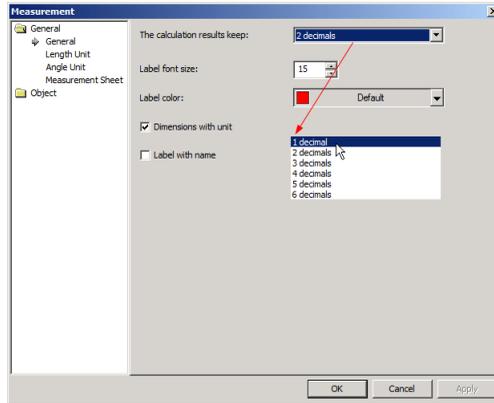


The calculation results keep: The combo box allows you to set the calculation results accuracy among 1 decimal and 6 decimals. Default is **2 decimals**;

Label font size: The **Label** is used to display the **Measurement Objects** dimensions. the **Label font size** is among 5 and 15; Default is 15;

Label Color: The Label Font color for the **Measurement Objects** dimensions; Default is red;

Dimensions with unit: If this box is checked, the **Dimension** will be displayed together with the unit for the **Measurement Objects**. If unchecked, only the **Dimension** will be displayed with the **Measurement Objects**.

Label with name: Checking **Label with name** will add a **Prefix** before the **Measurement Objects** number. The **Prefix** for the different **Object** is specified in the **Object** page's different subpage, where there is an item called: **Label: Prefix before number**. Here, we show an example for **Line Object** with and without **Prefix** as below:

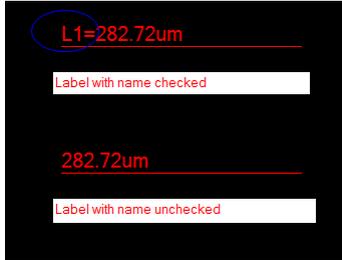

The defaults **Object** prefixes are shown below:

Object	Object Prefix
↗ Point	→ P
Line	→ L
Parallel	→ Pa
Two Parallels	→ Tp
Vertical	→ V
Angle	→ A
Rectangle	→ R
RoundRect	→ Rr
Circle	→ C
Arc	→ Ar
Annulus	→ An
Two Circles	→ Tc
Ellipse	→ E
Polygon	→ Py
Text	→

15.2.2 General>Length Unit

Length Unit page: One can select the **Length Unit** for the **Measurements** menu's **Objects** on the **Layer Measurement** operation.


Unit: The system unit defined by the ToupView. It cannot be deleted;

Current: The **Unit** selected;


Type: Type of unit. It can be **System** (Defined by ToupView) or **User** (Defined by the user) type;

Scale: Represents the ratio of "Meter by unit". For example, if the unit is μm , then "Meter by μm " should be 1000000, the **Scale** should be $1\text{e}+6$;

Help Manual

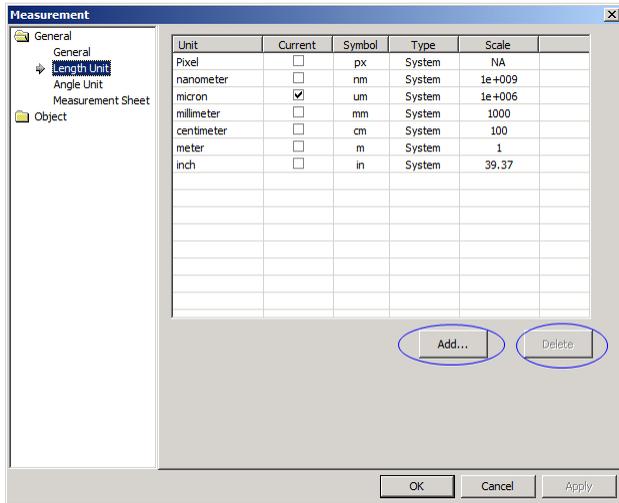


Add: User can also define his own Unit. Clicking **Add...** button and the **Add Unit** dialog will be brought up as below:

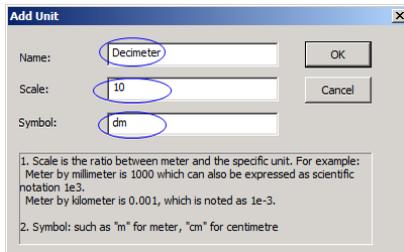

Here, one can enter the unit **Name** and **Symbol** in their fields. Here, we enter a name called **Decimeter** and its **Abbreviation** is **dm**, its **Scale** for **Meter** by **Decimeter** is 10. Click **OK** to end the **Add Unit** operation, or **Cancel** to cancel the **Add Unit** operation.

The final **Length Unit** list looks like the figure below:

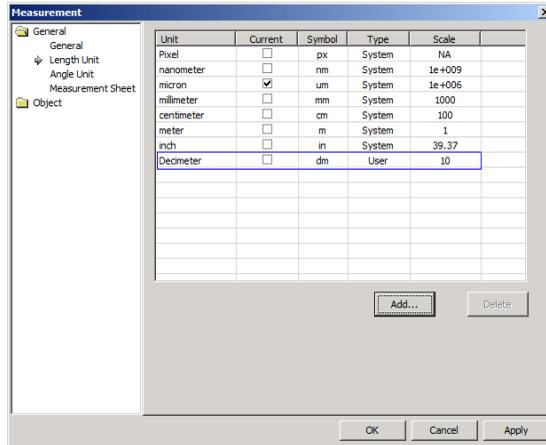

You can find the **Type** is now named **User**. This means that this **Unit** is not defined by the **System**, but by the **User**.

15.2.3 General>Angle Unit

On the **Angle Unit** page, one can select **Radian**, **Degree**, or **PI** as the **Angle Unit**.



15.2.4 General>Measurement Sheet


See the **View>Measurement Sheet** menu to understand its functions.

1. To modify the **Measurement Sheet**'s item order, click the item to highlight it, the **Up** or **Down** buttons will be enabled (If the item is in the third position, the **Up** button will be disabled, if the item is in the last position, the **Down** button will be disabled). Click the **Up** or **Down** buttons to modify the item position;

Note: The item **Index** and **Name** are always in grayed states, this means these two items cannot be modified (Both their hide/show states and their positions).

2. Checking/Unchecking the item will show/hide the item in the **Measurement Sheet**;

3. Clicking **Default** will return to the ToupView's default settings;
4. Item 1(**Index**) and item 2(**Name**) cannot be changed.

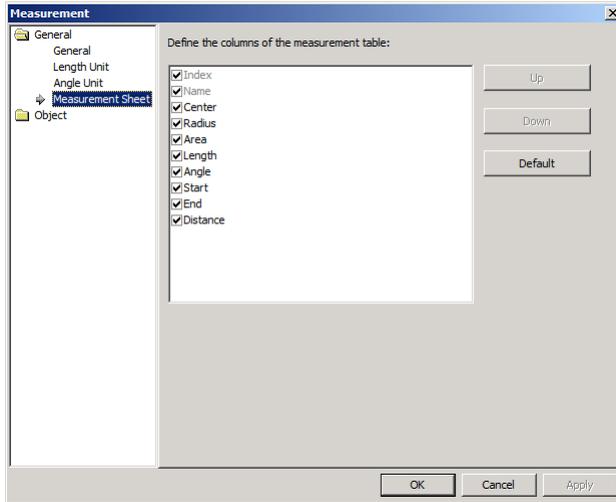

15.2.5 Object>Point

The **Object>Point's Appearance** describes the point **Size** and point **Color** and its **Prefix**.



Appearance: Size: The point will be represented by a cross. The size is the cross size in pixel. It is among 2-16. Default is 8;


Appearance: Color: Clicking the **Color** button's drop down arrow can select the point(or Cross) **Color**. Default is Blue (0,0,255);

Label: Prefix before the number: The **Label** is for all the **Layer Points**. Checking **Prefix before the number** or checking it on the **General>General** page's **Label with name** will

add a **Prefix** before the numeral for clarification. Unchecking it, the **Prefix Label** will never be displayed on any **Measurement Object**.

15.2.6 Object>Line

The **Line** page can set the **Line Width**, **Line Color**, **Line Style**, **Arrow** shape, and determine whether the objects are being labeled or not.

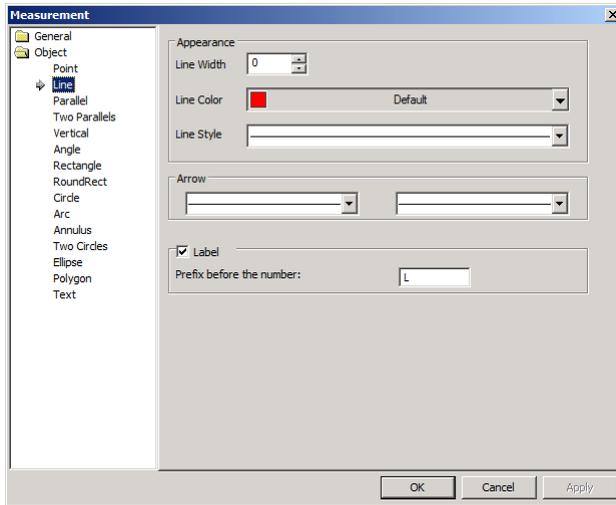

Appearance: Line Width: The **Line Width**, It is among 0 to 8. Default is 0(1 pixel);

Appearance: Line Color: Clicking the **Line Color** button's drop down arrow can select the **Line Color**. Default is Blue (255,0,0);

Appearance: Line Style: The **Line Styles** can be:



Appearance: Line Arrow: Different **Arrow** styles can be assigned separately to the **Line** start and end points. The **Arrow** styles can be:



Label: Prefix before the number: The **Label** is for the entire **Layer Lines**. Checking **Prefix before the number** or checking it on the **General>General** page's **Label with name** will add a **Prefix** before the numeral for clarification. Unchecking it, the **Prefix Label** will never be displayed on any **Measurement Object**.


6. If everything is ok, press **OK** to end the **Magnification** management dialog.

15.3 Calibrate●●●




The detailed calibration steps are as follows:

1. Run ToupView;
2. Connect the camera to the computer and microscope;
3. Start the camera (Here, it is **UCMOS00350KPA**);
4. Switch the microscope object to 10X and put the TS-M1 micrometer in the middle of the microscope field and try to find the ruler clearly. Set the **Unit** to **Pixel** and the video **Resolution** to the maximum one (640X480 for **UCMOS00350KPA**) and **Zoom** ratio to 100% (see );

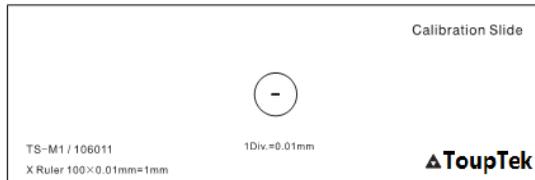

5. Choosing the **Options>Calibrate●●●** command or clicking “” on the toolbar, a red line with pixels number and 0.000um length is overlaid on the video window. At the same time, a dialog called **Calibrate** will be displayed over the video window;

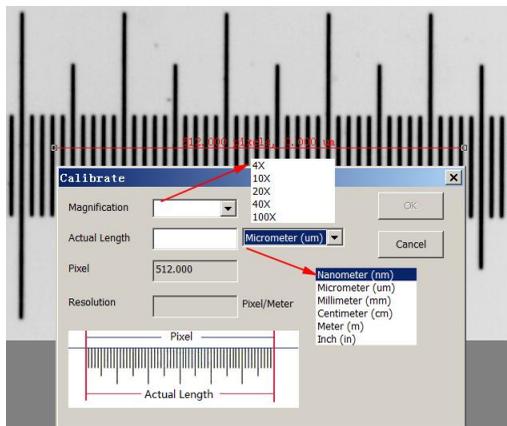


6. Try to align the two ends of the red line with the video ruler scale (Try to pull the

red line as long as possible to keep the calibration accuracy);

7. Enter or select the current microscope objective lens **Magnification** in the **Magnification** field. The current **Magnification** is 10X;


8. Read the overlaid (by the red line) micro-ruler actual length and fill it in the **Actual Length** field, the current **Actual Length** is 450um. This **Actual Length** will also be displayed in the middle of the red line just behind the pixel number. At the same time, the **Resolution** will be calculated and displayed simultaneously in the **Resolution** field.

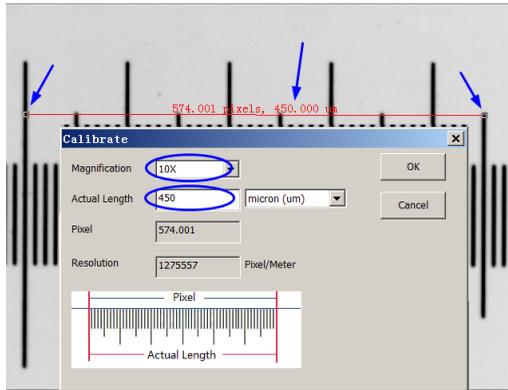


9. If everything is ok, click **Ok** to end the calibration. The **Magnification** 10X (This is the number you entered in the **Magnification** edit box) will be available in the

Magnification dropdown combobox on the video window toolbar 

10. If you wish to make the measurement under this **Magnification**, choose 10X in the above **Magnification** dropdown combobox first and then choose **Unit** in the **Unit** dropdown list box () which is just in the left of the **Magnification**'s dropdown combobox (User can also choose **Unit** through the **Option>Measurements...** menu, a dialog called **Measurement** will be brought up. Click **Length Unit** page to display the **Length Unit** page and checking **Unit** in the **Current** item.

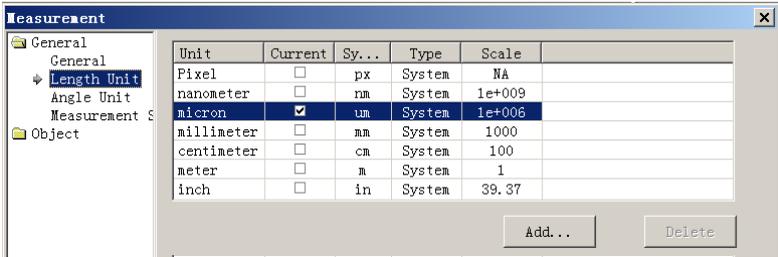

Now, you can measure the objects with the selected **Unit** at ease.

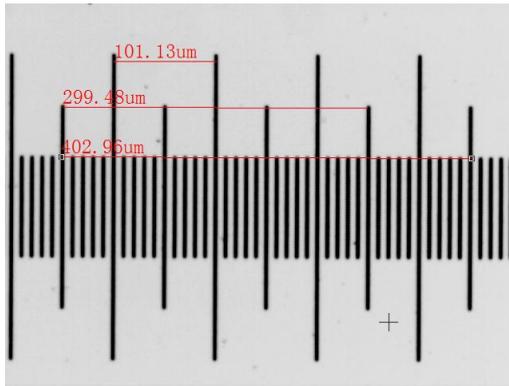


10. The selected **Resolution** can be saved for the future image **Measurement** operations. If the **Measurements** are performed on the video, the objects and **Resolution** can be saved in the image with **ToupView File Format (*.tft)** for the future applications;

11. The other microscope **Magnification** such as 4X, 40X, 100X can also be defined just as the above steps. Thus when you switch the microscope objective lens, you just need to select the **Magnification** () for example) on the toolbar to perform the **Measurement** operations.

15.4 Edit Dye List●●●

The new **Dye List** feature in ToupView allows you to select from a list of dyes, or add your custom dyes to the list. The list of dyes (**toupview.dye**) is stored under **C:\Program Files\ToupTek\ToupView** directory.

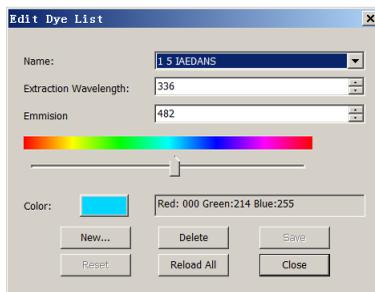

When you choose **Options>Edit Dye List** command, you will see the **Edit Dye List** dialog box:


The list includes any dye definitions found in the current location. When this feature is used to prompt for a specific dye, the list indicates the currently selected dye, or you may select a dye from the drop-down list.

Name: This list includes all dye definitions found in the current dye file. You may select a different dye from the drop-down list, and all the remaining controls will be updated to show the characteristics of the new dye.


Emissions Wavelength: This field displays the emissions wavelength for the dye you have chosen. You can adjust the emissions wavelength by entering a new value. The default color will change in response to changes in the emissions wavelength.

Excitation wavelength: The default wavelength for your dye is displayed here. You can adjust the wavelength by entering a new value.

Color: Use the slider to adjust the emissions wavelength for your dye, which also sets the default display color. (You can also use the **Color** button  to select a specific color using a standard Windows color selection dialog) The color will change in response to changes in the emissions wavelength. Here you can change the color associated with the selected dye, or define a custom color. The color default is linked to the selection of a wavelength for your dye. If you change the **Emissions Wavelength**, the color displayed will be the standard color for that wavelength. To set a custom color, first you should set the wavelength and then edit the color.

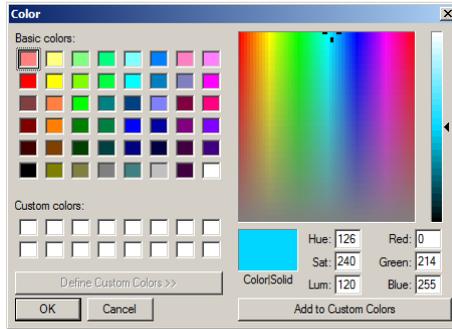

New: Clicking **New** will setup a new dye, this will invoke a dialog called **New Dye**. Input your **New Dye** name and click **Ok** to end the dialog and click **Cancel** to cancel the input. This will return to the **Edit Dye List** again and the new name will be the default on in the Name list box. Try to set the **Emissions wavelength**, **Excitation wavelength**, **Color** for your new dye.

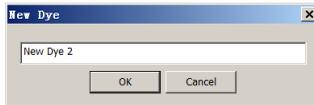


Save: Click **Save** on **Edit Dye List** to save your dye selection

Close: Click the **Close** button to end the **Edit Dye List** dialog:

15.5 Auto Correction●●●

Set the low and high ranges for the **Image>Adjust>Auto Level** and **Image>Adjust>Auto Contrast** menus. The default value is 0.5% for both, but ToupView suggests this value be smaller than 1%.

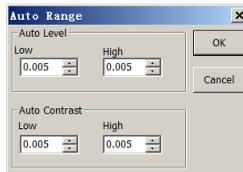


See **Image>Adjust>Auto Level** and **Image>Adjust>Auto Contrast** for details.


16 Windows

16.1 Close All


Choosing the **Closes All** command to close all of the pictures opened or captured inside ToupView frame. If you have made any modifications to the pictures or if you have captured some pictures from the camera, Choosing the **Close All** will let you finish the saving operations quickly.

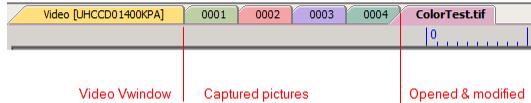

If the above image windows exist, choosing the **Windows>Close All** command will bring up a **Save Files** dialog as below:

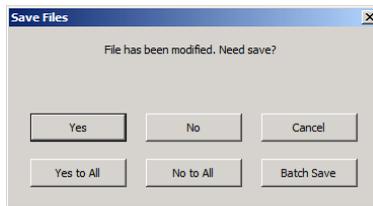

1. Clicking **Yes** on **Save Files** dialog will bring up a ToupView dialog that tell you the current active **Image has been modified**, And ToupView will ask you if you **Need save** it or not in a dialog as below:

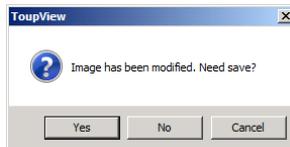

2. Clicking **Yes** on the **ToupView** dialog will bring up a **Save As** dialog to prompt you to enter the **File name** to save files, **No** will not save the current files or **Cancel** will cancel this file and let it still there and the process will step to the next image window. It will also bring up ToupView dialog for the new file. Repeat step 2 until all of the file saving operations have been finished;

3. Clicking **No** on **Save Files** dialog will bring up the **Save Files** dialog again if the

unprocessed file number is large then 1. If it is the final file, it will only bring up a **TouView** dialog as step 2. Click **Yes** to save the final file, or **No** to give up the file;

4. Clicking **Yes to All** on **Save Files** dialog will always bring up **Save As** dialog to let you enter the **File name** to save files one by one;

5. Clicking **No to All** on **Save Files** dialog will close all of the image windows without saving them;

6. Clicking **Batch Save** on **Save Files** dialog will transfer the save process to **File>Batch Save...** command. Please check the **File>Batch Save...** menu for details;

Note: If you have a lot of pictures opened and have made some modifications on them and want to close them quickly without saving anything, you may use this command and choose Step 5. All the images will be closed instantly, no warning will be given.

16.2 Windows...

Choosing the **Windows>Windows...** will open a dialog box that lets you manage the currently opened windows.

The dialog allows you to manage large lists of open windows by assigning them into groups. Let us say you have eight windows open, but want to close four of them scattered through the list.

1. Choose the **Windows>Windows...** command;
2. Select the windows you want to close. Hold **Ctrl** key to select more than one at a time. Here, 4 windows are selected;
3. Click the button that says **Close Window(s)**. Just those four will be closed, leaving the other four files or captures available for editing;

You can use this command to switch to the checked window on a list by hitting **Activate** (or "bring up" from an icon multiple windows).

17 Help

17.1 Help Contents

Choose the **Help>Help Contents** command to load ToupView help file. The help file is in Portable Document Format (PDF).

PDF is a file format used to represent documents in a manner independent of application software, hardware, and operating systems. Each PDF file encapsulates a complete description of a fixed-layout flat document, including the text, fonts, graphics, and other information needed to display it. In 1991, Adobe Systems co-founder John Warnock outlined a system called "Camelot" that evolved into PDF.

While the PDF specification has been available free of charge since at least 2001, PDF was originally a proprietary format controlled by Adobe. It was officially released as an open standard on July 1, 2008, and published by the International Organization for Standardization as ISO 32000-1:2008. In 2008, Adobe published a Public Patent License to ISO 32000-1 granting royalty-free rights for all patents owned by Adobe that are necessary to make, use, sell and distribute PDF compliant implementations.

If your computer still does not install the PDF reader, please try to download from:

<http://www.adobe.com/downloads/>

17.2 About

Display the related information about ToupView, including ToupView version, compatibility, built date and its developer's www etc.

Clicking on the **URL** link will direct you to the address of the camera supplier. If you have any problem with the camera or ToupView, please feel free to contact your supplier.

